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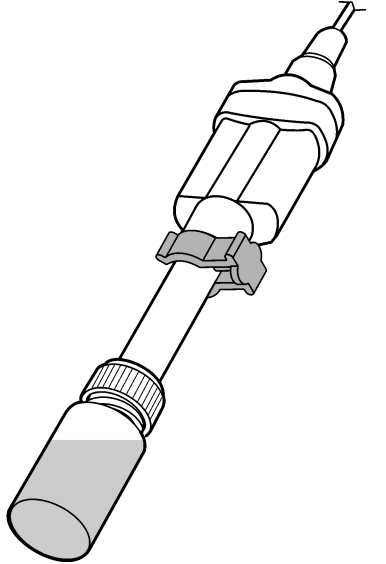
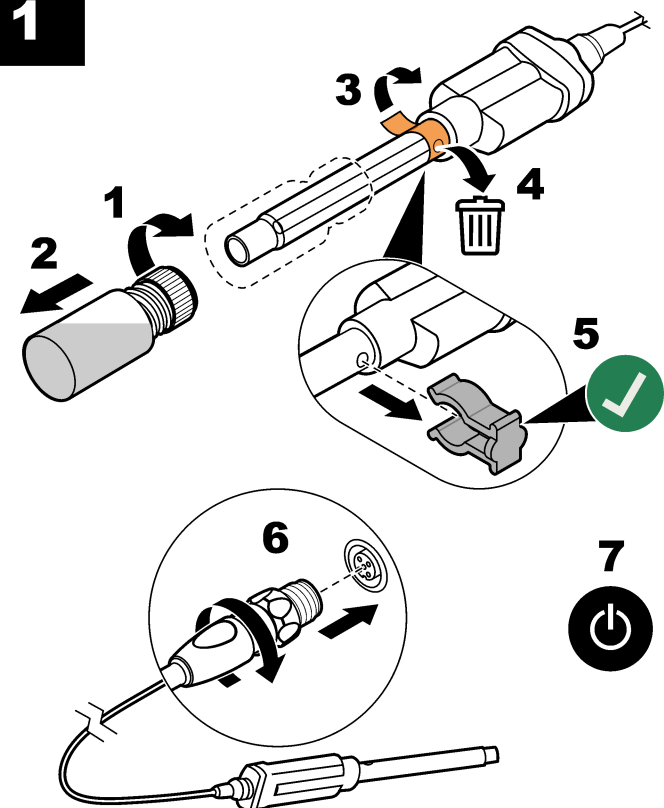
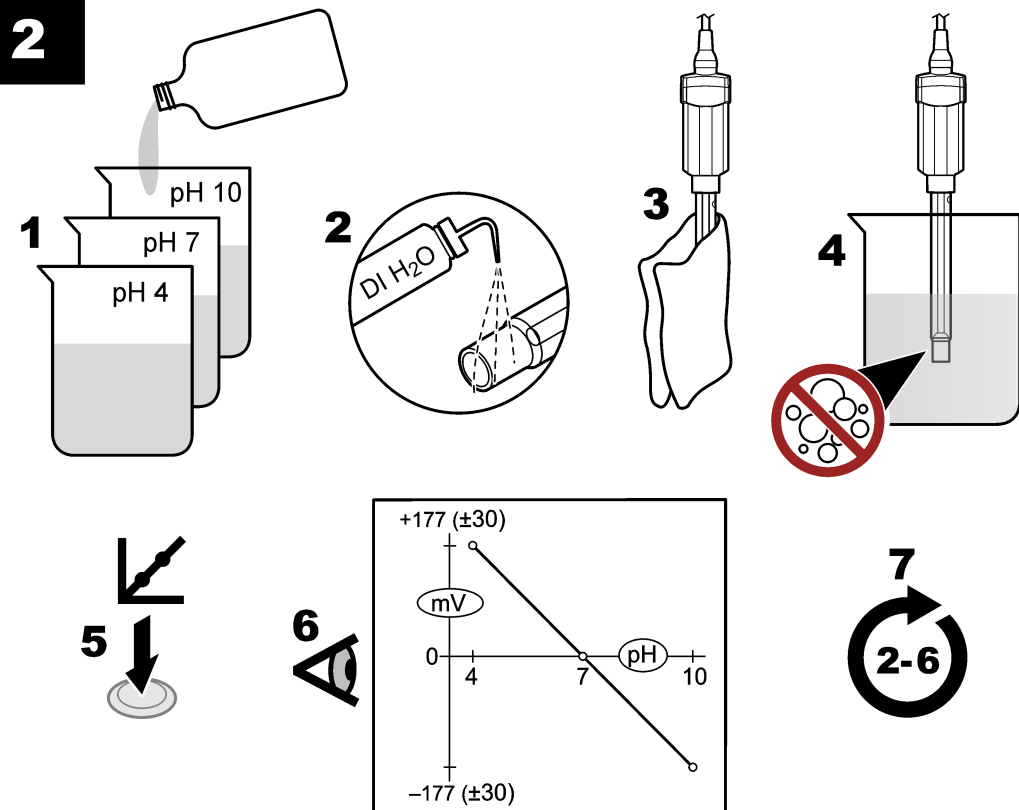
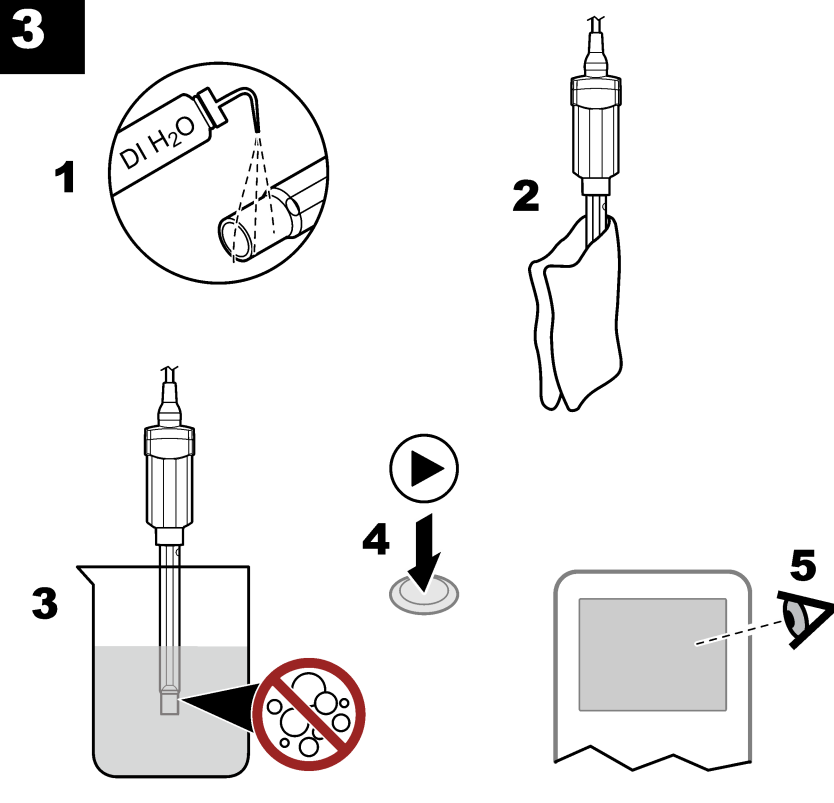
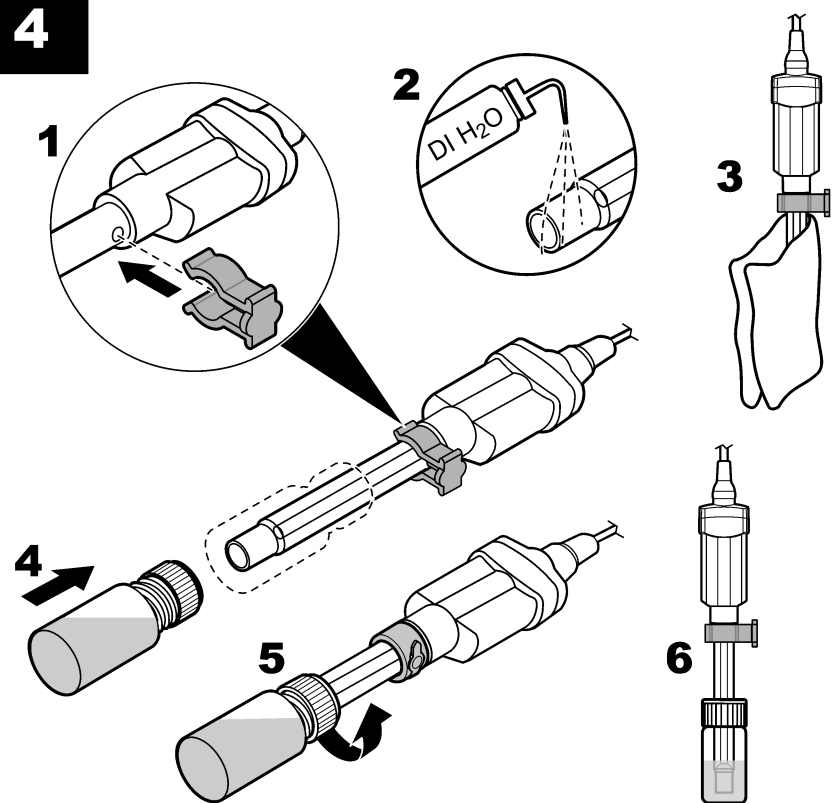
www.hach.com

**1****2****3****4****EN****PHC729****Introduction**

The probe is a pH electrode with refillable electrolyte, a glass body and an integrated temperature sensor. The probe measures pH on flat surfaces or where there is a minimal sample depth.

**Warning**

**⚠ Chemical exposure hazard. Obey laboratory safety procedures and wear all of the personal protective equipment appropriate to the chemicals that are handled. Refer to the current safety data sheets (MSDS/SDS) for safety protocols.**

**User manual**

A user manual with all instructions is on the manufacturer's website.

**Items to collect:**

Compatible meter, pH standard solutions, beakers, DI water, no-lint cloth, filling solution

**1 Connect the probe****2 Calibration**

*Air bubbles under the probe tip when submerged can cause slow stabilization or error in measurement. If bubbles are present, gently shake the probe until bubbles are removed.*

**3 Measurement****4 Storage****Technical information**

pH range	pH 0 to 12
Slope	-59 mV/pH (85 to 115% at 25 °C)
Temperature	-10 to 100 °C (14 to 212 °F)
Minimum sample depth	<1 mm (<0.04 in.)
Filling solution	Saturated KCl
Storage solution	Saturated KCl

**Maintenance and storage**

For maintenance and storage details, refer to the expanded version of this manual on the manufacturer's website.

**Warranty**

1 year for manufacturing faults. This warranty does not cover improper use or wear.

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