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



PHC735

PHC735

© Hach Company/Hach Lange GmbH, 2014, 2021. All rights reserved. Printed in U.S.A. 09/2021, Edition 2

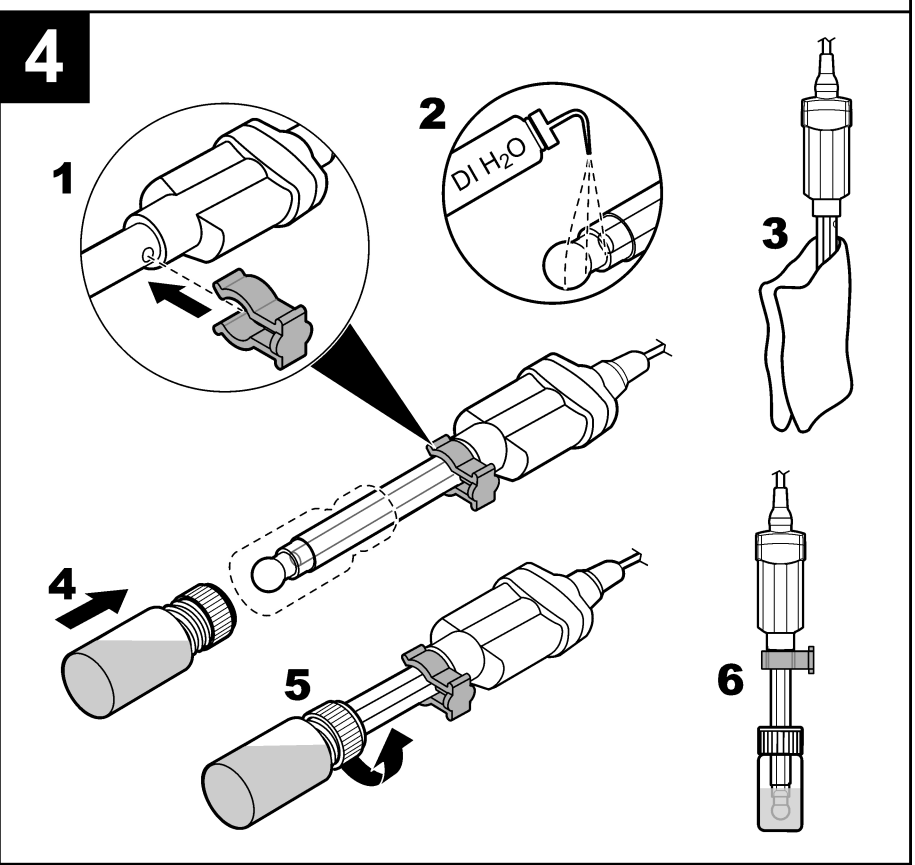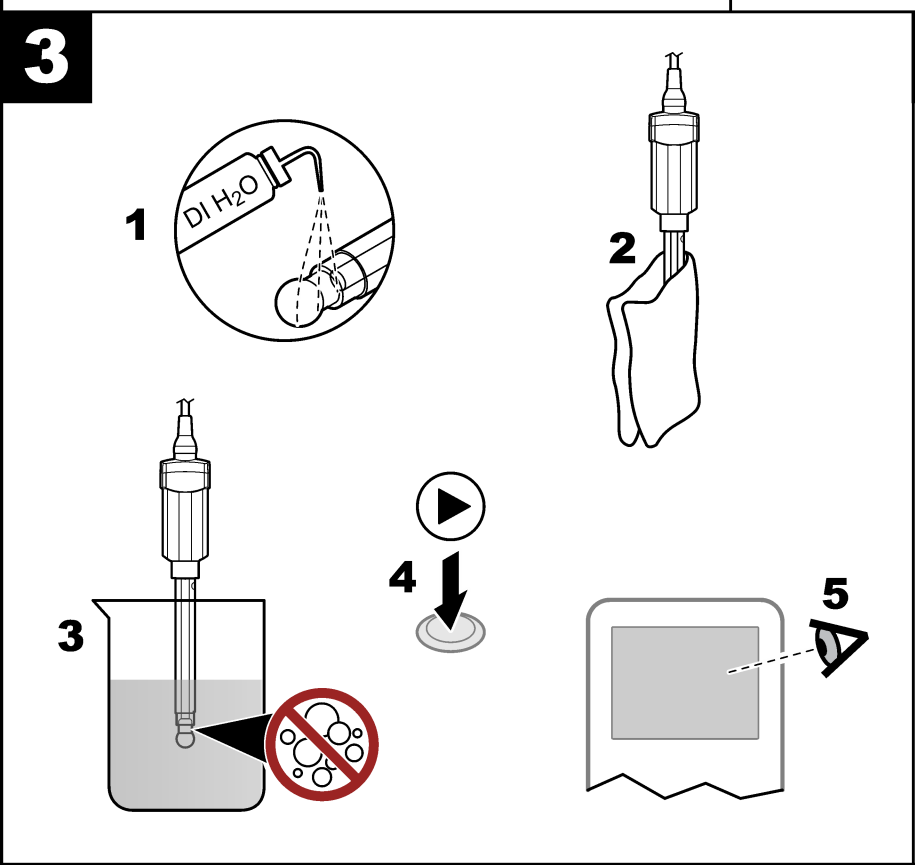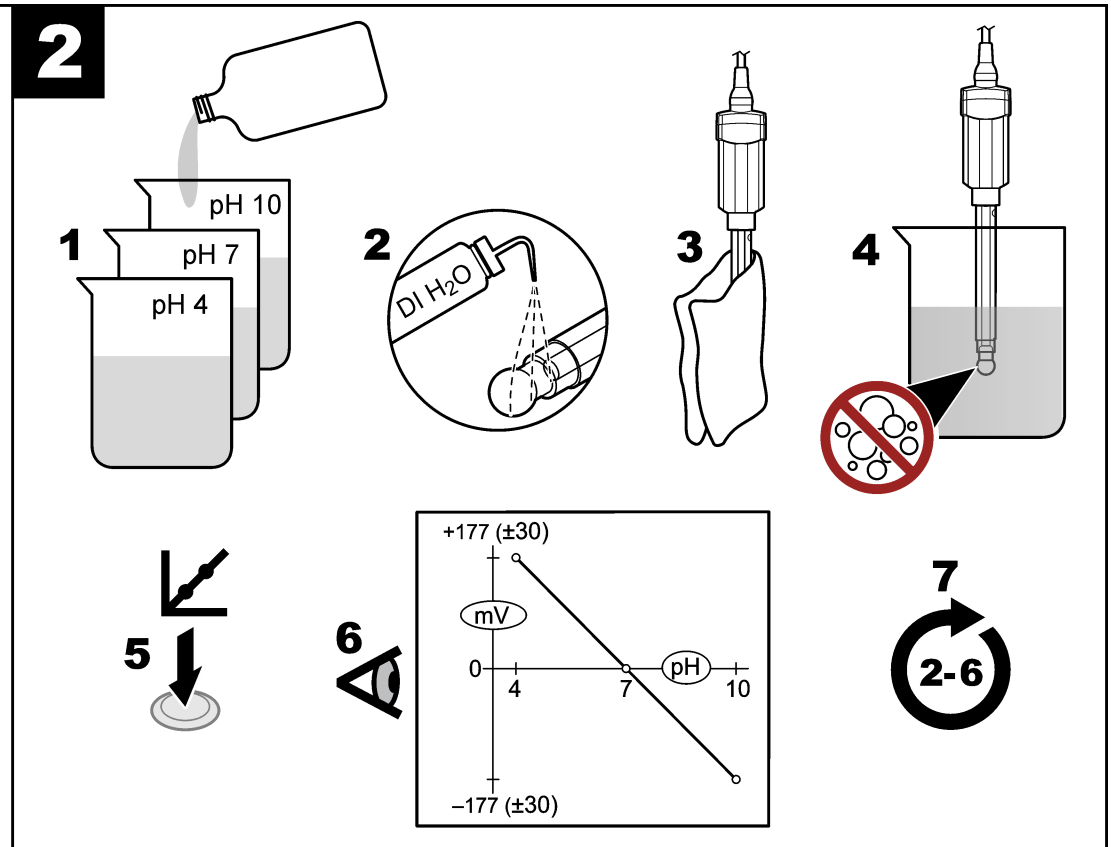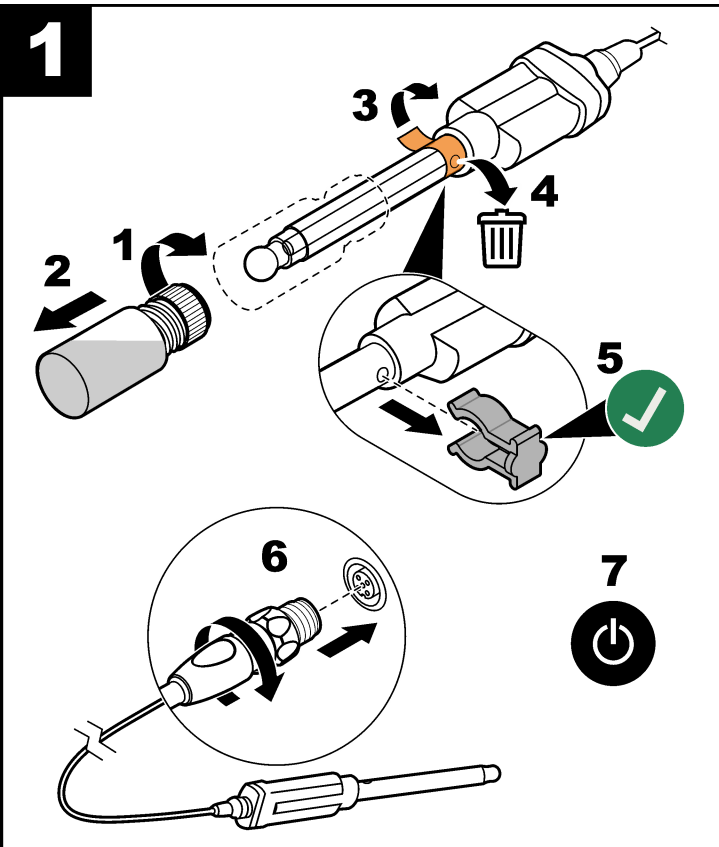PHC735 Introduction

The probe is a pH electrode with refillable electrolyte, a glass body and an integrated temperature sensor. The probe has a high-flow annular ring junction that measures pH in wastewater or "dirty" applications.

Warning
 ⚠ *Chemical exposure hazard. Obey laboratory safety procedures and wear all of the personal protective equipment appropriate to the chemicals that are handled. Refer to the current safety data sheets (MSDS/SDS) for safety protocols.*

User manual
 A user manual with all instructions is on the manufacturer's website.

Items to collect:
 Compatible meter, pH standard solutions, beakers, DI water, no-lint cloth, filling solution

- 1 Connect the probe**
- 2 Calibration**
Air bubbles under the probe tip when submerged can cause slow stabilization or error in measurement. If bubbles are present, gently shake the probe until bubbles are removed.
- 3 Measurement**
- 4 Storage**

Technical information

pH range	pH 0 to 14
Slope	-59 mV/pH (85 to 115% at 25 °C)
Temperature	-10 to 100 °C (14 to 212 °F)
Minimum sample depth	14 mm (0.55 in.)
Filling solution	Saturated KCl
Storage solution	Saturated KCl

Maintenance and storage
 For maintenance and storage details, refer to the expanded version of this manual on the manufacturer's website.

Warranty
 1 year for manufacturing faults. This warranty does not cover improper use or wear.

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