PHC745

Introdução

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



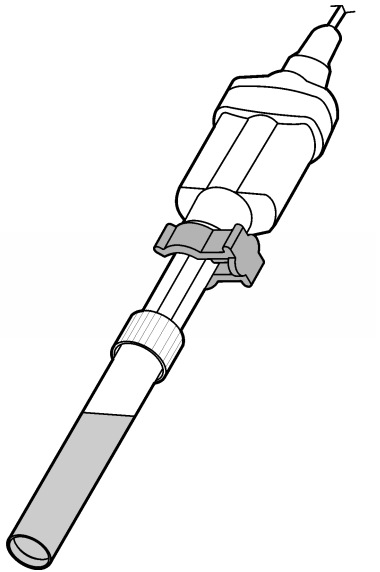
www.hach.com

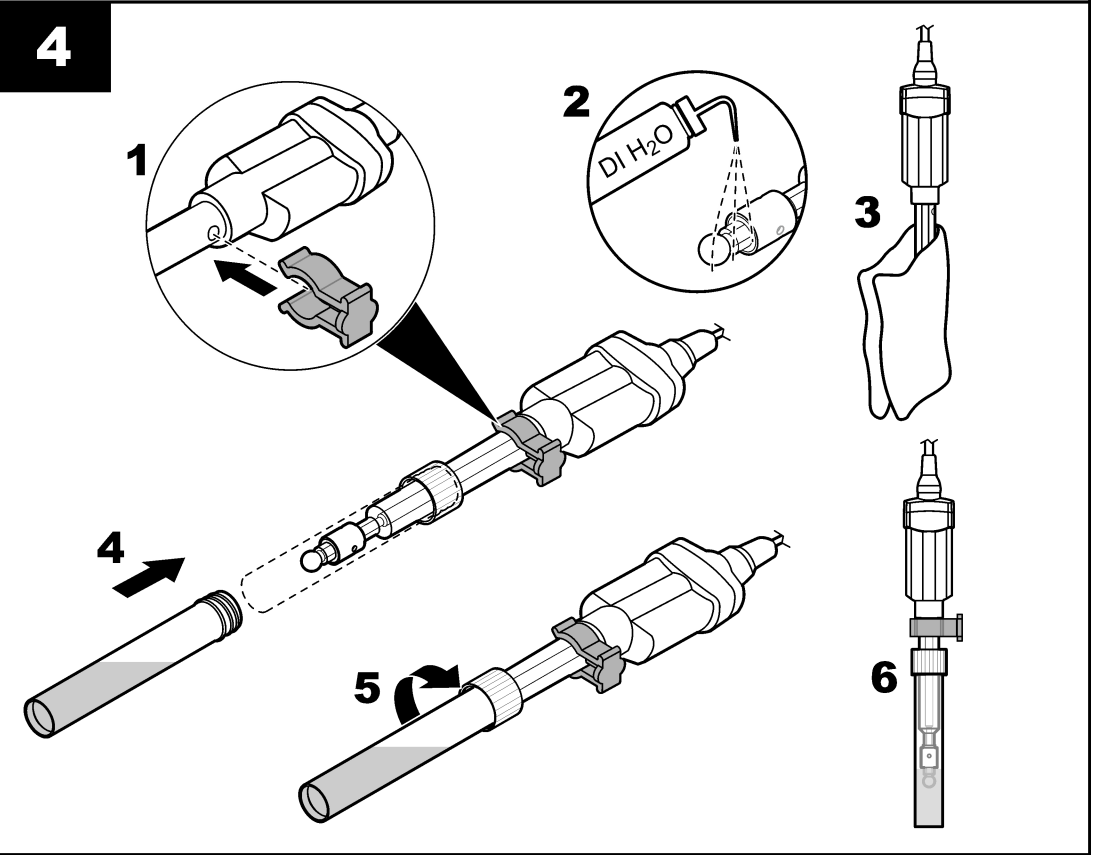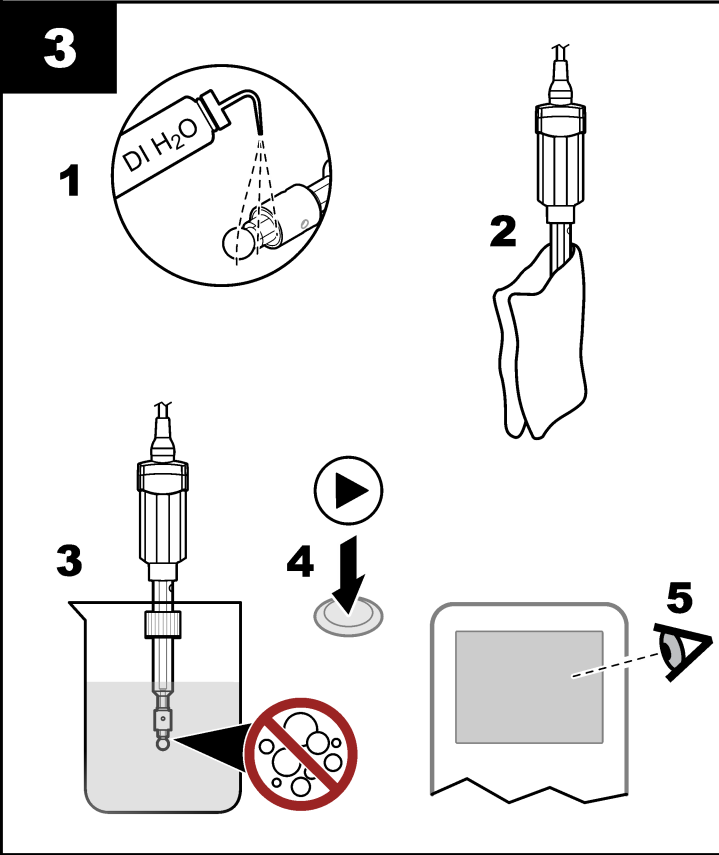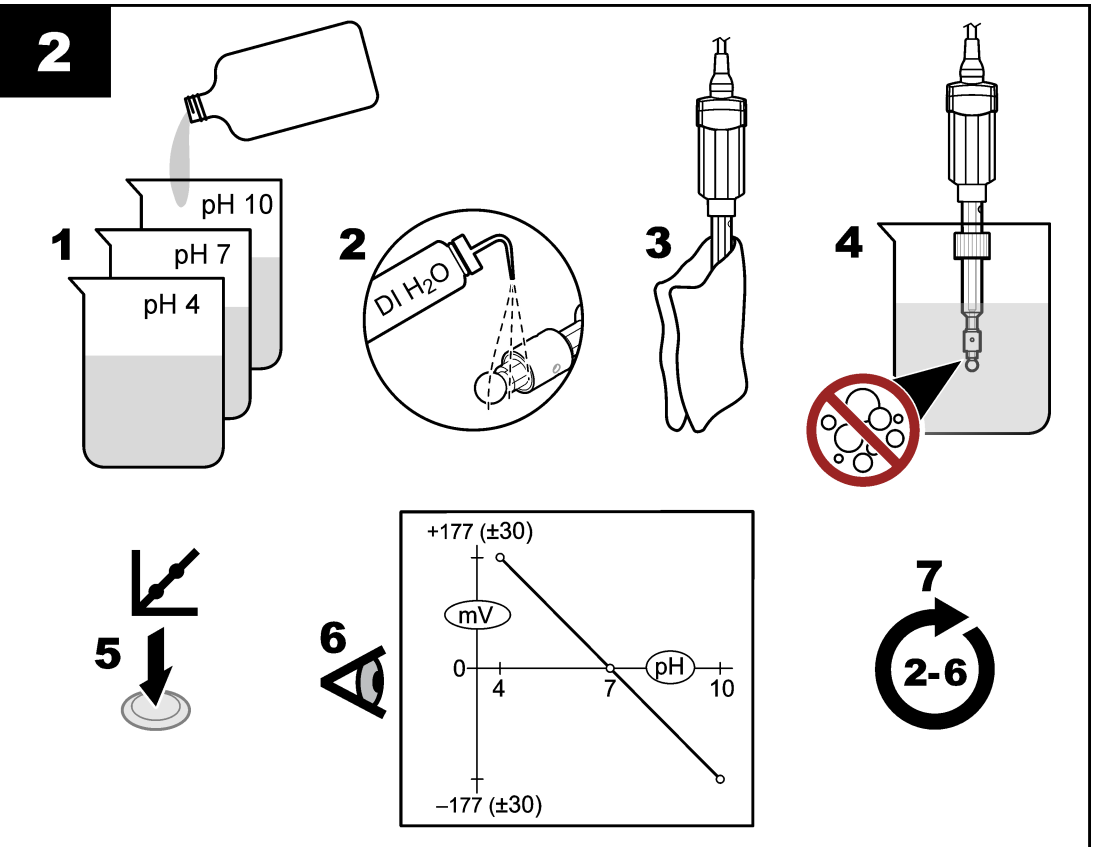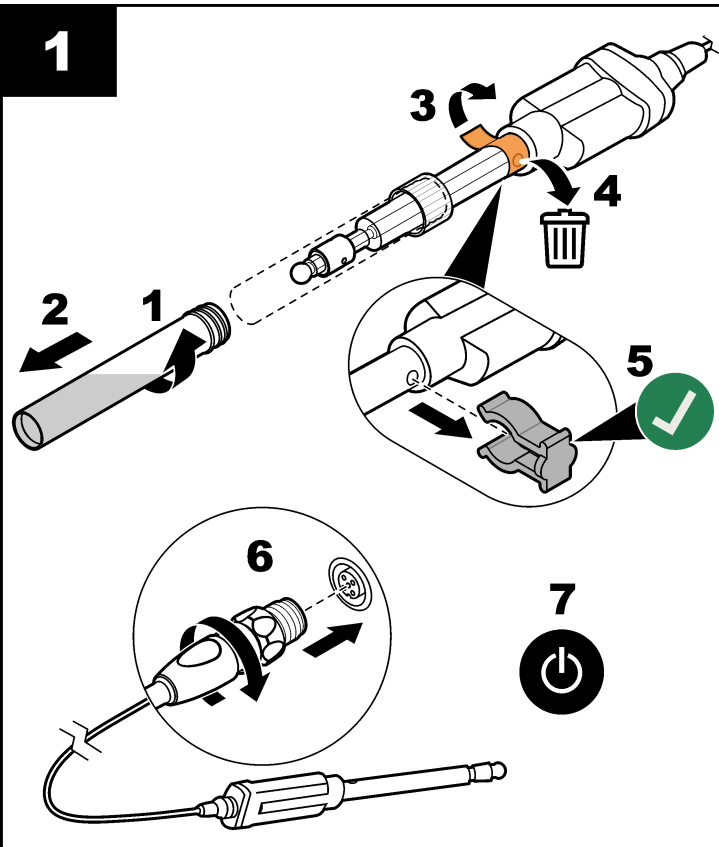PHC745 Introduction

The probe is a pH electrode with refillable electrolyte, a glass body and an integrated temperature sensor. The probe has a glass sleeve junction that measures pH in wastewater, emulsions and viscous liquids.

Warning
 ⚠ *Chemical exposure hazard. Obey laboratory safety procedures and wear all of the personal protective equipment appropriate to the chemicals that are handled. Refer to the current safety data sheets (MSDS/SDS) for safety protocols.*

User manual
 A user manual with all instructions is on the manufacturer's website.

Items to collect:
 Compatible meter, pH standard solutions, beakers, DI water, no-lint cloth, filling solution

1 Connect the probe
 The sleeve junction must be open to let the electrolyte touch the sample.

2 Calibration
Air bubbles under the probe tip when submerged can cause slow stabilization or error in measurement. If bubbles are present, gently shake the probe until bubbles are removed.

3 Measurement

4 Storage
 Always close the sleeve junction when the probe is not in use.

Technical information

pH range	pH 0 to 14
Slope	-59 mV/pH (85 to 115% at 25 °C)
Temperature	-10 to 100 °C (14 to 212 °F)
Minimum sample depth	18 mm (0.71 in.)
Filling solution	Saturated KCl
Storage solution	Saturated KCl

Maintenance and storage
 For maintenance and storage details, refer to the expanded version of this manual on the manufacturer's website.

Warranty
 1 year for manufacturing faults. This warranty does not cover improper use or wear.

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