Sonda PHC101

Introdução

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




PHC101

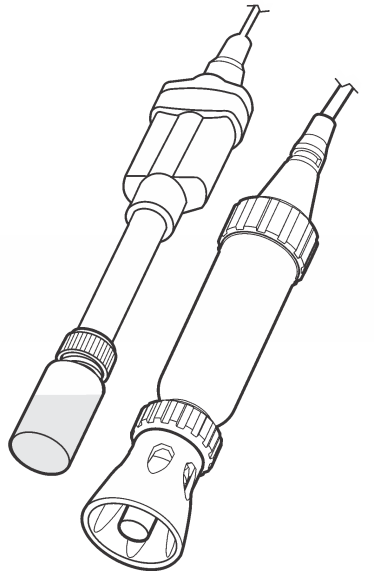www.hach.com

© Hach Company/Hach Lange GmbH, 2010, 2013, 2021. All rights reserved. Printed in U.S.A. 09/2021, Edition 5

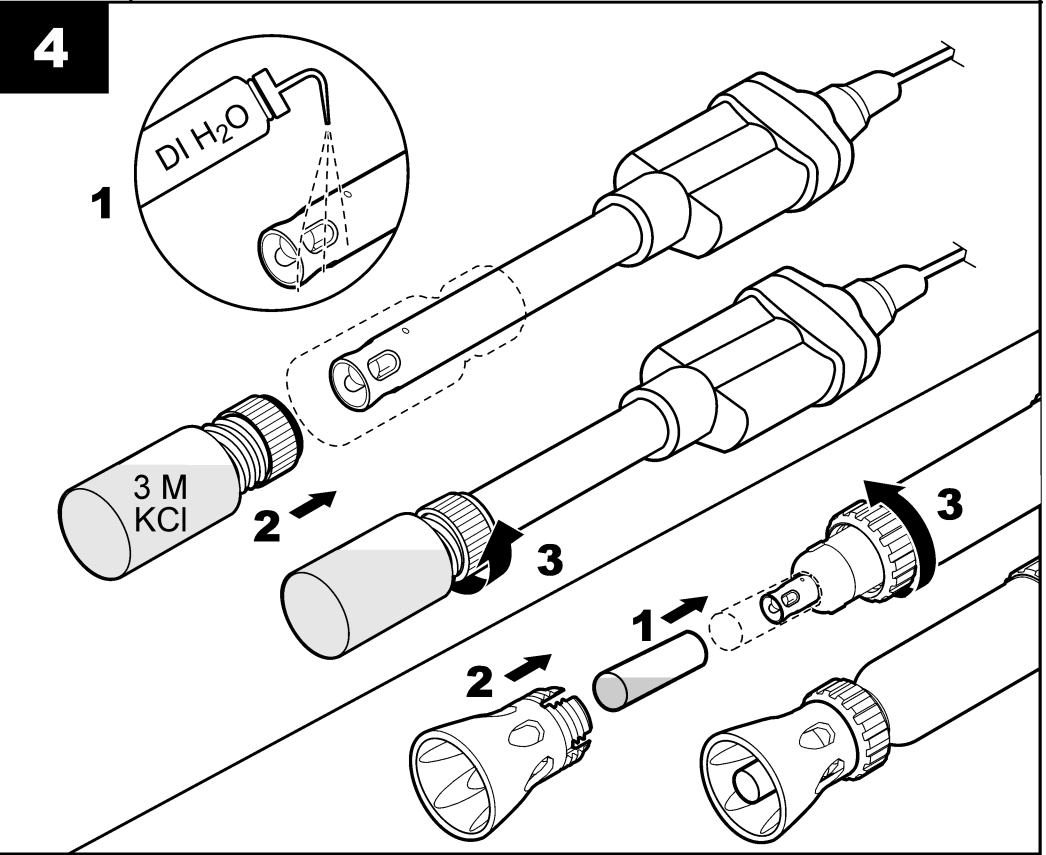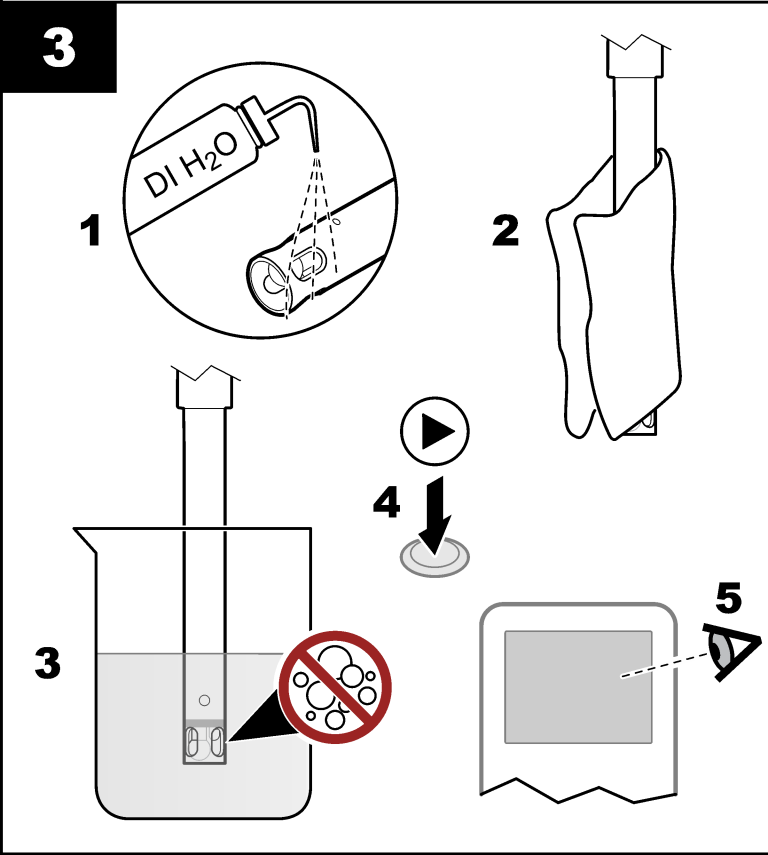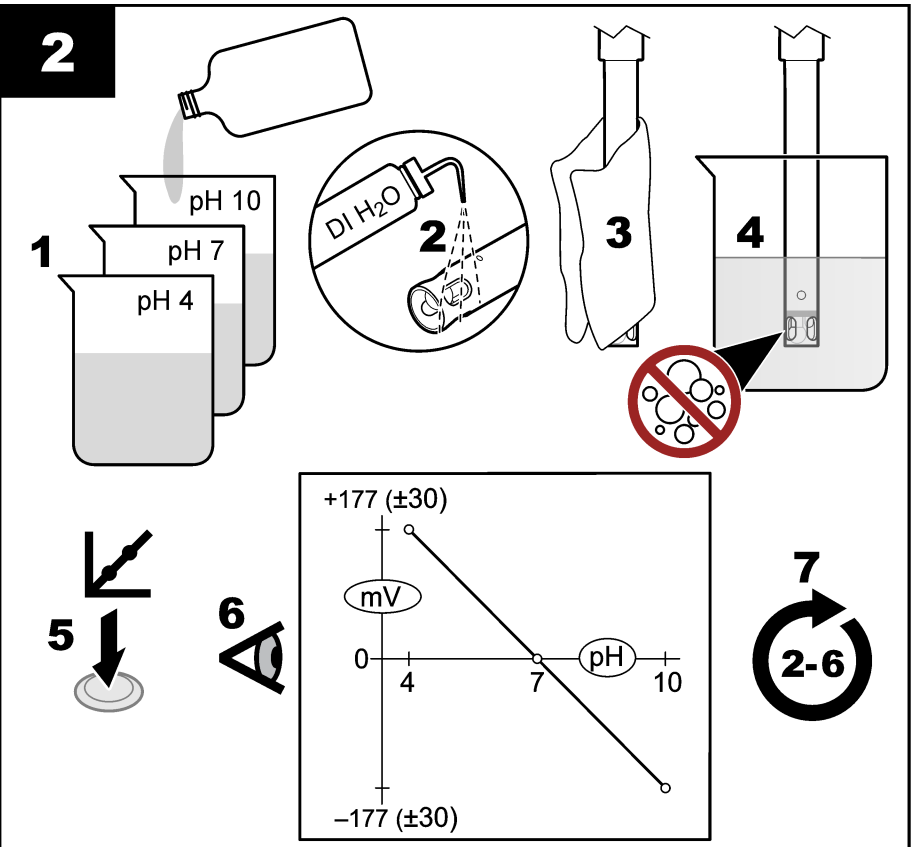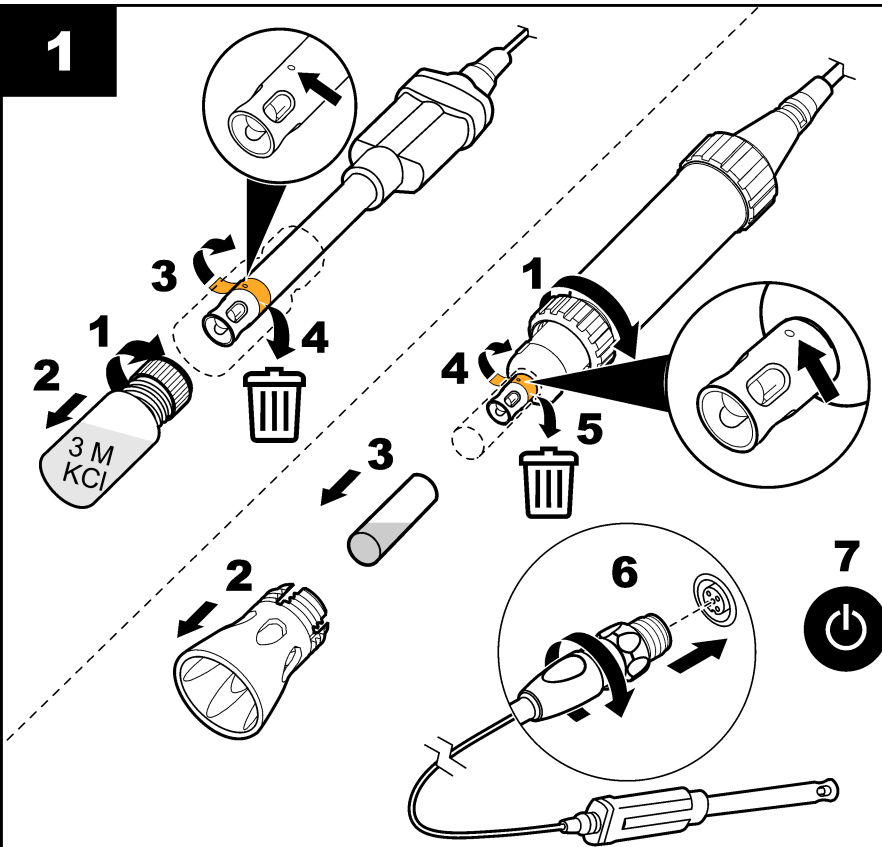EN

PHC101 probe Introduction

The probe is a non-refillable, gel-filled combination pH probe with a built-in temperature sensor for drinking water and general aqueous applications.

Warning
 ⚠ *Chemical exposure hazard. Obey laboratory safety procedures and wear all of the personal protective equipment appropriate to the chemicals that are handled. Refer to the current safety data sheets (MSDS/SDS) for safety protocols.*

User manual
 A user manual with all instructions is on the manufacturer's website.

Collect the following items:
 Compatible meter; pH standard solutions; beakers; DI water; lint-free cloth.

- 1 Connect the probe**
- 2 Calibration**
Air bubbles under the probe tip when submerged can cause slow stabilization or error in measurement. If bubbles are present, gently shake the probe until bubbles are removed.
- 3 Measurement (direct method)**
 ⚠ *Keep the shroud installed on the rugged probe during sample measurements to prevent damage to the sensor.*
- 4 Storage**

Technical information

pH range	pH 2 to 14
Slope	-59 mV/pH (90 to 110% at 25 °C (77 °F) per Nernstian theoretical value)
Operating temperature	0 to 50 °C (32 to 122 °F)
Temperature accuracy	±0.3 °C (±0.54 °F)
Minimum sample depth	20 mm (0.79 in.)
Storage solution	Hach pH electrode storage solution

Warranty

6 months on the probe. This warranty covers manufacturing defects, but not improper use or wear.

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