INFORMATION

EN

Safety information

In conformity with European local and national regulations, European electrical equipment users must now return old or end-of-life equipment to the manufacturer for disposal at no charge to the user.

The internal gel and solution do not contain dangerous compounds.

Prepare the electrode for use

Carefully remove the protection cap. Rinse the sensor with clean water. Make sure that the junction opening does not have an obstruction. If necessary, use a scraper to carefully remove the obstruction. Gently shake the sensor downward in the same manner as a clinical thermometer to remove any trapped air bubbles.

Clean the sensor

Clean the sensor when there is a buildup of deposits or when performance has decreased: Rinse with clean water. Soak the sensor for some minutes in a cleaning solution (Hach item no. S16M001 or 2965249). Then rinse with clean water. Soak the sensor for some minutes in a 3M KCl solution. Then rinse with clean water. Be sure to calibrate the sensor before use.

Storage

Keep the sensor in a vertical position for storage. Put the probe in the protector with KCl 3 M solution. Do not keep the sensor dry or in distilled water. The shelf life decreases during storage. Make sure to install the sensor before the use-by date shown on the box label.

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P.O. Box 389, Loveland, CO 80539-0389 U.S.A.
(800) 227-4224 (U.S.A. only)