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

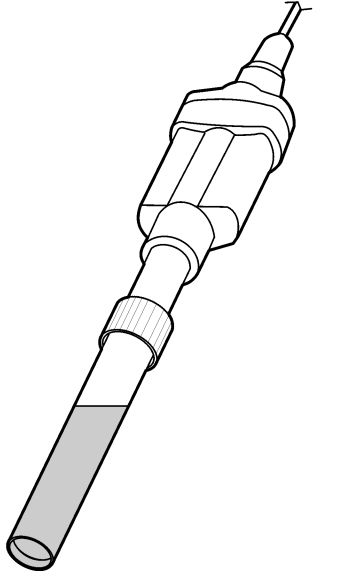
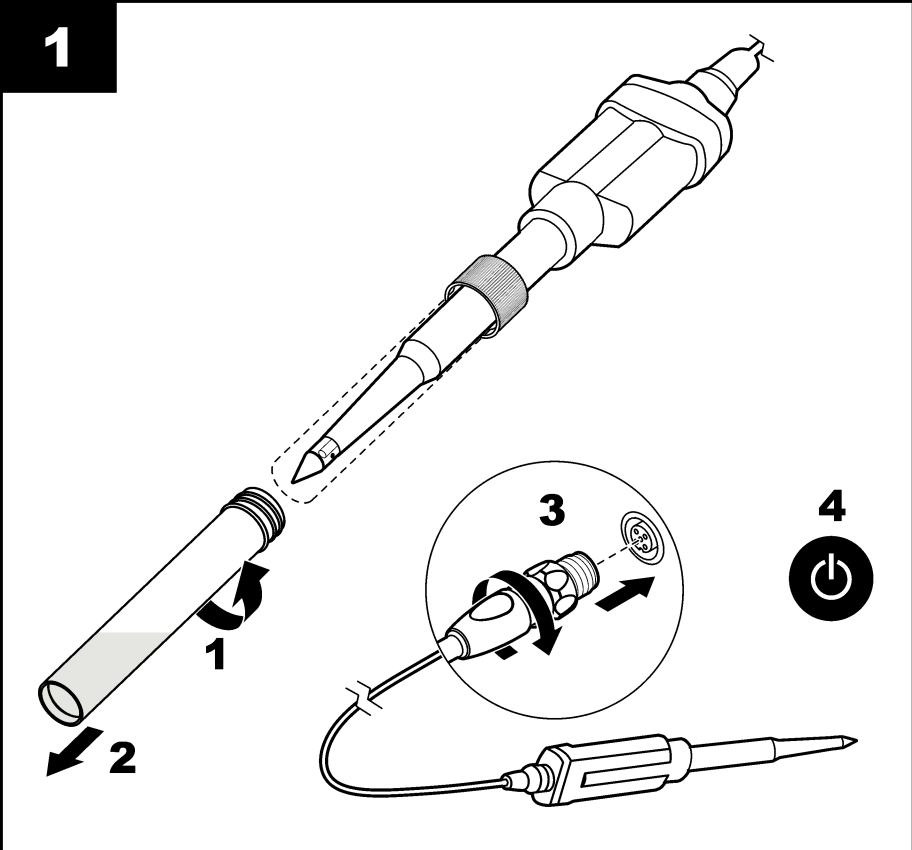
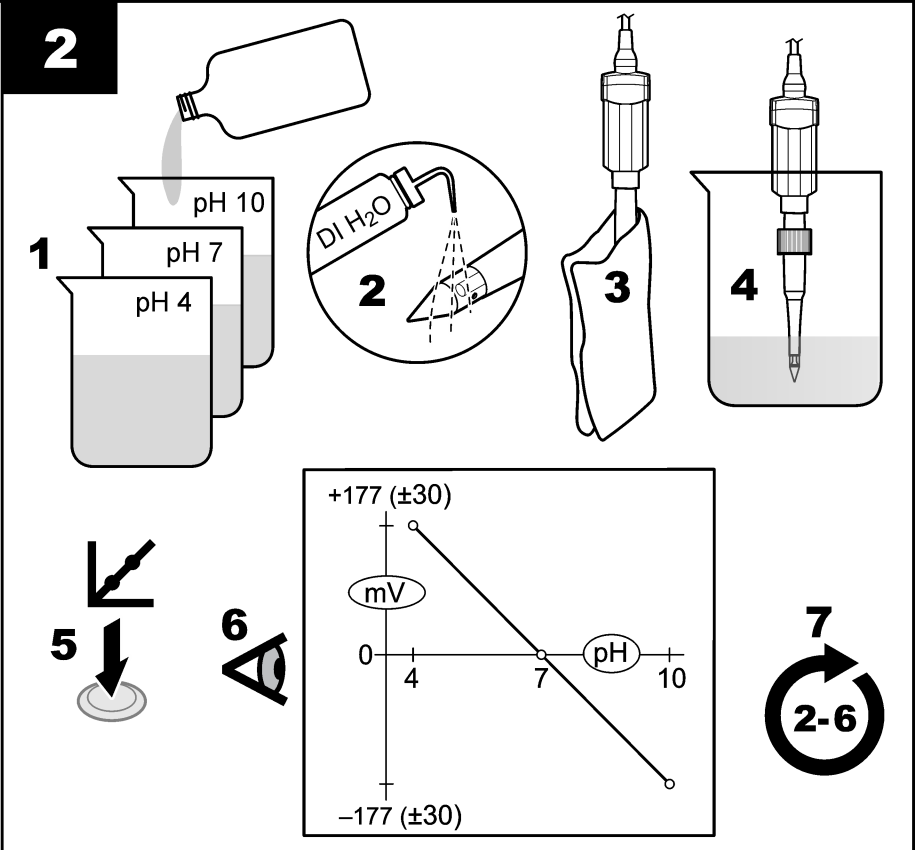
3

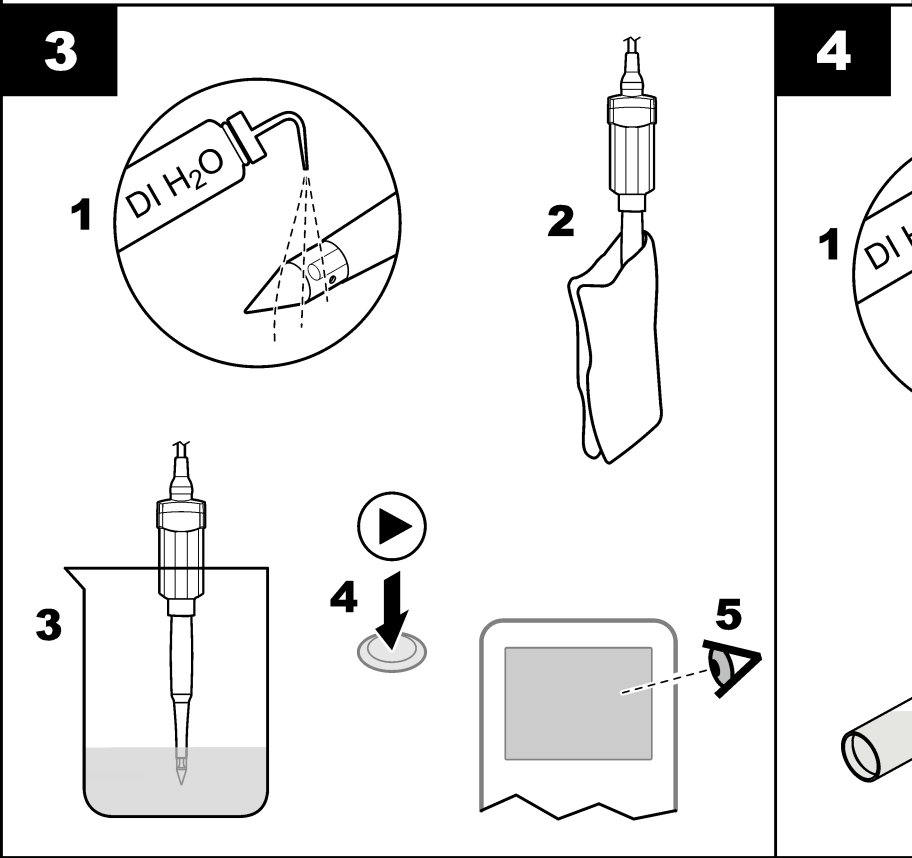4

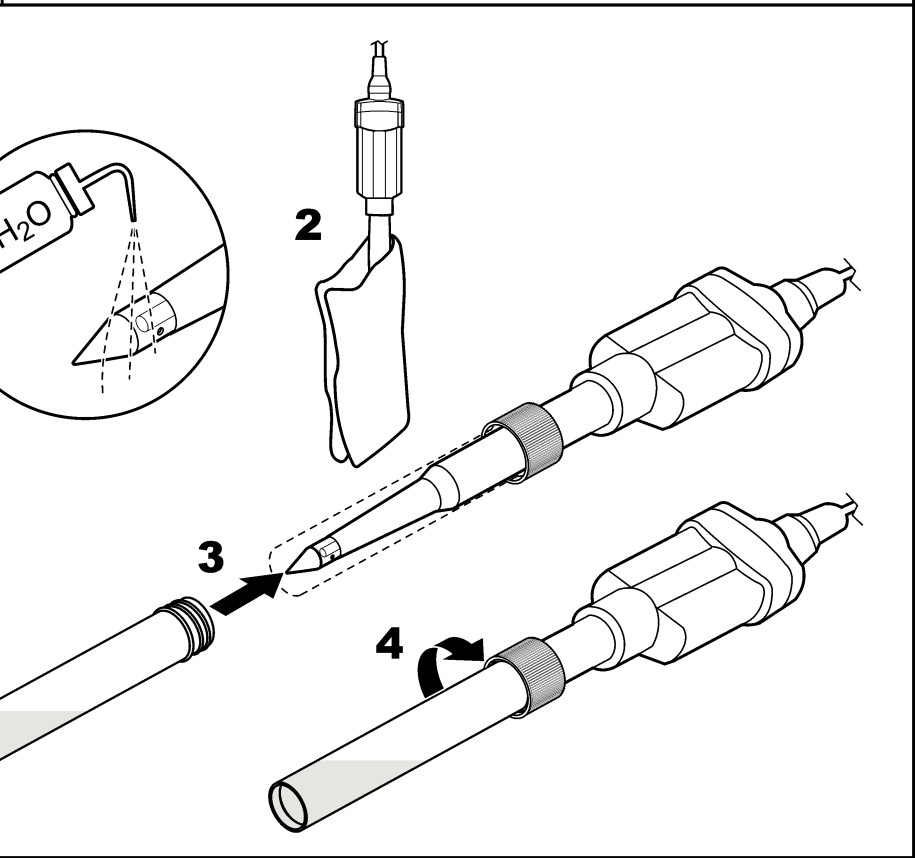PHC108
Introduction
The probe is a pH electrode with non-refillable electrolyte and an integrated temperature sensor. The probe has a spear-shaped glass sensor that measures pH in semi-solid media applications.
Warning
 **Chemical exposure hazard.** Obey laboratory safety procedures and wear all of the personal protective equipment appropriate to the chemicals that are handled. Refer to the current safety data sheets (MSDS/SDS) for safety protocols.
User manual
A user manual with all instructions is on the manufacturer's website.
Items to collect:
Compatible meter; pH standard solutions; beakers; DI water; lint-free cloth.

1 Connect the probe
2 Calibration
3 Measurement
Do not use the probe in applications where must apply pressure to introduce the probe into the sample (e.g., solids). For these applications, it is necessary to first open a path with a stainless steel perforator.
4 Storage

Technical information	
pH range	pH 2 to 12
Slope	-59 mV/pH (85 to 115% at 25 °C)
Temperature	0 to 60 °C (32 to 140 °F)
Minimum sample depth	15 mm (0.59 in.)
Storage solution	Hach pH electrode storage solution or 3 M KCl

Maintenance and storage
For maintenance and storage details, refer to the expanded version of this manual on the manufacturer's website.
Warranty
1 year for manufacturing faults. This warranty does not cover improper use or wear.

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