
DOC022.97.80339

QuikChem[®] 8500 Series 2

04/2015, Edition 2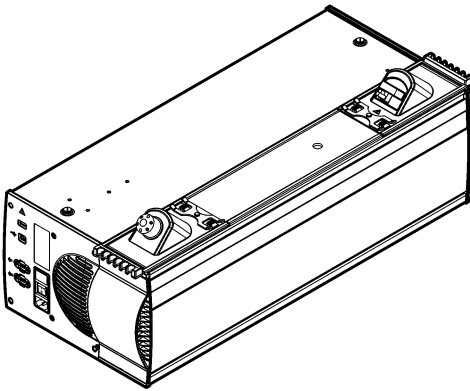

Installation
Installation
Instalación
安装

English 3

Français 21

Español 40

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Table of contents

[Specifications](#) on page 3

[General information](#) on page 4

[Installation](#) on page 7

Additional information

Additional information is available on the manufacturer's website.

Specifications

Specifications are subject to change without notice.

Specification	Details
Dimensions bench space (W x D x H)	2.72 x 0.83 x 0.58 m (9 x 2.75 x 2 ft) (recommended)
Location	Laboratory (indoor use only)
Pollution degree/installation category	II
Protection class	I
Operating temperature	5 to 40 °C (41 to 104 °F), maximum 80% relative humidity non-condensing for temperatures up to 31 °C and decreases linearly to 50% relative humidity at 40 °C
Storage temperature	-20 to 70 °C (-4 to 158 °F)
Altitude limit	Up to 3000 m (10,000 feet) above sea level
Certifications	CE: IEC/EN 61010-1 per LV Directive (2006/95/EC), IEC/EN 61326 per EMC: Directive (2004/108/EC)
	cETL: CSA C22.2 No 61010-1
	ETLus: UL 61010-1
Warranty	1 year
System unit (5 channels)	
Dimensions (W x D x H)	70.5 x 71.2 x 26.7 cm (27.7 x 28.0 x 10.5 in.)
Weight	46.5 kg (102.51 lb)
Power requirements	100–115 V ±10%, 230 V ±10%; 50–60 Hz, 800 W
Fuse rating	T, 8 A, 100–115 V, T, 4 A, 230 V
Lamp	Tungsten/Halogen, 12 V, 50 W, MR16
Sample processing module (SPM)	
Dimensions (W x D x H)	69 x 9 x 26 cm (27.2 x 3.5 x 10.2 in.)
Weight	4 kg (8.8 lb)
Power requirements	115 VAC, 1.2 A, 60 Hz or 230 VAC, 0.8 A, 50 Hz
Fuse rating	T, 1.6 A, 250 V, 5 mm x 20 mm
Pump (12 position)	
Dimensions (W x D x H)	15 x 26 x 26.7 cm (5.9 x 10.2 x 10.5 in.)
Weight	5.6 kg (12.3 lb)

Specification	Details
Power requirements	100 or 120 VAC, 220 or 240 VAC; 50–60 Hz, 15 VA
Fuse rating	T, ½ A, 250 V
ASX-520 sampler	
Dimensions (W x D x H)	52.5 x 46 x 25 cm (20.6 x 18.1 x 9.8 in.)
Weight	10 kg (22 lb)
Power requirements	100–240 VAC; 50–60 Hz, 1.5 A
Fuse rating	Not user-replaceable
ASX-260 sampler	
Dimensions (W x D x H)	30 x 50.8 x 61 cm (13 x 20 x 24 in.)
Weight	8.4 kg (18.5 lb)
Power requirements	100–240 VAC; 50–60 Hz, 1.5 A
Fuse rating	Not user-replaceable
PDS200 dilutor (optional)	
Dimensions (W x D x H)	12.7 x 22.5 x 23.4 cm (5 x 8.9 x 9.2 in.)
Weight	4 kg (8.8 lb)
Power requirements	110–240 VAC; 50–60 Hz, 0.5 A
Fuse rating	T, ½ A, 250 V

General information

In no event will the manufacturer be liable for direct, indirect, special, incidental or consequential damages resulting from any defect or omission in this manual. The manufacturer reserves the right to make changes in this manual and the products it describes at any time, without notice or obligation. Revised editions are found on the manufacturer's website.

Safety information

NOTICE

The manufacturer is not responsible for any damages due to misapplication or misuse of this product including, without limitation, direct, incidental and consequential damages, and disclaims such damages to the full extent permitted under applicable law. The user is solely responsible to identify critical application risks and install appropriate mechanisms to protect processes during a possible equipment malfunction.

Please read this entire manual before unpacking, setting up or operating this equipment. Pay attention to all danger and caution statements. Failure to do so could result in serious injury to the operator or damage to the equipment.

Make sure that the protection provided by this equipment is not impaired. Do not use or install this equipment in any manner other than that specified in this manual.

Use of hazard information

▲ DANGER

Indicates a potentially or imminently hazardous situation which, if not avoided, will result in death or serious injury.

▲ WARNING

Indicates a potentially or imminently hazardous situation which, if not avoided, could result in death or serious injury.

⚠ CAUTION

Indicates a potentially hazardous situation that may result in minor or moderate injury.

NOTICE

Indicates a situation which, if not avoided, may cause damage to the instrument. Information that requires special emphasis.

Precautionary labels

Read all labels and tags attached to the instrument. Personal injury or damage to the instrument could occur if not observed. A symbol on the instrument is referenced in the manual with a precautionary statement.

	This symbol, if noted on the instrument, references the instruction manual for operation and/or safety information.
	Electrical equipment marked with this symbol may not be disposed of in European public disposal systems after 12 August of 2005. In conformity with European local and national regulations (EU Directive 2002/96/EC), European electrical equipment users must now return old or end-of-life equipment to the Producer for disposal at no charge to the user. <i>Note: For return for recycling, please contact the equipment producer or supplier for instructions on how to return end-of-life equipment, producer-supplied electrical accessories, and all auxiliary items for proper disposal.</i>
	This symbol indicates that a risk of electrical shock and/or electrocution exists.
	This symbol indicates the need for protective eye wear.
	This symbol indicates that the marked item can be hot and should not be touched without care.
	This symbol, when noted on the product, identifies the location of a fuse or current limiting device.
	This symbol indicates that the object is heavy.

Certification

Canadian Radio Interference-Causing Equipment Regulation, IECS-003, Class A:

Supporting test records reside with the manufacturer.

This Class A digital apparatus meets all requirements of the Canadian Interference-Causing Equipment Regulations.

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FCC Part 15, Class "A" Limits

Supporting test records reside with the manufacturer. The device complies with Part 15 of the FCC Rules. Operation is subject to the following conditions:

1. The equipment may not cause harmful interference.
2. The equipment must accept any interference received, including interference that may cause undesired operation.

Changes or modifications to this equipment not expressly approved by the party responsible for compliance could void the user's authority to operate the equipment. This equipment has been tested and found to comply with the limits for a Class A digital device, pursuant to Part 15 of the FCC rules. These limits are designed to provide reasonable protection against harmful interference when the equipment is operated in a commercial environment. This equipment generates, uses and can radiate radio frequency energy and, if not installed and used in accordance with the instruction manual, may cause harmful interference to radio communications. Operation of this equipment in a residential area is likely to cause harmful interference, in which case the user will be required to correct the interference at their expense. The following techniques can be used to reduce interference problems:

1. Disconnect the equipment from its power source to verify that it is or is not the source of the interference.
2. If the equipment is connected to the same outlet as the device experiencing interference, connect the equipment to a different outlet.
3. Move the equipment away from the device receiving the interference.
4. Reposition the receiving antenna for the device receiving the interference.
5. Try combinations of the above.

System overview

▲ DANGER	
	<p>Chemical or biological hazards. If this instrument is used to monitor a treatment process and/or chemical feed system for which there are regulatory limits and monitoring requirements related to public health, public safety, food or beverage manufacture or processing, it is the responsibility of the user of this instrument to know and abide by any applicable regulation and to have sufficient and appropriate mechanisms in place for compliance with applicable regulations in the event of malfunction of the instrument.</p>
▲ WARNING	
	<p>Fire hazard. The user is responsible to make sure that sufficient precautions are taken when the equipment is used with methods that use flammable liquids. Make sure to obey correct user precautions and safety protocols. This includes, but is not limited to, spill and leak controls, proper ventilation, no unattended use, and that the instrument is never left unattended while power is applied.</p>

The QuikChem[®] 8500 Series 2 FIA (Flow Injection Analysis) System Unit is configured to complete a flow injection analysis with a photometric detection. Alternate configurations, such as potentiometric and amperometric detection methods are described in the analytical method where they are necessary. For complete information on the software, refer to the online help system in the software.


The QuikChem[®] 8500 Series 2 analyzer is modular in design and includes the modules that follow. Refer to [Figure 1](#).

- **QuikChem[®] FIA System Unit**—The System Unit includes the core unit and the sample processing modules (SPM). The core unit supplies power, light and electronic signals to and from each SPM. An SPM or a channel has an injection valve, a manifold, a photometric detector and a heating module (optional). A System Unit can have up to five channels that operate simultaneously. A single computer controls two System Units in order to operate up to eight channels at once. A chemical reaction module or a manifold is installed on a channel for the quantitative analysis of an analyte. There are 450 QuikChem methods available to find over 40 analytes. There are many methods in several matrices.
- **Sampler**—The QuikChem[®] 8500 Series 2 analyzer operates with either the ASX-260 or the ASX-520 series autosamplers.
- **Pump (12 or 16 cartridges)**—The multichannel pump is used to pump reagents to the manifold. It also aspirates sample aliquots and moves them from the sampler to the valve. These portions are then loaded into the sample loop and moved to the manifold by the carrier stream to mix with the reagents.

- Computer with Omnion software loaded
- Printer (optional)
- Dilutor (optional)—The PDS200 uses positive displacement to pull diluent water to mix with non-flammable samples for analysis.

Make sure that all components have been received. If any items are missing or damaged, contact the manufacturer or a sales representative immediately.

Figure 1 System overview

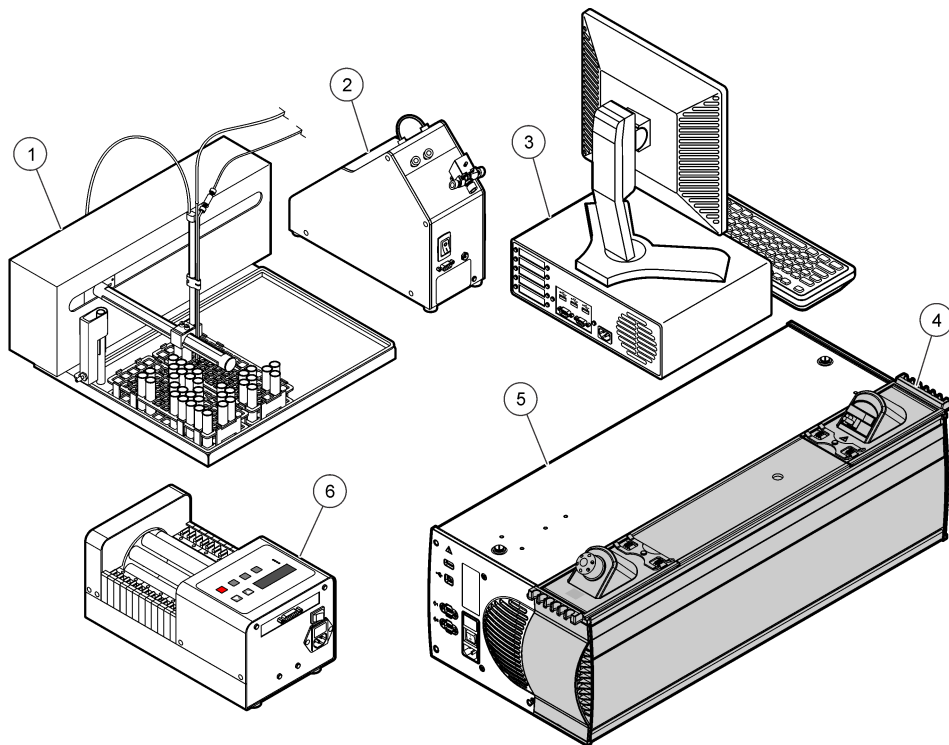1 Sampler	4 Channel (1–5 on 1 system, up to 8 on 2 systems)
2 Dilutor (optional)	5 QuikChem® FIA System Unit
3 Computer with Omnion software loaded	6 Pump

Installation

▲ DANGER

Multiple hazards. Only qualified personnel must conduct the tasks described in this section of the document.

▲ DANGER

Personal injury hazard. Make sure that the equipment is stable and use assistance to install and move.

Attach the Z-drive to the sampler


Refer to the instructions that come with the sampler.

Attach the probe to the sampler

When the probe is installed, make sure that the probe is flush with the hole in the guide plate. Refer to [Figure 2](#).

Note: Always move the sliders with the wheel on the sampler to prevent damage to the tubing.

Figure 2 Probe installation

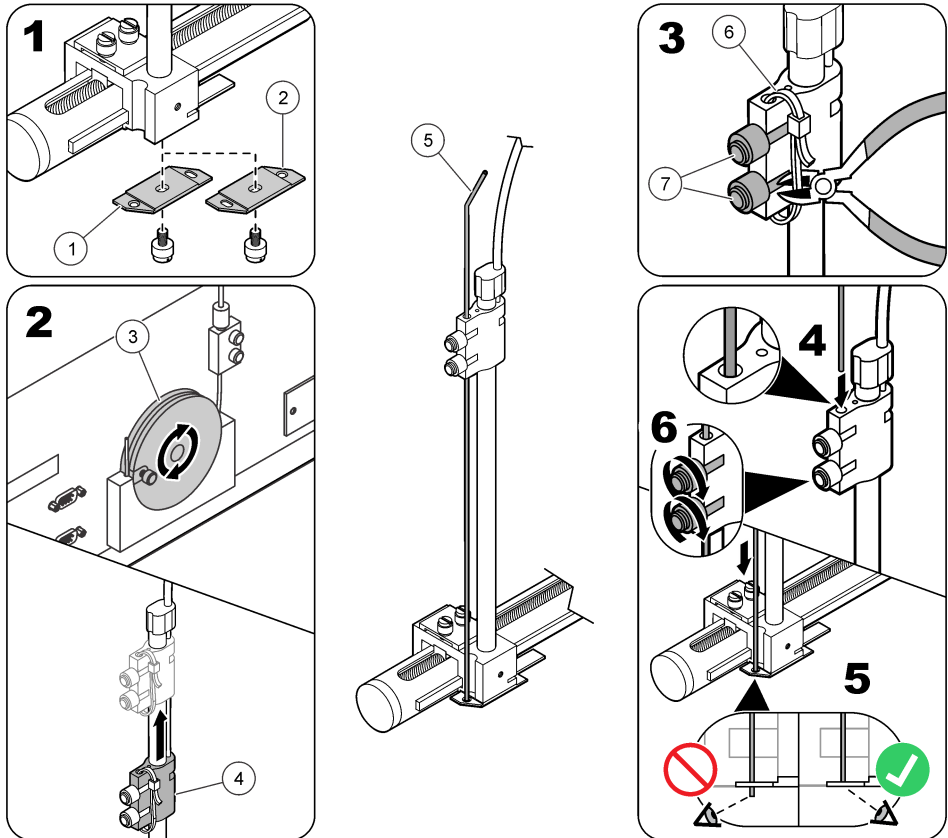1 Probe guide plate (single probe installation)	4 Probe slider	7 Slider knobs
2 Probe guide plate (dual probe installation)	5 Probe	
3 Slider wheel	6 Tie wrap	

Install additional channels

If additional channels were ordered, attach the channels that are shipped separately. Use the documentation that is shipped with the channels.

Electrical installation

⚠ DANGER

Electrocution hazard. Always remove power to the instrument before making electrical connections.

⚠ DANGER

Electrocution hazard. Do not connect AC power directly to a DC powered instrument.

⚠ DANGER

Electrocution hazard. If this equipment is used outdoors or in potentially wet locations, a Ground Fault Circuit Interrupt (GFCI/GFI) device must be used for connecting the equipment to its main power source.

⚠ DANGER

Electrocution hazard. Protective Earth Ground (PE) connection is required.

⚠ DANGER

Electrocution and fire hazards. For instruments with a power cord, make sure that there is easy access to the local power disconnect.

⚠ WARNING

Electrical shock hazard. Externally connected equipment must have an applicable country safety standard assessment.


NOTICE

Make sure that the equipment is connected to the instrument in accordance with local, regional and national requirements.

Connect the interfaces

Refer to [Figure 3](#).


Figure 3 System interface connections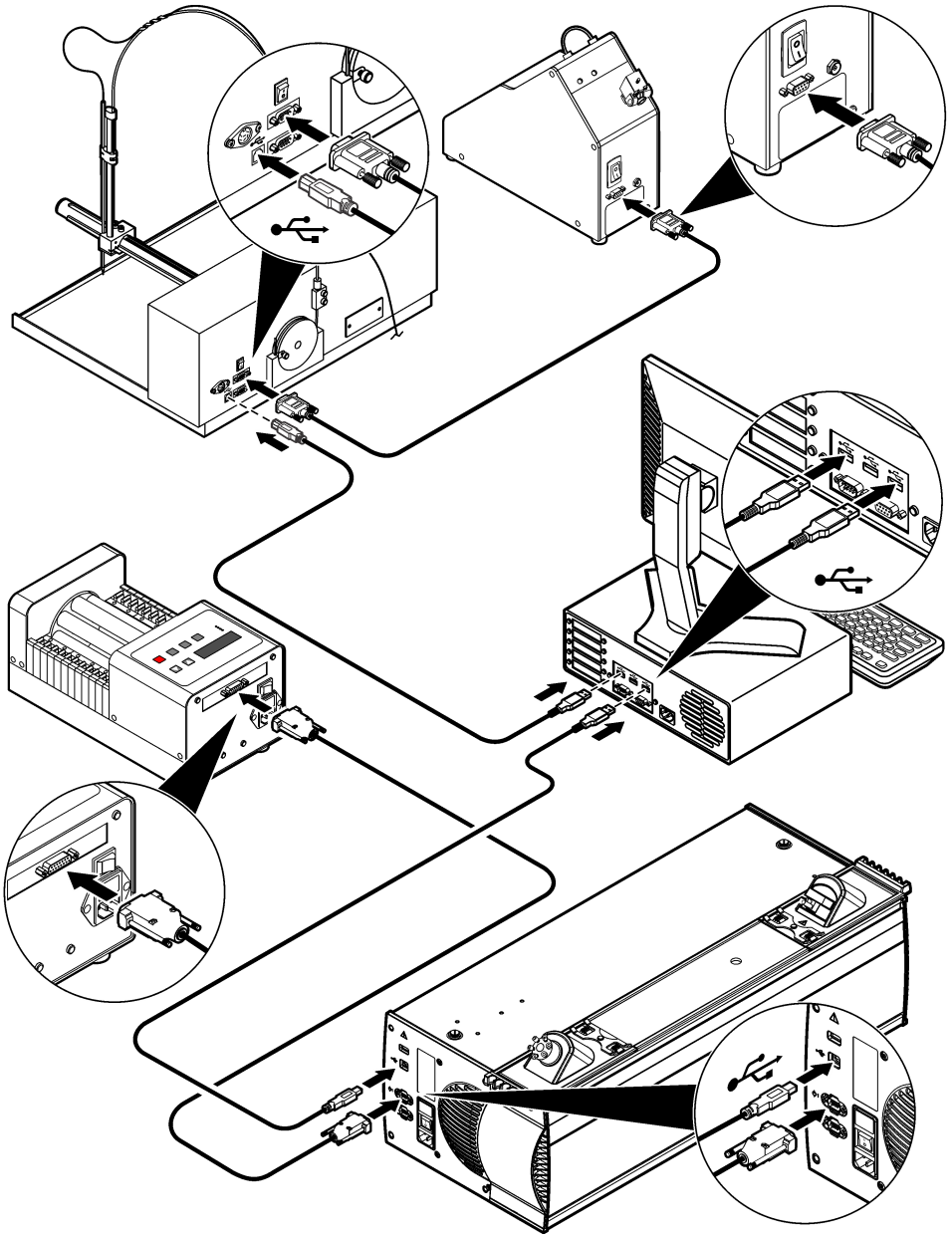

Connect to power

Do not set the power to on for these instruments or the power strip. Refer to [Figure 4](#).

Figure 4 System power connections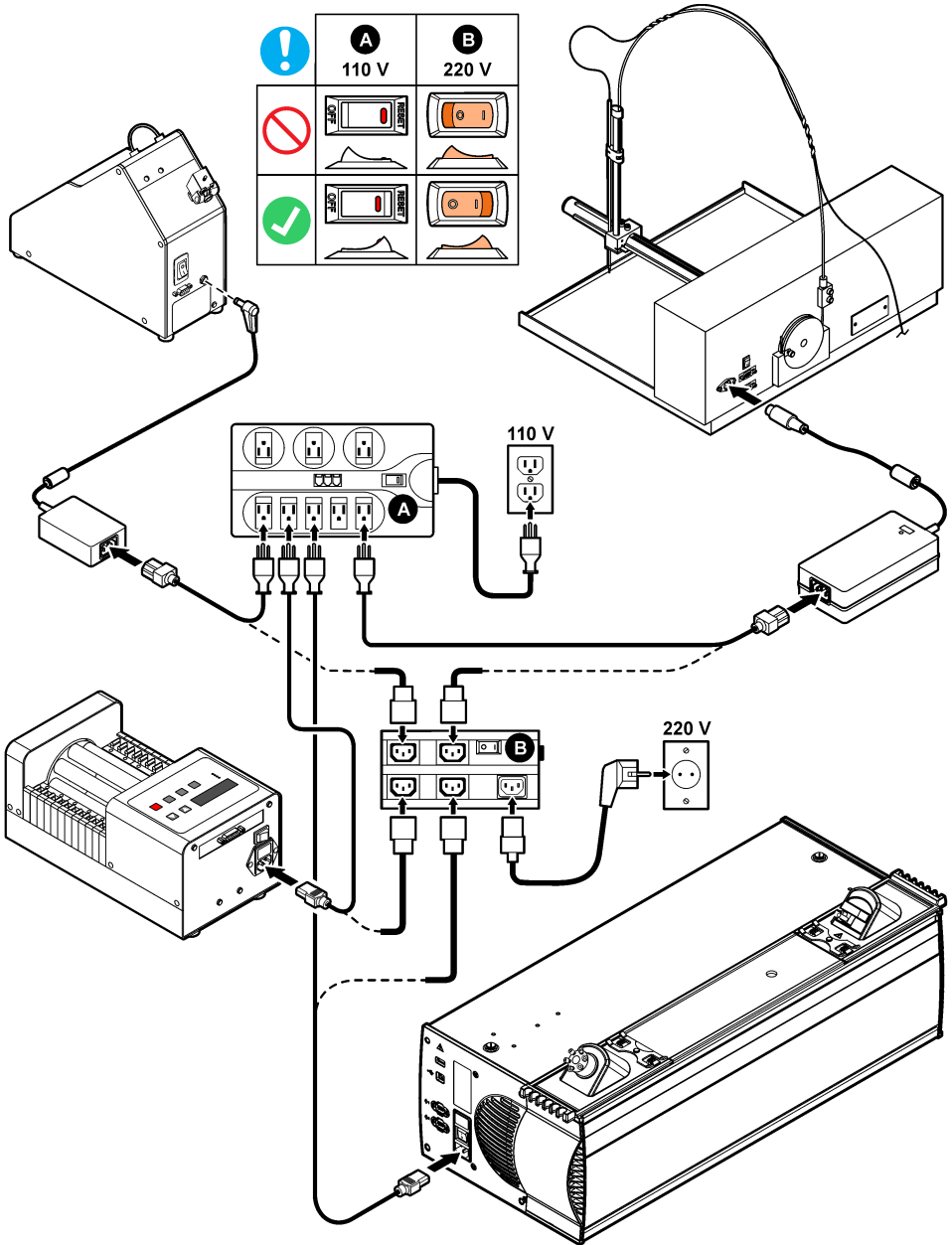


Plumbing

The plumbing includes tubing attachment to the pump, sampler and dilutor. Make sure to use the specified tubing size. Refer to the QuikChem method for details.

Install the pump tubing


1. Remove all of the cartridges from the pump. Refer to [Figure 5](#).
2. Attach the tubing to the cartridge. Refer to [Figure 6](#).
3. Attach the cartridge to the pump. Set the tension lever one click to the left. Refer to [Figure 7](#).

Figure 5 Cartridge removal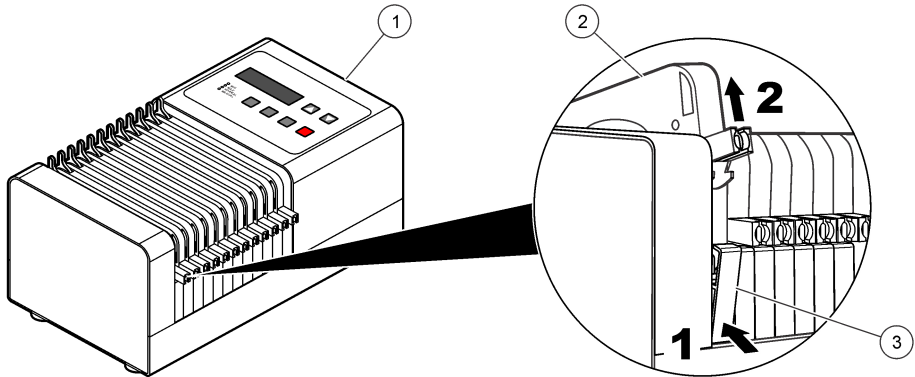


1 Pump	2 Pump cartridge	3 Cartridge holder
--------	------------------	--------------------

Figure 6 Attach cartridge and tabs to the tubing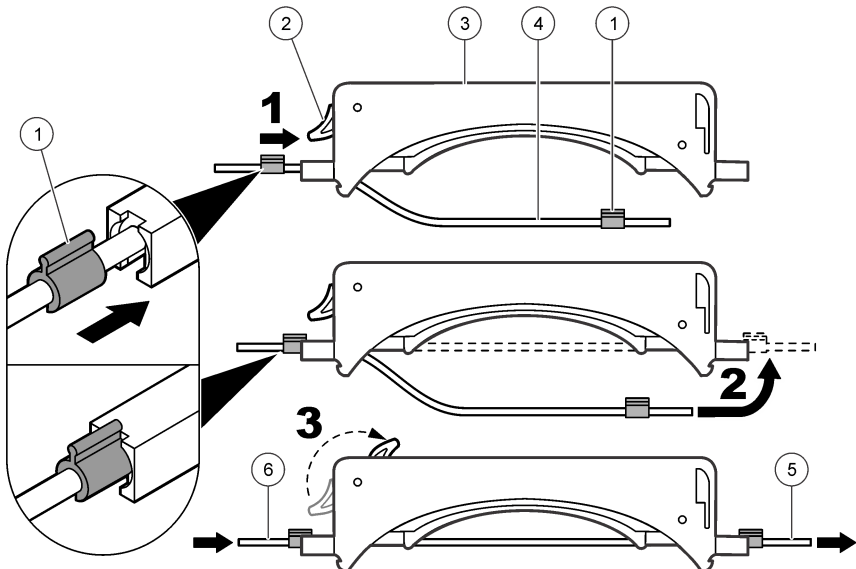

1 Color-coded tab	3 Pump cartridge	5 To injection valve, wash station or manifold
2 Tension lever	4 Pump tubing	6 From reagent, sample, wash solution, buffer or other chemicals (varies with method)

Figure 7 Cartridge installation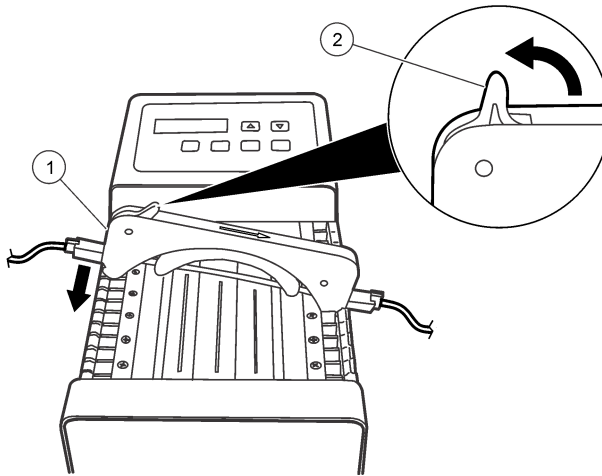


1 Pump cartridge

2 Tension lever at 12 o'clock position

Install the sampler tubing

To install the sampler tubing, refer to [Figure 8](#).

Figure 8 Sampler tubing installation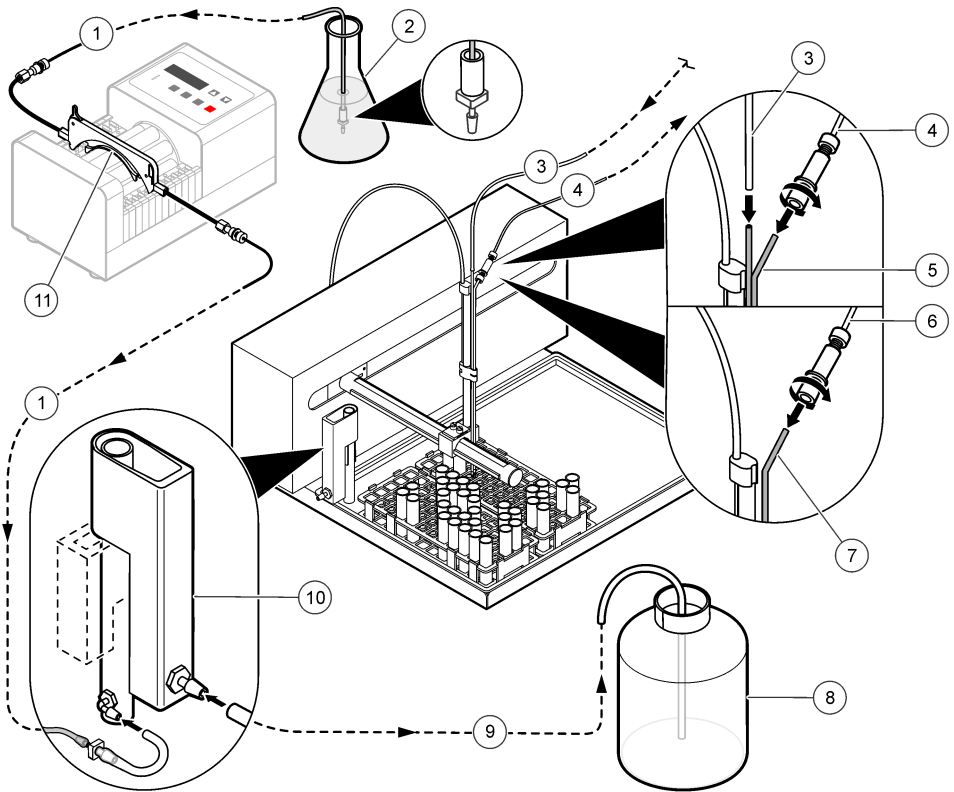


1 Rinse solution tube assembly	5 Dual probe (for dilutor option)	9 Rinse station waste tube
2 Rinse solution	6 Sample tubing to pump and injection valve	10 Sampler rinse station
3 Dilution tube from dilutor	7 Sample probe	11 Pump cartridge
4 Sample tube to dilutor	8 Waste container	

Install the dilutor tubing

To install the dilutor tubing, refer to [Figure 9](#).

Figure 9 Dilutor tubing connections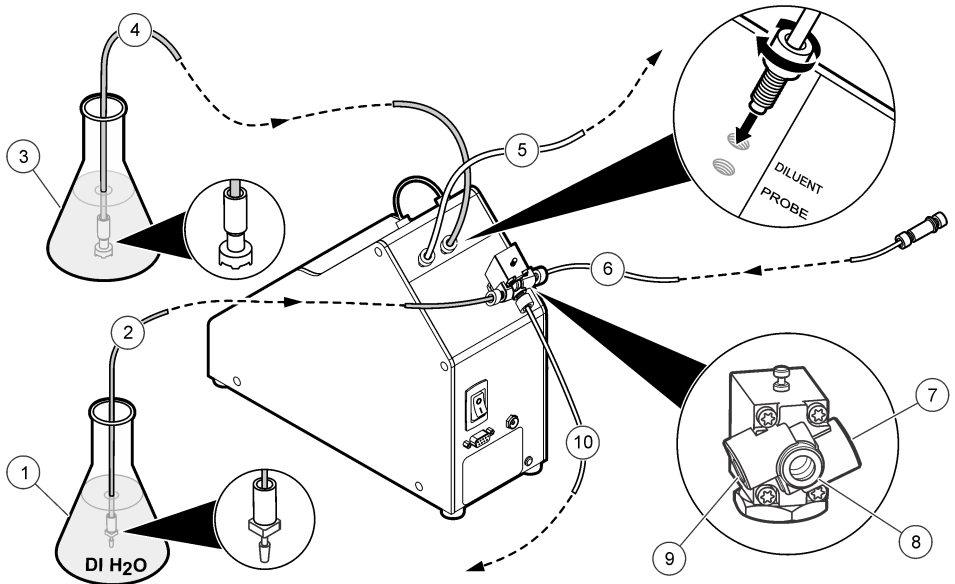


1 DI water	5 Tube to sampler probe	9 Solenoid valve port 1
2 Tube to solenoid valve	6 Sample tube from sample probe	10 Sample tube to pump and injection valve
3 Diluent solution	7 Solenoid valve port 2	
4 Tube to dilutor	8 Solenoid valve port 3	

Install the heater mandrel

If the instrument has a heater module, a heater mandrel must be installed. To install a heater mandrel:

1. Make sure that the heater mandrel has the applicable wrapping.
2. Remove the channel cover. Refer to [Figure 10](#).
3. Install the mandrel, the tubing and the thermocouple jack.
4. Install the manifold board.

Figure 10 Heater mandrel installation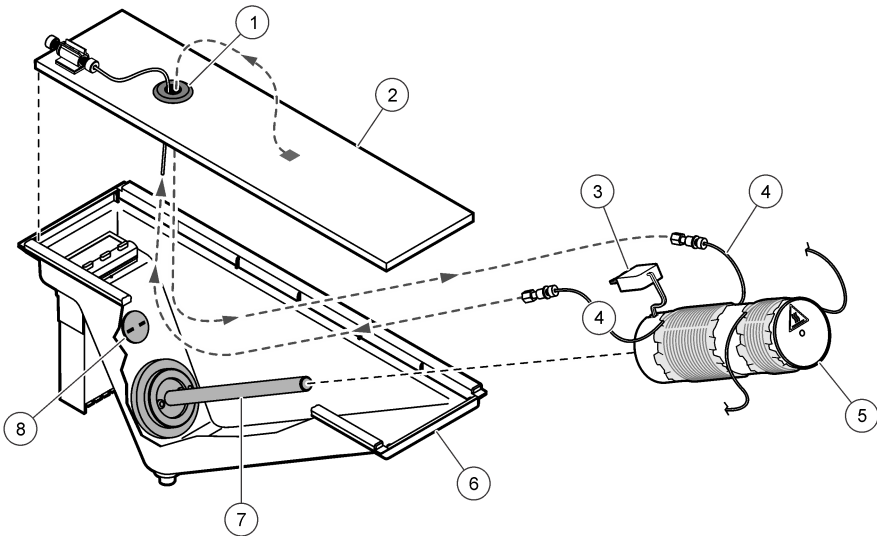


1 Access hole for heater tubing	4 Tubing to manifold	7 Heater element
2 Manifold board	5 Heater mandrel	8 Thermocouple jack
3 Thermocouple plug	6 Drip tub	

Install the external manifold plumbing

Connect the tubing to the manifold. Refer to [Figure 11](#).

Figure 11 External manifold plumbing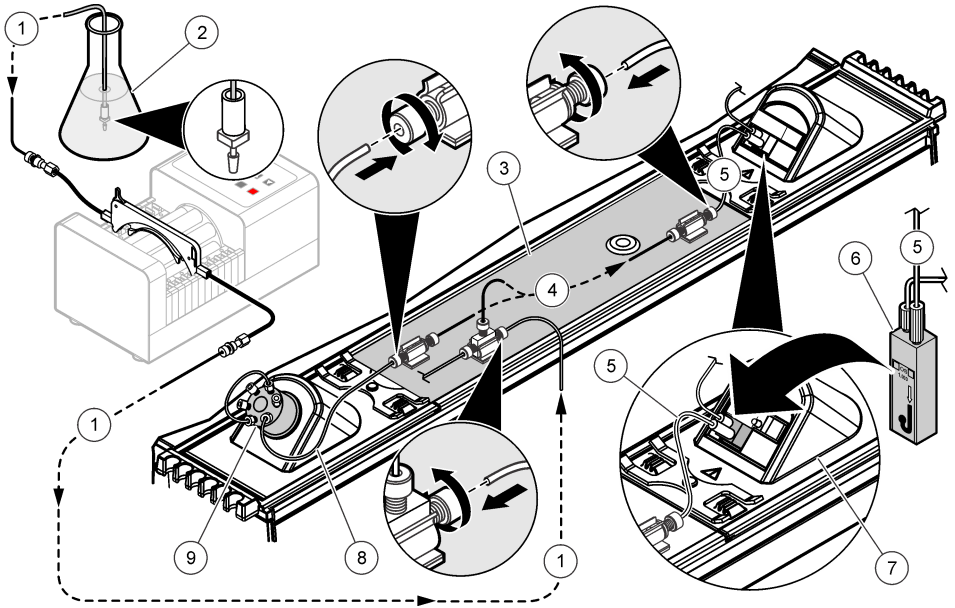


1 Transmission solution tubing to manifold	4 Manifold plumbing (varies with method)	7 Detector
2 Reagent solution (varies with method)	5 Tubing from manifold to flow cell	8 Tubing from valve port 3 to manifold
3 Manifold board	6 Flow cell	9 Injection valve

Install the channel plumbing

Connect the tubing to the channel ([Figure 12](#)).

Figure 12 Channel plumbing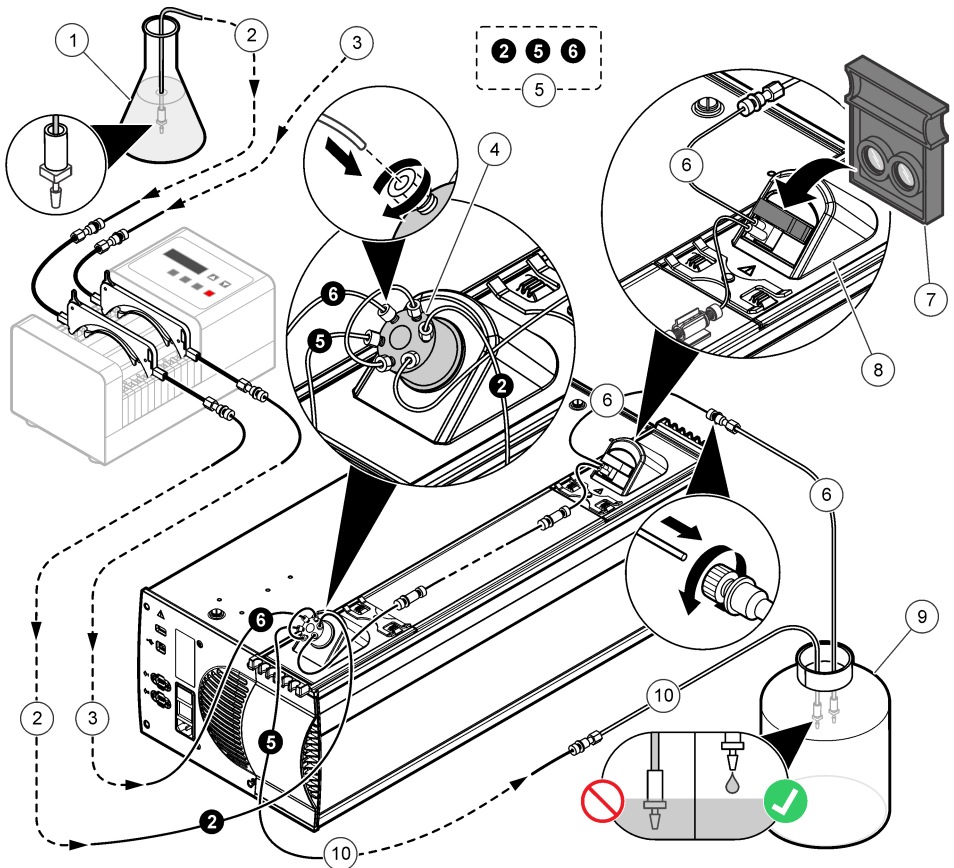


1 Carrier solution (varies with QuikChem method)	5 Injection valve port numbers	9 Waste container
2 Carrier tubing to injection valve	6 Waste tubing from flow cell	10 Waste tubing from injection valve
3 Sample tubing to injection valve	7 Interference filter	
4 Injection valve	8 Detector	

Install the plumbing to multiple channels

When multiple channels are used, use 15 cm of 0.8 mm Teflon® tubing to connect port 5 on the first valve to port 6 on the next valve (Figure 13).

Note: Do not put the tubing in the waste solution, or back pressure and flow restrictions can occur. Make sure that the waste solution always drips in the container.

Figure 13 Multi-channel plumbing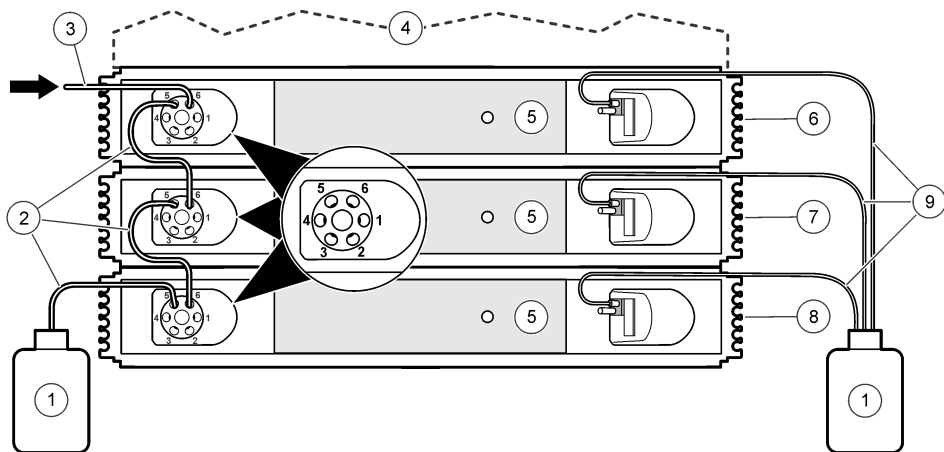

1 Waste container	4 System unit	7 Channel 2
2 Injection valve waste tubing	5 Manifolds	8 Channel 3
3 Sample tubing in	6 Channel 1	9 Flow cell waste tubing


Non-photometric detectors

Other types of detectors are necessary for some methods, such as fluoride, flame photometers, conductivity or pH. The regular flow cell or interference filters are not necessary for non-photometric detectors, such as an ISE detector for fluoride determination. Refer to the instructions in the QuikChem method for fluidic connections.

Set the power to on

1. Set the power to on for each instrument. Refer to [Figure 14](#).
2. Set the power to on for the power strip.

Figure 14 Power switches and buttons

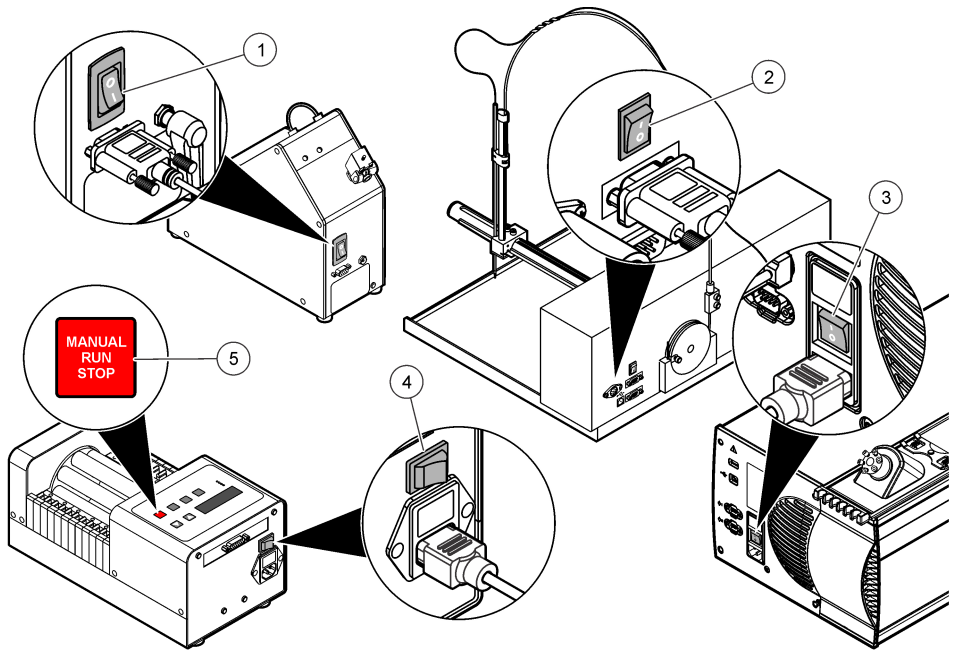1 Dilutor power switch	4 Pump power switch
2 Sampler power button	5 Pump start/stop button
3 System unit power switch	

Do the initial pump test

1. Make sure that the Omnion software is not on at this time. Push the **NORMAL RUN** button. Make sure that the pump operates at a normal speed of 35. The NORMAL LED illuminates. If the display shows a different speed, push the **ARROW** button to change the setting to 35.
2. Push the **MIN** button. The MIN LED illuminates. The display shows a speed of 4.
3. Hold down the **MAX** button until the MAX LED illuminates. The display shows a speed of 999. The pump goes back to the original speed when the **MAX** button is released.
4. To stop the pump, push the **MANUAL RUN/STOP** button. No LEDs illuminate, but the display shows "35" (the pump speed).
5. Push the **MANUAL RUN/STOP** button again. The MANUAL LED illuminates. The pump operates at a normal speed of 35.
6. Push the **MANUAL RUN/STOP** button to stop the pump. To do a step from above again, set the pump power to off and then to on. Start with step 1.
7. If there is a failure on a step from above, contact technical support.

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P.O. Box 389, Loveland, CO 80539-0389 U.S.A.
(800) 227-4224 (U.S.A. only)