TitraLab[®]
SAC850 & SAC950
Sample Changers

User's Guide

RADIOMETER ANALYTICAL SAS
72 rue d'Alsace, 69627 Villeurbanne Cedex, France
E-mail: radiometer@analytical.com Web: www.radiometer-analytical.com
Tel. : +33 (0)4 78 03 38 38 - Fax : +33 (0)4 78 68 88 12

Safety information

Please read this entire manual before unpacking, setting up, or operating this equipment. Pay attention to all danger and caution statements. Failure to do so could result in serious injury to the operator or damage to the equipment. To ensure that the protection provided by this equipment is not impaired, do not use and do not install this equipment in any manner other than that specified in this manual.

Precautionary labels

Read all labels and tags attached to the instrument. Personal injury or damage to the instrument could occur if not observed. A symbol, if noted on the instrument, will be included with a danger or caution statement in the manual.

	<p>This symbol, if noted on the instrument, references the instruction manual for operation and/or safety information.</p>
	<p>Electrical equipment marked with this symbol may not be disposed of in European public disposal systems after 12 August of 2005. In conformity with European local and national regulations (EU Directive 2002/96/EC), European electrical equipment users must now return old or end-of life equipment to the Producer for disposal at no charge to the user.</p> <p>Note: For return for recycling, please contact the equipment producer or supplier for instructions on how to return end-of-life equipment, producer-supplied electrical accessories and all auxiliary items for proper disposal.</p>

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Contents

1. Introduction.....	25
2. Presentation.....	29
SAC850	29
SAC950	31
RFID technology on turntables	34
3. Setting up.....	37
Mounting the turntable.....	38
Placing beakers	39
Mounting the propeller stirrer.....	41
Mounting the electrodes and addition tips	42
Connections.....	43
Assembling option accessories	47
Dynamic rinse module.....	47
Reagent addition module	56
Beaker cover module	62
4. Use.....	65
Switching on	65
Running sequence with a workstation	66
5. Troubleshooting	73
Autotest and manual mode errors	73
Errors displayed by the workstation.....	74

6. Maintenance	79
Cleaning	79
Transporting the instrument	79
Servicing	79
Replacing the fuses	80
Adjusting the beaker detector position	81
Adjusting the mechanics reference positions	83
7. Accessories	85
8. Technical specifications	93

1. Introduction

The SAC850 and SAC950 Sample Changers are designed for automatic analysis in various types of trays.

Each tray can hold up to 70 samples depending on the sample volume and the application to be performed.

The SAC850 Basic model is dedicated for small sample series automation. The SAC850 is fitted with a single radius tray and can hold 14 to 25 samples.

The SAC950 Extended model is dedicated for large sample series automation. The SAC950 is fitted with a 1, 2 or 3-radii tray and can hold 14 to 70 samples.

The SAC850 and SAC950 Sample Changers are ideal for handling samples in TitraLab® or MeterLab® compatible systems.

The aim of this User's Guide is to provide you with detailed information about setting up and operating your sample changer.

For information on how to enter data into your workstation, please refer to the instructions delivered with the workstation.

Warning !

The sample changer has been developed to meet the requirements of volumetric titration applications. It is therefore aimed at experienced users who have the knowledge required to operate the instrument and implement the security instructions enclosed. Please remember that the sample changer must not, under any circumstances, be used to perform tests on living beings.

We accept no responsibility for using the sample changer and its peripheral devices under conditions that are not specified in this User's Guide.

Compliance with FCC rules, part 15

Information to the user



NOTE: This equipment has been tested and found to comply with the limits for a Class B digital device, pursuant to part 15 of the FCC Rules. These limits are designed to provide reasonable protection against harmful interference in a residential installation. This equipment generates, uses and can radiate radio frequency energy and, if not installed and used in accordance with the instruction, may cause harmful interference to radio communications. However, there is no guarantee that interference will not occur in a particular installation. If this equipment does cause harmful interference to radio or television reception which can be determined by turning the equipment off and on, the user is encouraged to try to correct interference by one or more of the following measures:

- Reorient or relocate the receiving antenna.
- Increase the separation between the equipment and receiver.
- Connect the equipment into an outlet on circuit different from that to which the receiver is connected.
- Consult the dealer or an experienced radio/TV technician for help.
- Use a shielded RS232 serial cord and a shielded USB cord fitted with a ferrite.

Warning!!

Changes and modifications that are not expressly approved by the party responsible for compliance could void the user's authority to operate the equipment.

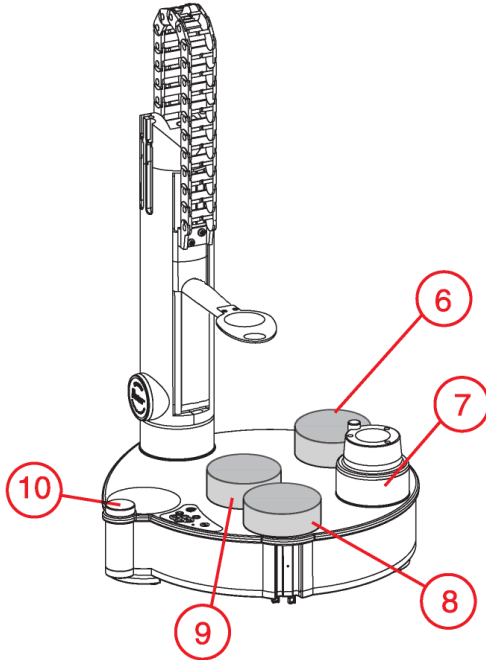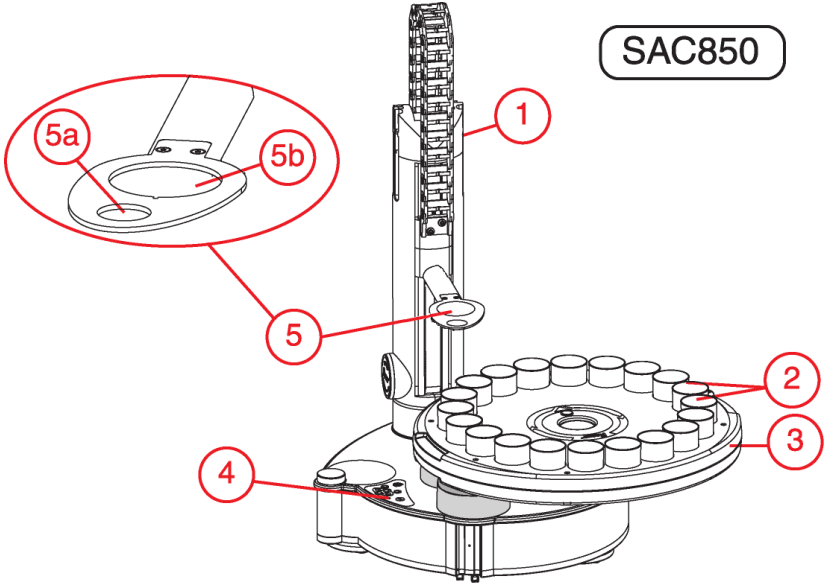SAC850


2. Presentation

SAC850

- ① 4 locations for spare electrodes and antidiffusion delivery tips (2 locations on each side).
- ② Sample/Rinse/Park beakers (1 Park, 0 to 2 Rinses, remaining positions dedicated for samples).
- ③ Mono radius turntable. Capacities depend on the beaker size and diameter : see chapter "Accessories" - Table "Turntables and beakers (SAC850 and SAC950)".
- ④ Keypad.
- ⑤ Electrodes/Stirrer/Beaker presence detector holder.
Beaker presence detector location **(5a)**.
Electrodes/stirrer head holder location **(5b)**.
- ⑥ Optional peristaltic pump (optional).
- ⑦ Turntable holder.
- ⑧ Optional shower rinse pump/emptying pump module. Comprises 2 diaphragm pumps (rinsing and emptying).
- ⑨ Magnetic stirrer assembly comprising also a RFID reader device for turntable automatic recognition.
- ⑩ Status lamp:
Blue, blinking : Free for operation. Manual operations possible.
Purple: Sample changer controlled by the workstation (analysis in progress). Only the keyboard **Stop** key is available.
Red: An error has been detected.

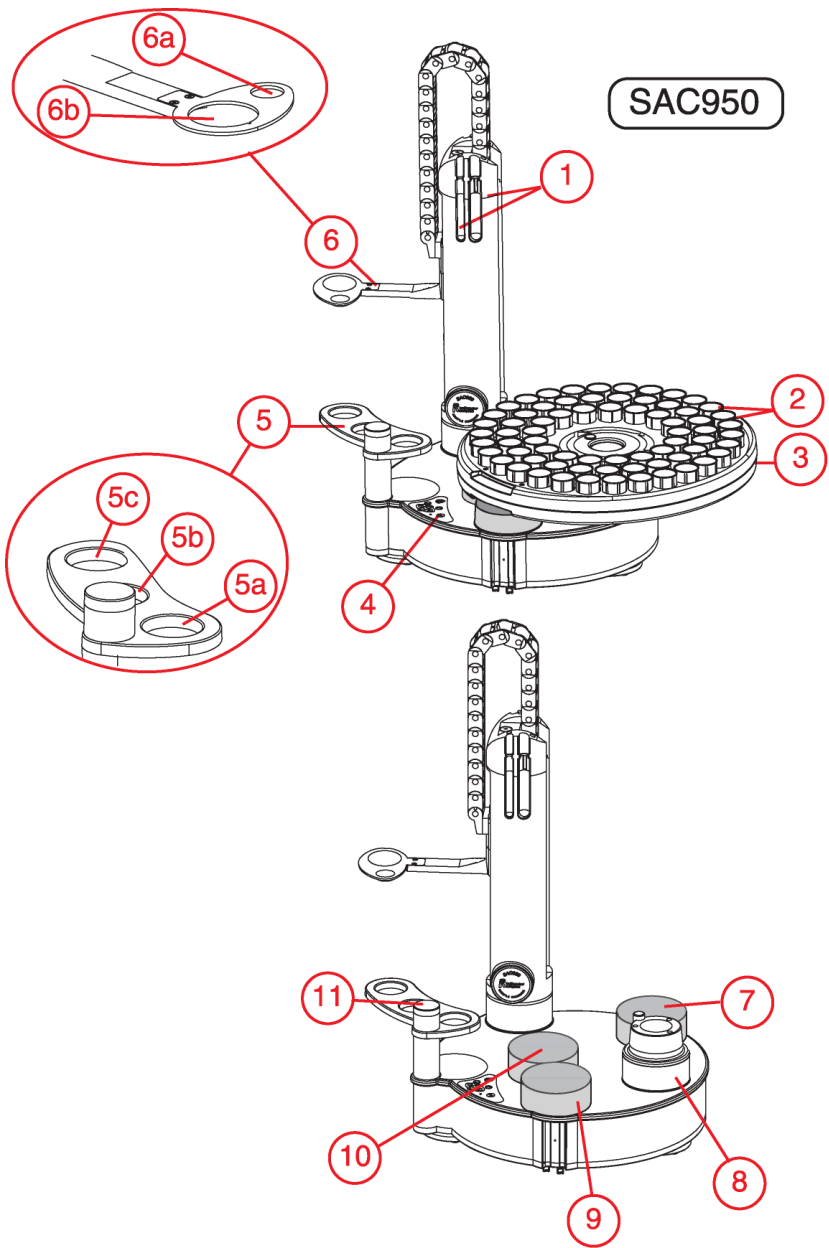
SAC950

SAC950

- ① 4 locations for spare electrodes and antidiffusion delivery tips (2 locations on each side).
- ② Sample beakers.
- ③ 1, 2 or 3 radii turntable. Capacities depend on the beaker size and diameter : see chapter "Accessories" - Tables "Turntables and beakers".
- ④ Keypad.
- ⑤ Beaker extension with Park (5a), Rinse 1 (5b) and Rinse 2 (5c) beaker locations.
- ⑥ Electrodes/Stirrer/Beaker presence detector holder.
Beaker presence detector location (6a).
Electrodes/stirrer head holder location (6b).
- ⑦ Optional peristaltic pump.
- ⑧ Turntable holder.
- ⑨ Optional shower rinse pump/emptying pump module. Comprises 2 diaphragm pumps (rinsing and emptying).
- ⑩ Magnetic stirrer assembly comprising also a RFID reader device for turntable automatic recognition.
- ⑪ Status lamp:
Blue, blinking : Free for operation.
Manual operations on keyboard are possible
Purple: The sample changer is controlled by the workstation unit (analysis in progress). Only the keyboard **Stop** key is available.
Red: An error has been detected.

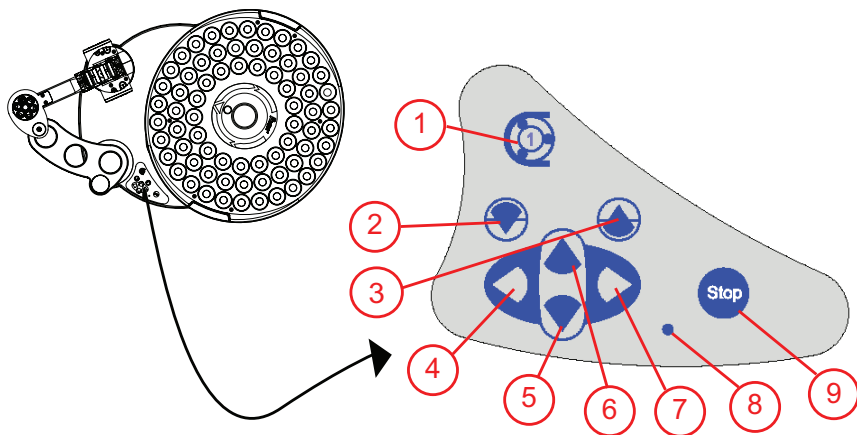Keypad

- ① Starts adding reagent with peristaltic pump (*). Stops addition when the key is latched down.
- ② Starts shower rinsing using a membrane pump of the Dynamic rinsing module (**). Stops action when the key is latched down.
- ③ Starts emptying a beaker using the second membrane pump of the Dynamic rinsing module (**). Stops action when the key is latched down.

(*) If a Reagent addition module is installed (see chapter "Setting up") and if the electrode holder is at its downmost position.

(**) If a Dynamic rinsing module is installed (see chapter "Setting up") and if the electrode holder is at its downmost position.

④ Moves to the previous beaker. If the electrodes dip into the beaker (or are above the beaker) before pressing the key, at the end of the operation, they will dip into the previous beaker (or will be above the previous beaker). The beakers are ranked as follows:

Rinse beaker 2 (R2)
Rinse beaker 1 (R1)
Park beaker
Beaker no.1 of the turntable
.....
Beaker no.n-1 of the turntable
Beaker no.n (last beaker) of the turntable
Rinse beaker 2 (R2)
.....

⑤ Moves the arm (electrode head holder) downwards.

⑥ Moves the arm (electrode head holder) upwards or moves the electrodes above the Park beaker if the electrodes are already on service position (upmost position).

⑦ Moves to the next beaker. Same observations as for the "Moves to the previous beaker" key (④).

⑧ Blue pilot lamp. Lights on when a key is pressed and lights off when the key is latched down. This lamp blinks when the key is not available (example: during an analysis, when a key other than **Stop** is pressed).

⑨ **Stop** key. Stops the action in progress while a sequence is running. Removes an error status (red status pilot lamp). The **Stop** key is the only active key while a sequence is running.

(***) Beaker no.n = last beaker of a SAC950 turntable or last sample beaker of a SAC850 turntable.

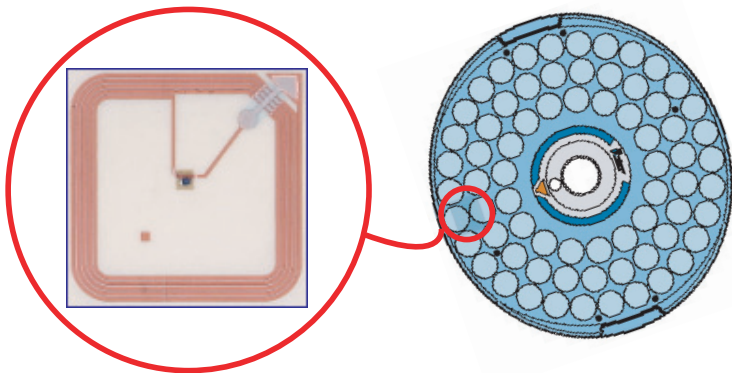
RFID technology on turntables

What's RFID means and how does it work?


Radio Frequency IDentification (RFID) is a generic term that is used to describe a system that transmits the identity (in the form of a unique serial number) of an object or person wirelessly, using radio waves. It's grouped under the broad category of automatic identification technologies.

It's the combination of a "tag" and a "reader". A typical RFID tag consists of a microchip attached to a radio antenna mounted on a substrate. The chip can store data. The reader is used to retrieve the data stored on an RFID tag. A typical reader is a device that has one or more antennas that emit radio waves and receive signals back from the tag. The reader then passes the information in digital form to a computer system.

In our assembly an RFID tag is attached inside the turntable at the first beaker level position. This tag is clearly visible.

RFID tag placed on the turntable

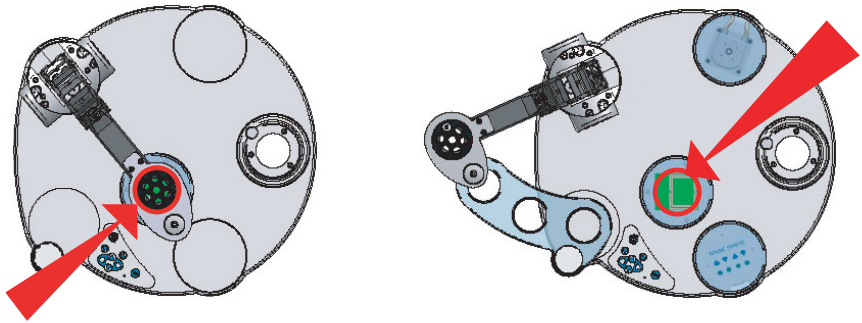*RFID reader mounted inside the Magnetic stirrer assembly
SAC850 on the left and SAC950 on the right*

The RFID reader is mounted inside the Magnetic stirrer assembly (see figure above). When an analysis is started and a turntable is in place, the sample changer rotates the tray (and the arm in the case of a SAC950) so that the tag and the reader are face to face for a moment.

Which data contains a turntable RFID tag?

The turntable identification data (type of turntable, number of beakers, beaker location and beaker size/height).

Benefits of using RFID technology on turntables

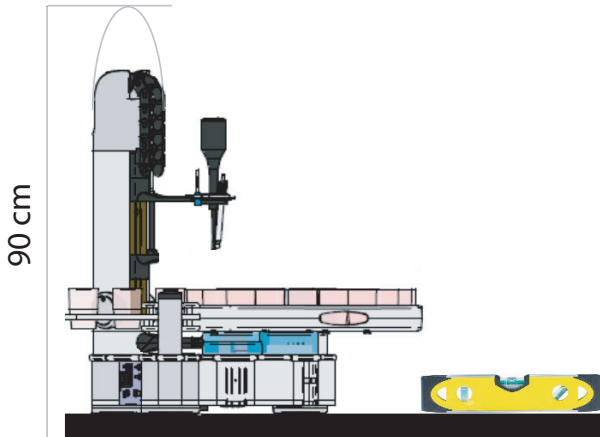
The Sample Changer knows at any moment if a turntable is installed and which turntable is used.

3. Setting up

Important: Never, at any time during setup or use of the instrument, manually move the arm or turntable even if blocked. Manual movement may cause damage to the motor axis, which in turn can cause excessive play resulting in positioning errors.

Note: Unless otherwise indicated, the following setting up instructions are common to both SAC850 and SAC950 Sample Changers.

Set up the instrument on a flat and horizontal surface and in a properly ventilated place. Use a spirit level to check the surface is flat.

The power supply connectors on the rear panel must remain accessible so that you can quickly disconnect the cables in case of emergency. The room temperature should not exceed 40°C.

The relative humidity must be between 20 and 80%.


Minimum space requirements

(SAC850): 90 cm x 70 cm x 60 cm (36" x 28" x 24")


(SAC950): 90 cm x 80 cm x 60 cm (36" x 32" x 24")

70 cm

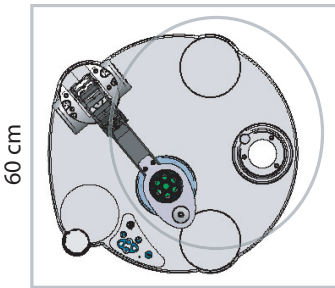80 cm


SAC850

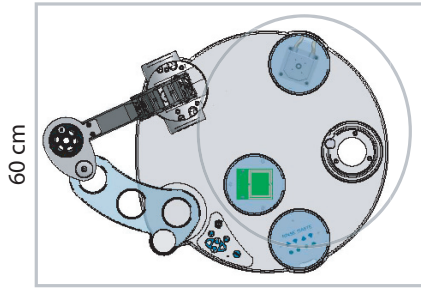SAC950

Mounting the turntable

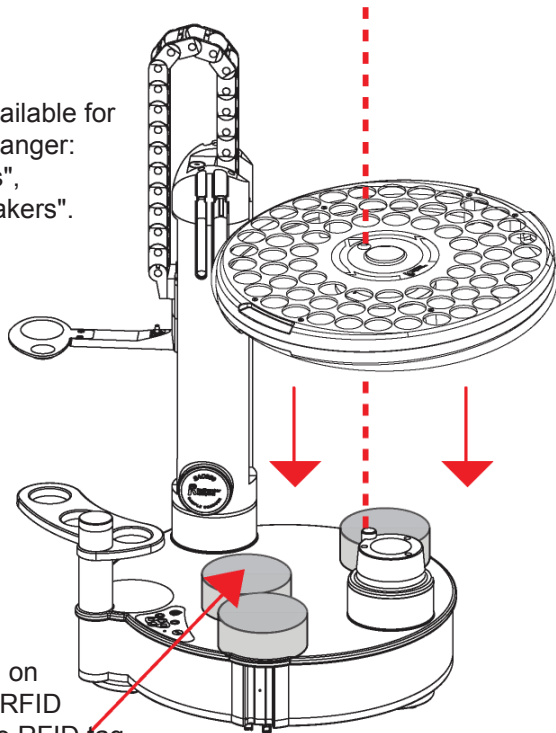Several turntables are available for each model of sample changer: see chapter "Accessories", table "Turntables and beakers".

The turntable is identified on starting an analysis. The RFID reader reads the turntable RFID tag

Placing beakers

Park beaker

At the end of a sample series, the electrodes are dipped in the park beaker or left above the park beaker. The park beaker can also be used to rinse dynamically the electrodes.

Positioning of the park beaker (see figures below):

SAC850: Last position of the turntable

SAC950: On the electrode reconditioning extension


Static rinse beakers

Between 2 sample runs, the electrodes are dipped once or twice into a rinse solution under stirring. 1 or 2 static rinse beakers can be used.

Positioning of the static rinse beaker(s)

SAC850: Last but one (**R1**) and previous (**R2**) positions of the turntable

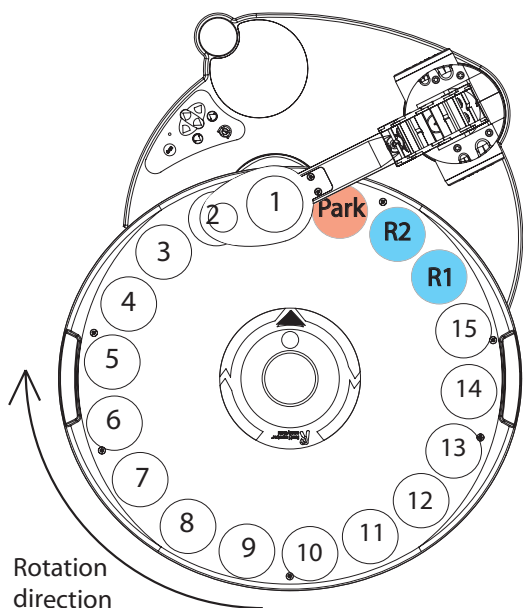
SAC950: On the electrode reconditioning extension, positions (**R1**) and (**R2**)

R1: Static rinse beaker no. 1
R2: Static rinse beaker no. 2

Other positions: Sample beakers (15 free position if 2 static rinses programmed)

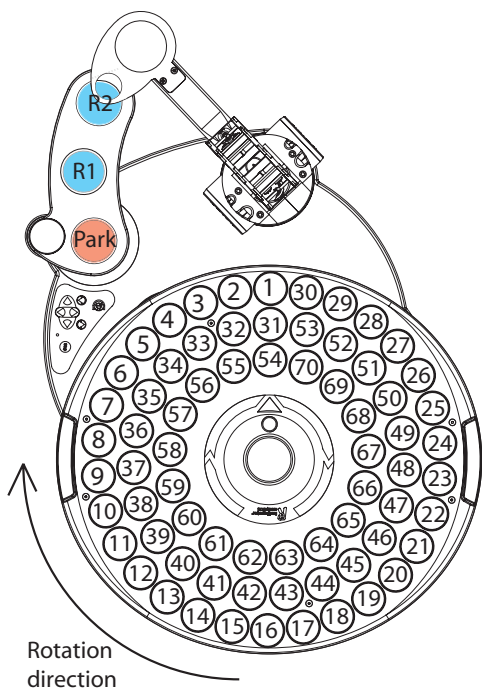Position of the beakers on a SAC850
Example of a 18-position turntable

R1: Static rinse beaker no. 1
R2: Static rinse beaker no. 2

Positions 1 to 70:
 Sample beakers

*Position of the beakers on a SAC950
 Example of a 70-position turntable*

SAC850 users:

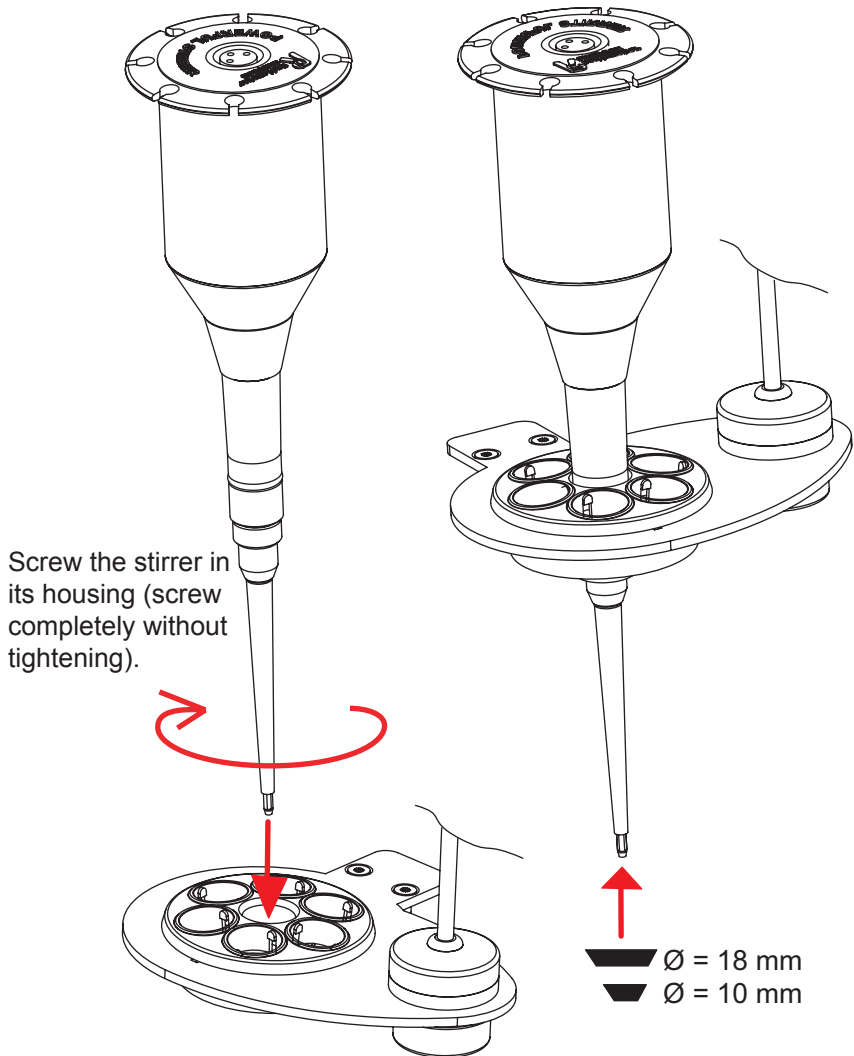
Park, static rinse and sample beaker sizes depend on the type of turntable used. These beakers are supplied with some of the turntable kits (see chapter "Accessories").

SAC950 users:

Park and static rinse beakers are 40-100 ml PP titration vessel, supplied with the SAC950, part no.: X91T055 (10 pcs).


Sample beaker sizes depend on the type of turntable used. These beakers are supplied with some of the turntable kits (see chapter "Accessories").

Mounting the propeller stirrer


The propeller stirrer is supplied with 2 propellers 10 & 18 mm. Fit the right propeller according to the size of the beaker (see chapter "Accessories", table "Turntables and beakers").

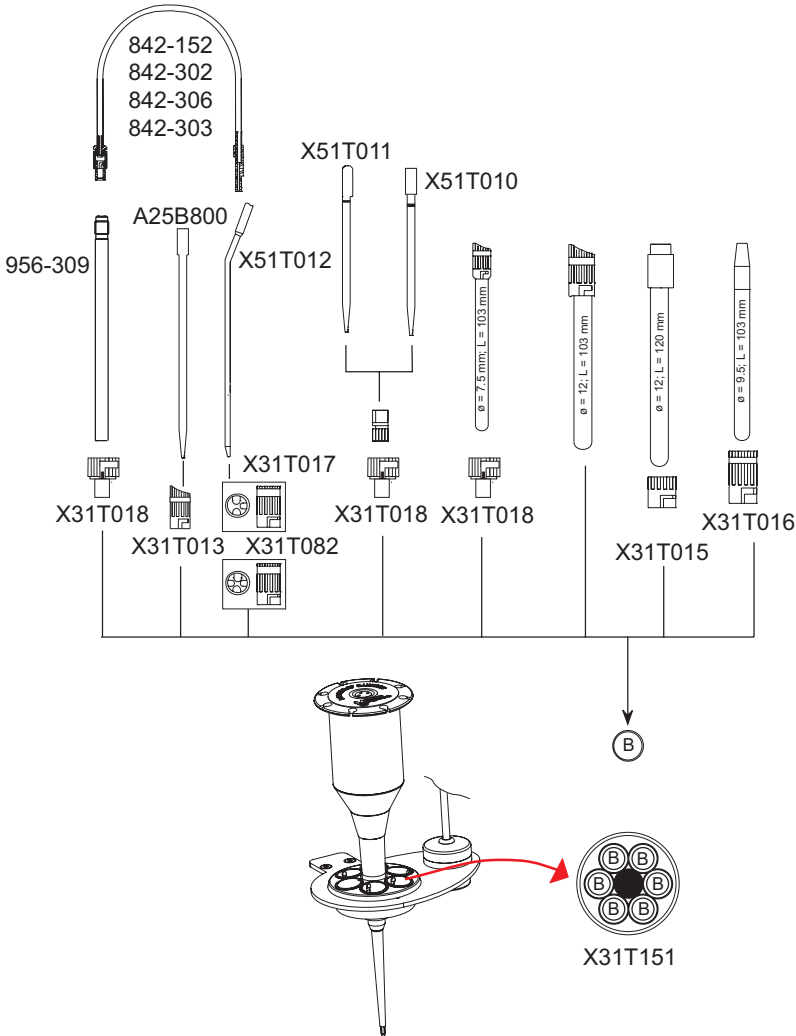Mounting the electrodes and addition tips


A set of adapters for bayonet electrode heads, part no. X91T101 is supplied with the sample changer.

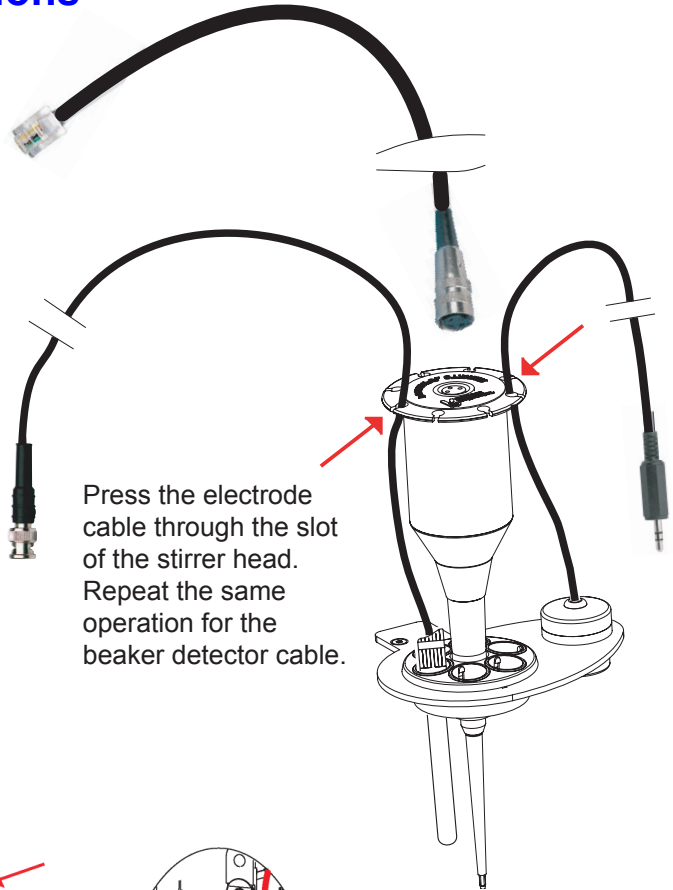Connections

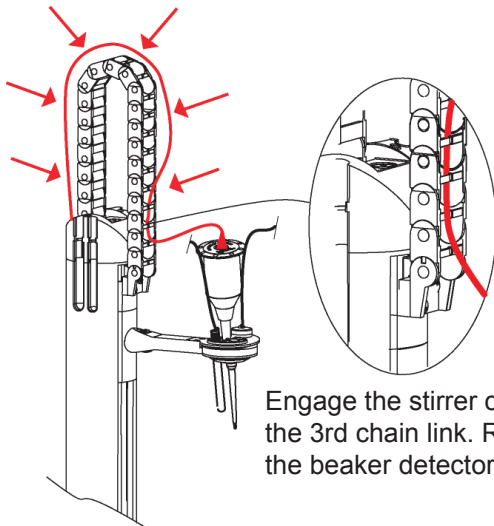
Press the electrode cable through the slot of the stirrer head. Repeat the same operation for the beaker detector cable.

Engage the stirrer cable in the cable carrier from the 3rd chain link. Repeat the same operation for the beaker detector and electrode cables.

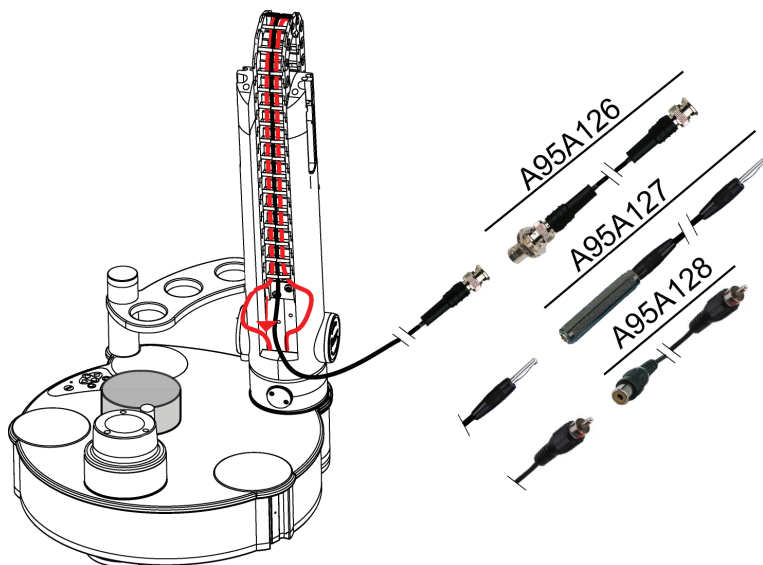
Insert the stirrer and beaker detector cables in the opening of the arm then slip these cables under the sample changer in order to present the plugs at the connection panel of the instrument. Slightly level up the instrument to facilitate this operation.

Connect the electrode the corresponding socket of the workstation unit.


- Use cable A95A126:
Electrodes with BNC plugs or screw cap electrodes + cable A94L114
- Use cable A95A127:
Electrodes with banana plugs or screw cap electrodes + cable A94L111
- Use cable A95A128:
Temperature sensors with Cinch plug

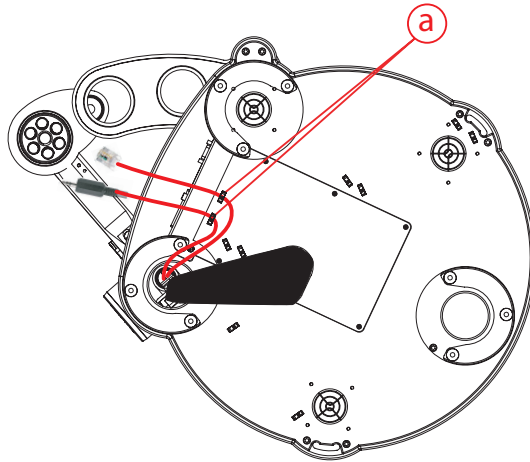Delicately tip the sample changer onto the side.
While doing this operation, do not pick up the instrument by the electrode holder or by the arm.
Slot the stirrer and beaker detector cables in holders **(a)**.

Put the sample changer back in place. While doing this operation, take care to not pinch the stirrer and beaker detector cables.

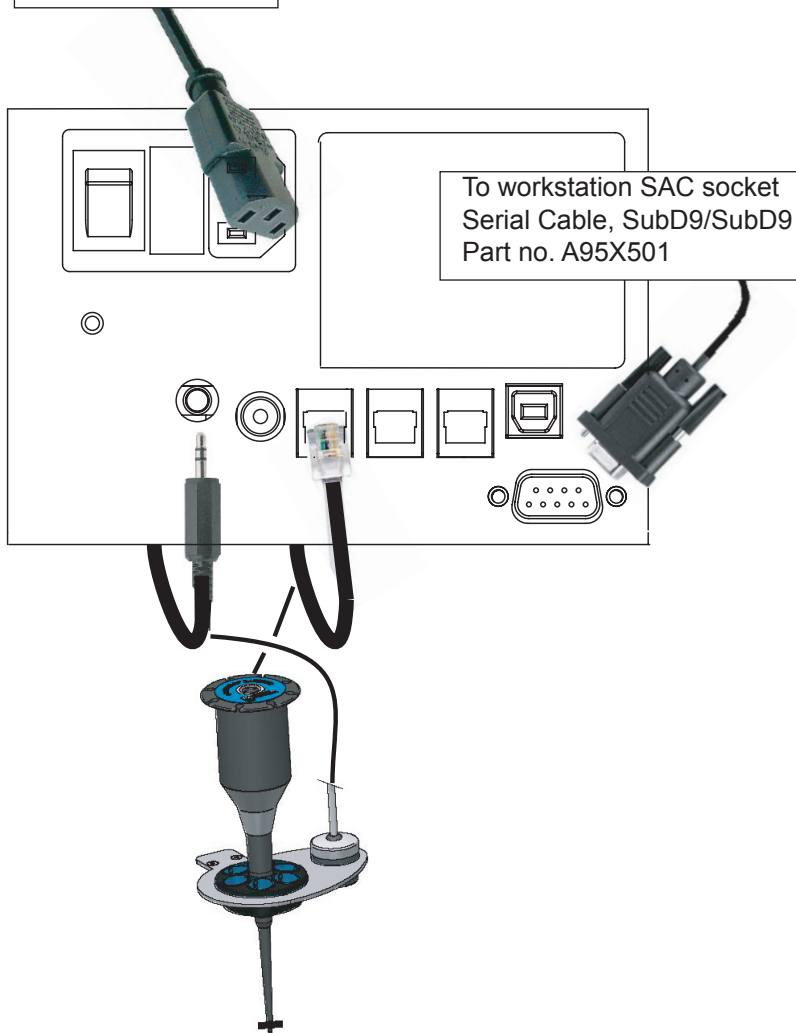To a mains socket.
Mains cable,

To workstation SAC socket
Serial Cable, SubD9/SubD9
Part no. A95X501

Assembling option accessories

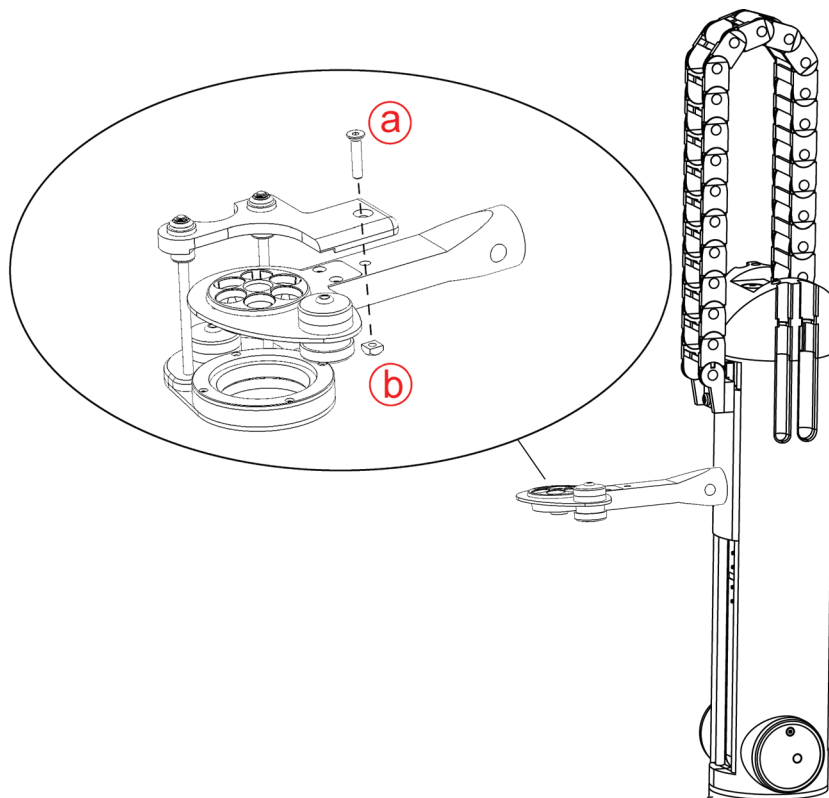
Dynamic rinse module

The Dynamic rinse module is used for shower rinsing and emptying operations. It consists of a pump assembly (two membrane pumps, one for rinsing and the other for emptying), a shower head, a suction tip and silicon tubings.

Switch on the sample changer and move the arm at its upmost position: press key  (see chapter "Presentation - Keypad", key (6)).



Switch off the sample changer.

Fix the module to the arm using the bolt (a), nut (b) and Allen key supplied.


Fix the shower head skirt **(g)** or **(h)** to the base of the module. The use of a **(g)** or **(h)** shower head skirt depends on the diameter of the beaker used (see table below).


Beakers	Use of parts (g) , (h)
400 ml tall form, d = 70 mm	(g)
250 ml low form, d = 70 mm	(g)
250 ml tall form, d = 60 mm	(g)
180 ml PP, d = 54 mm	(g)
150 ml tall form, d = 60 mm	(g)
150 ml low form, d = 62 mm	(g)
40-100 ml PP, d = 60 mm	(g)
100 ml low form, d = 48 mm	(h)
50 ml low form, d = 44 mm	(h)
22-45 ml, d = 44 mm	(h)
10-22 ml, d = 44 mm	(h)

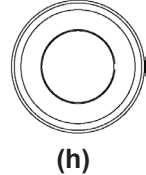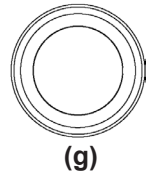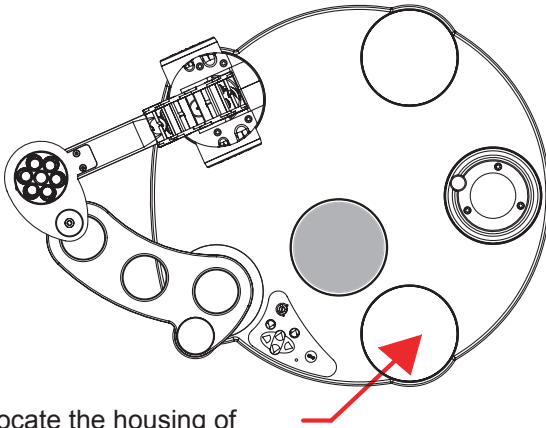If not already done,
remove the turntable.


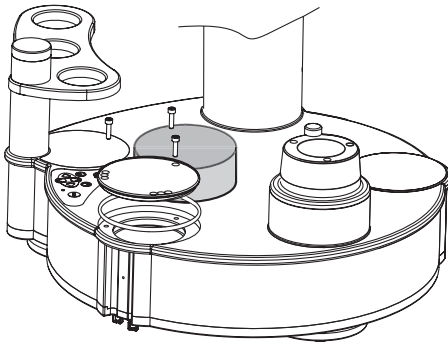
Locate the housing of
the pump assembly.


Remove the protection sticker.
Unscrew the 3 screws then
remove the cover plate. Use
the Allen key supplied.

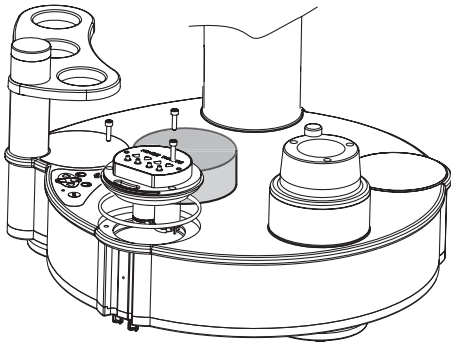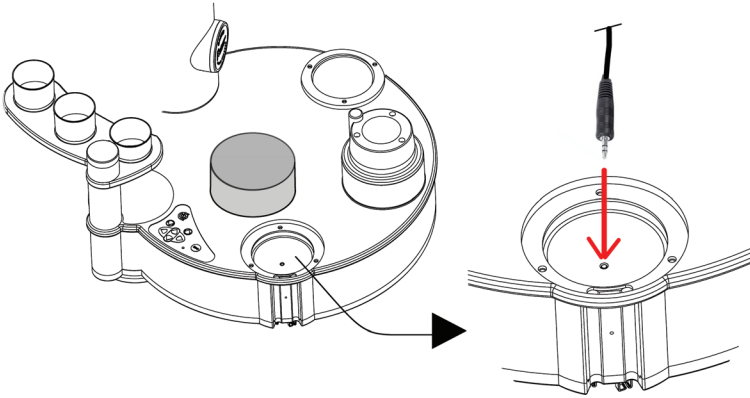
Connect the pump cable Jack plug to the corresponding socket.

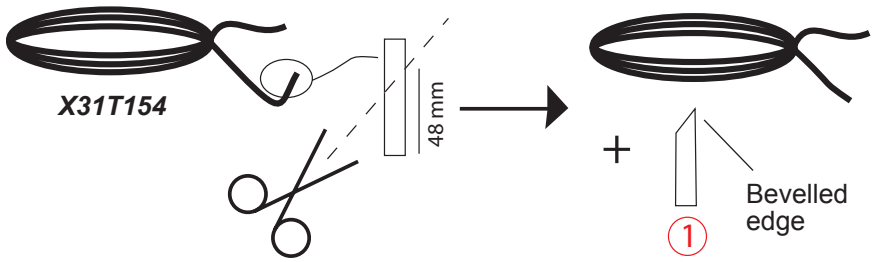
Insert the pump assembly and fix it in place with the 3 screws. Use the Allen key supplied.

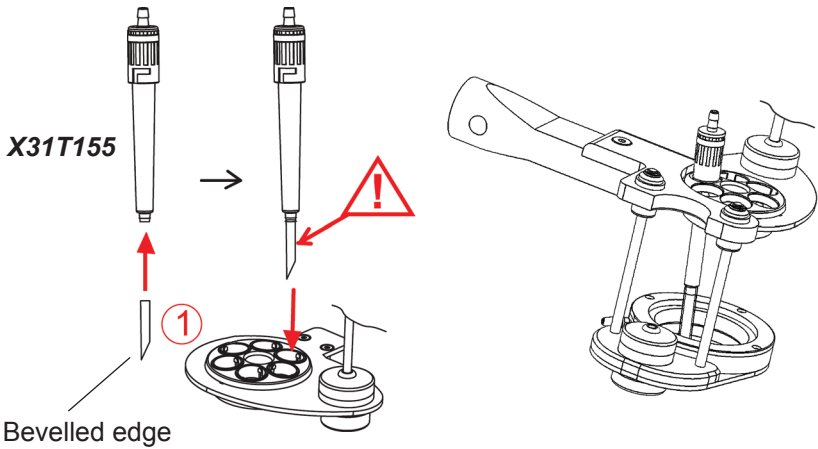
Cut a piece ① of 48 mm from the silicon tubing X31T154.

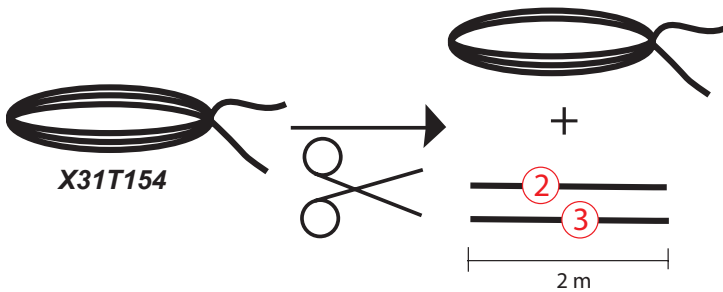Pay attention to the stirrer rod and propeller when setting up the suction tip X31T155. If necessary, direct the end of the tubing ①.

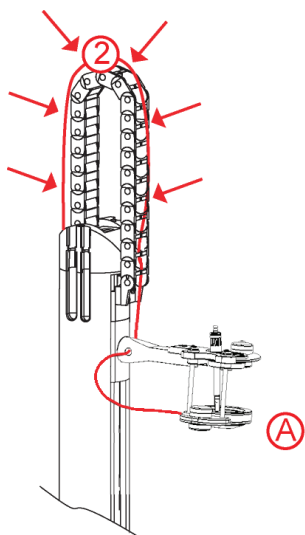Cut 2 pieces ② and ③ of 2 m each from the silicon tubing X31T154.


Mount one end of tubing (2) on the metal rod of the shower head.


Engage the tubing in the cable carrier by simple pressure on each link.

Important

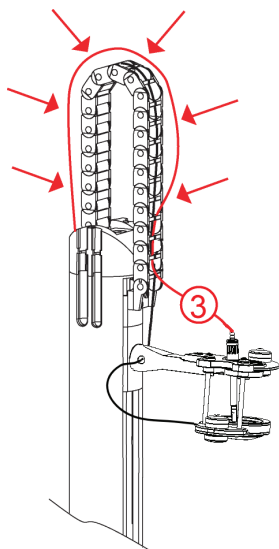
It is important to check at this stage that the tubing slides well when moving up part (A), this particularly after having set up all tubings and wires. To move up part (A), push up under this part.

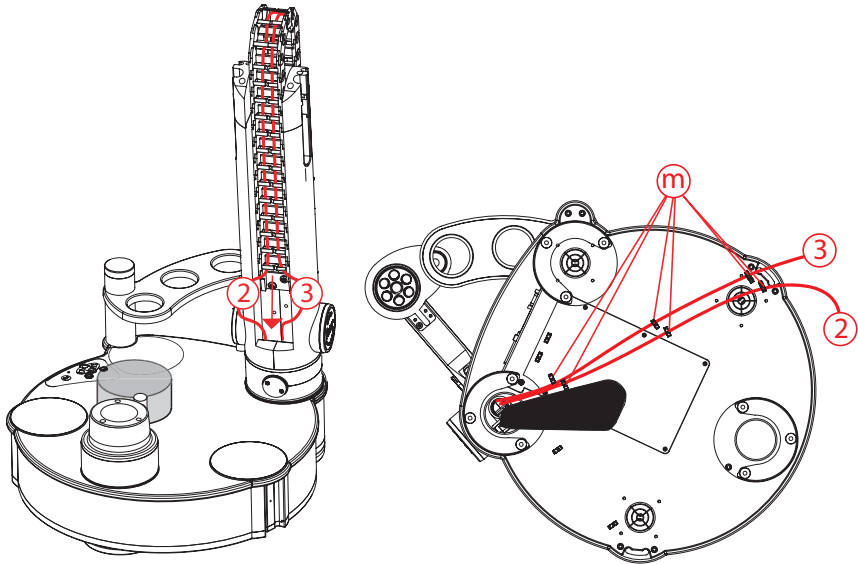
Mount one end of tubing (3) on the head of the suction tip X31T155. Engage the tubing in the cable carrier from the 3rd chain link.

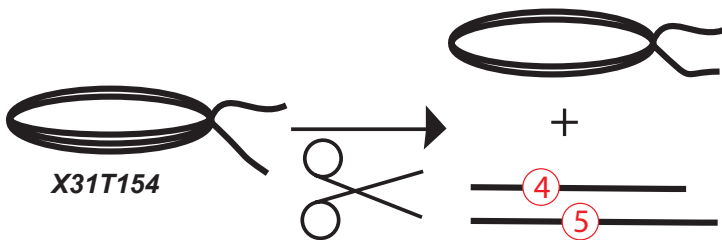
Insert tubings ② and ③ in the opening of the arm then slip these tubings under the sample changer in order to present them at the rinsing/draining pumps. The tubings must pass on the side of the black cable guide part. Delicately tip the sample changer onto the side to facilitate this operation. Slot the tubings in holders (m).

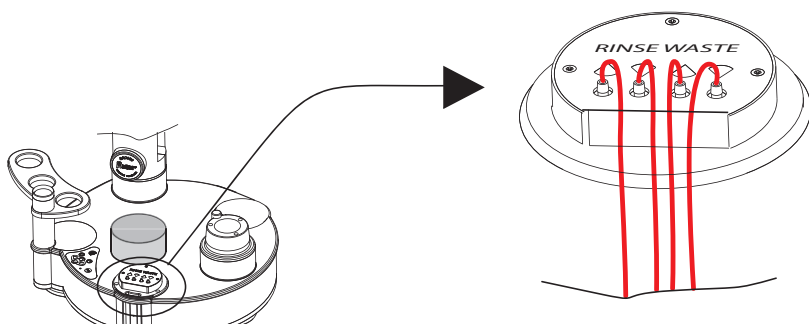Cut 2 pieces ④ and ⑤ from the silicon tubing X31T154. ④ is the pump/rinsing solvent can tubing and ⑤ is the pump/waste tubing. Determine the correct length of these tubings according to your needs.

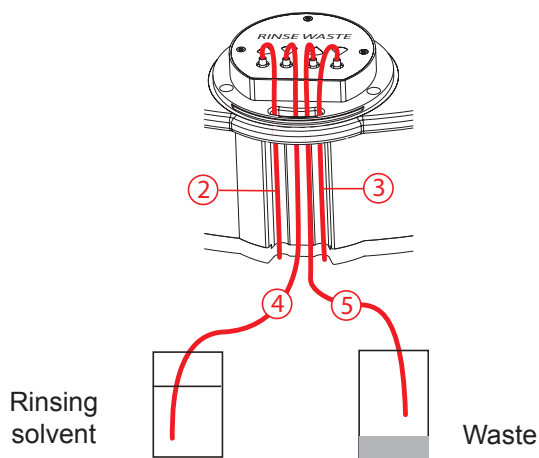
Adjust the length of the tubings to preserve a small loop after connection.

Mount tubing ② on the outlet of the rinsing pump and tubing ③ on the inlet draining pump.

Mount one end of tubing ④ on the inlet of rinsing pump and dip the other end in a can filled with rinsing solvent (demineralised water for example).

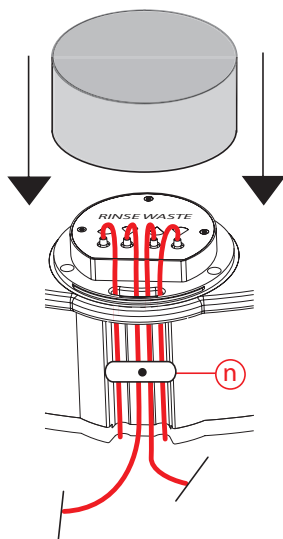Mount one end of tubing ⑤ on the outlet of draining pump and dip the other end into a waste.

Cover the pump assembly with the protection lid provided.

Fix the part (n) using the screw and Allen key 2.5 mm provided.

This part maintains the tubings in place.

Warning!

The pump body is made of polypropylene. So use the pump only with reagents compatible with polypropylene.

You can aspire precipitates with the pump to the only condition that this precipitate is resulting from a chemical reaction and that the particle size does not exceed 125 μ m.

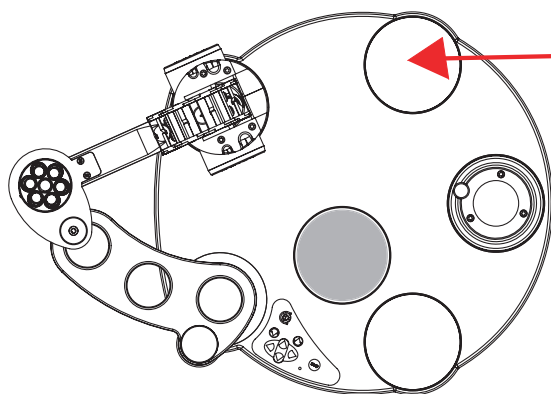
Moreover, it is recommended to carry out several cycles of rinsing in the event of samples charged in particles.

Reagent addition module


A reagent addition module can be mounted on the SAC850 or SAC950 sample changer. This peristaltic pump is used to add reagent as be part of a sample preparation. Volumes are added with an accuracy around 1%.

A reagent addition module consists of a peristaltic pump assembly, a reagent addition tip, silicon and Maprene[®] tubings.

If not already done, remove the turntable.



Locate the housing of peristaltic pump.


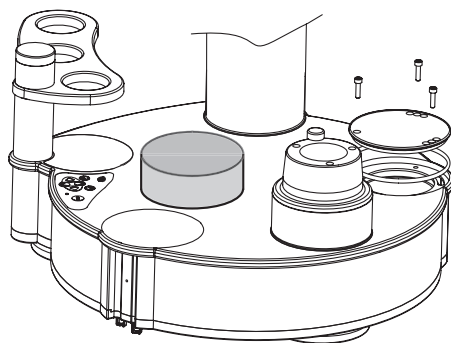
Remove the protection sticker, Unscrew the 3 screws then remove the cover plate. Use the Allen key supplied.

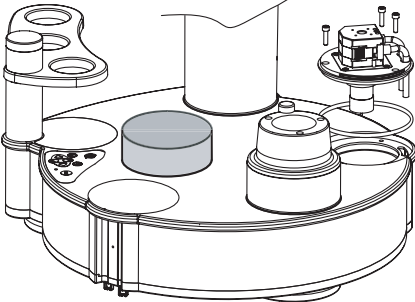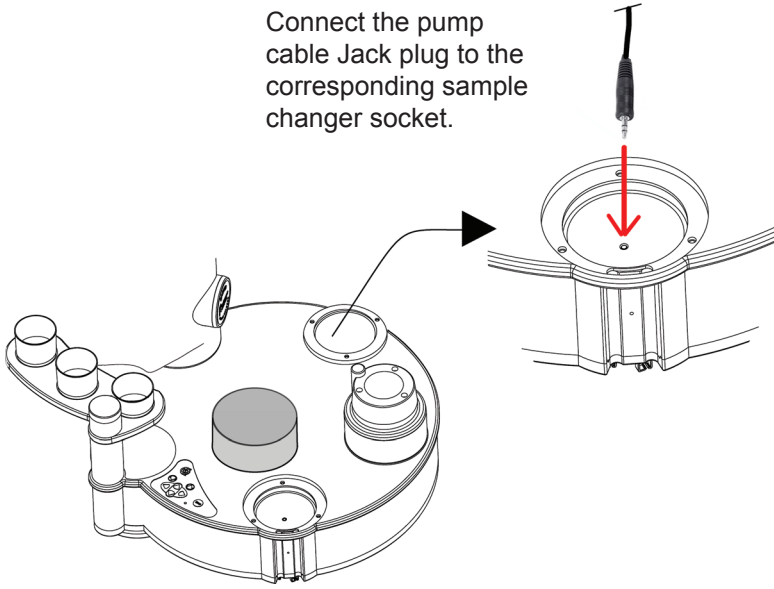
Connect the pump cable Jack plug to the corresponding sample changer socket.

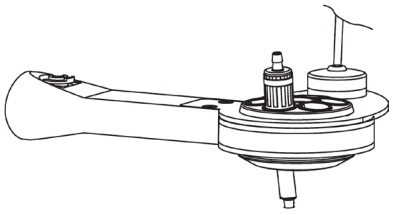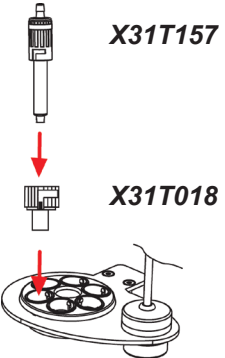
Insert the peristaltic pump assembly and fix it in place with the 3 screws. Use the Allen key supplied.

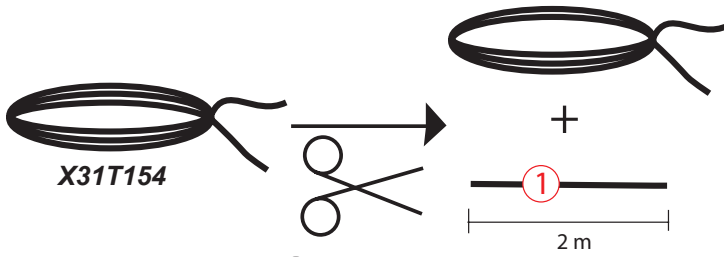
Set up the addition tip, part no. X31T157.

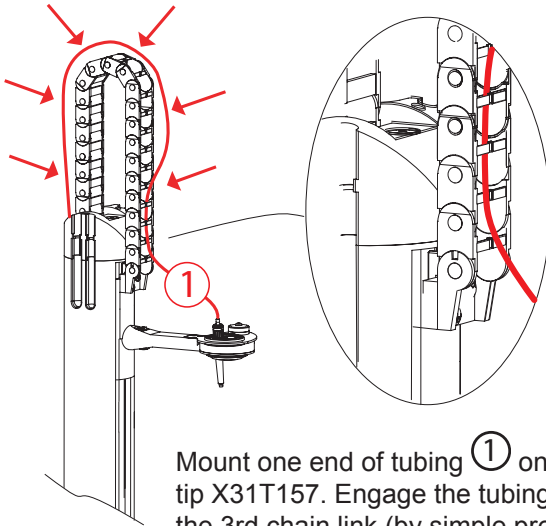Cut a piece of 2 m long ① in the silicon tubing X31T154.


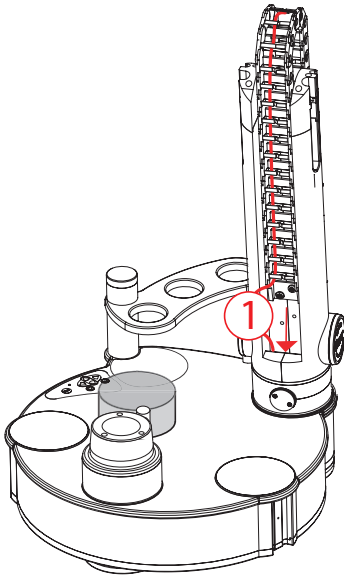
Mount one end of tubing ① on the head of the addition tip X31T157. Engage the tubing in the cable carrier from the 3rd chain link (by simple pressure on each link).

Insert tubing ① in the opening of the arm then slip this tubing under the sample changer in order to present it at the peristaltic pump. The tubing must pass on the side of the black cable guide part.


Delicately tip the sample changer onto the side to facilitate this operation.

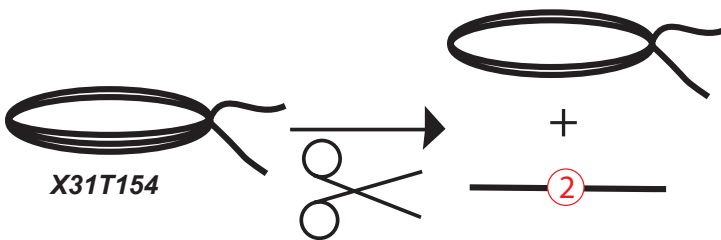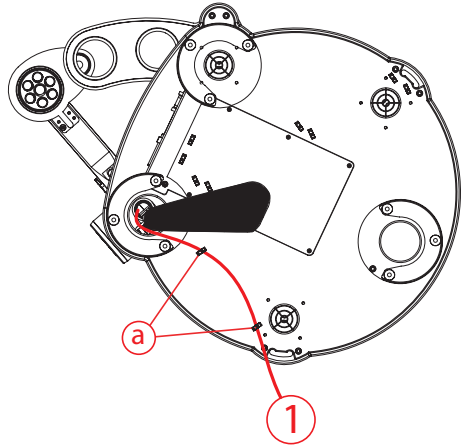Slot the tubings in holders (a).


Cut a piece ② from the silicon tubing X31T154.

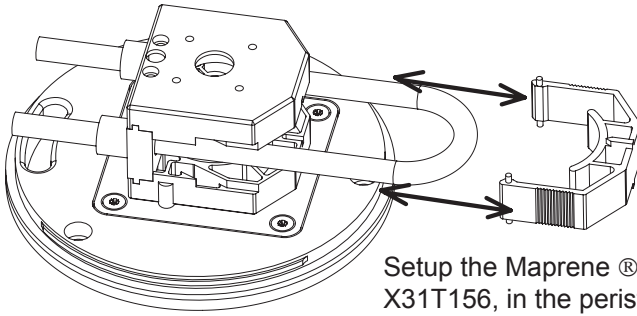② is the pump/reagent flask tubing. Determine the correct length of this tubing according to your needs.

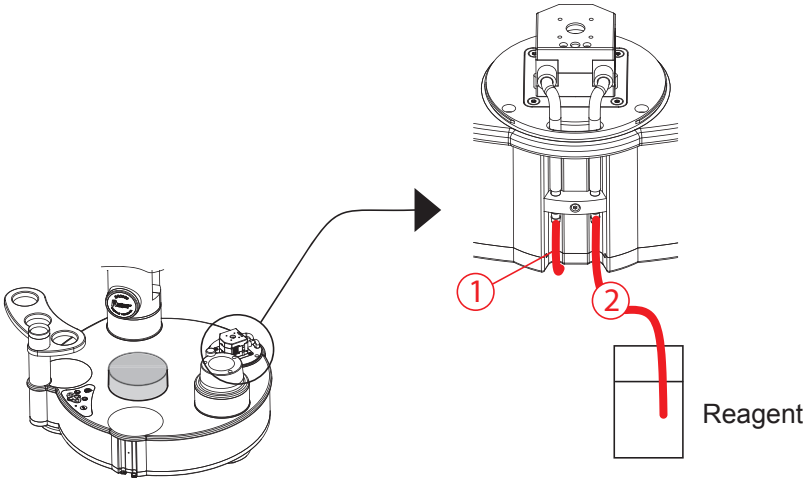
Setup the Maprene[®] tubing, part no. X31T156, in the peristaltic pump.


Mount tubing ① on the outlet of the peristaltic pump.

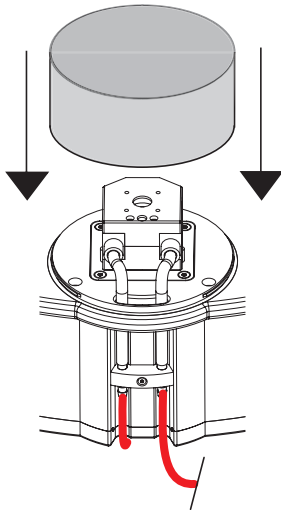Mount one end of tubing ② on the inlet of peristaltic pump and dip the other end in the reagent flask.

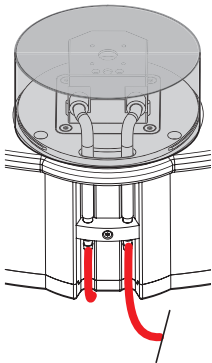Cover the pump with the protection lid provided.

Beaker cover module

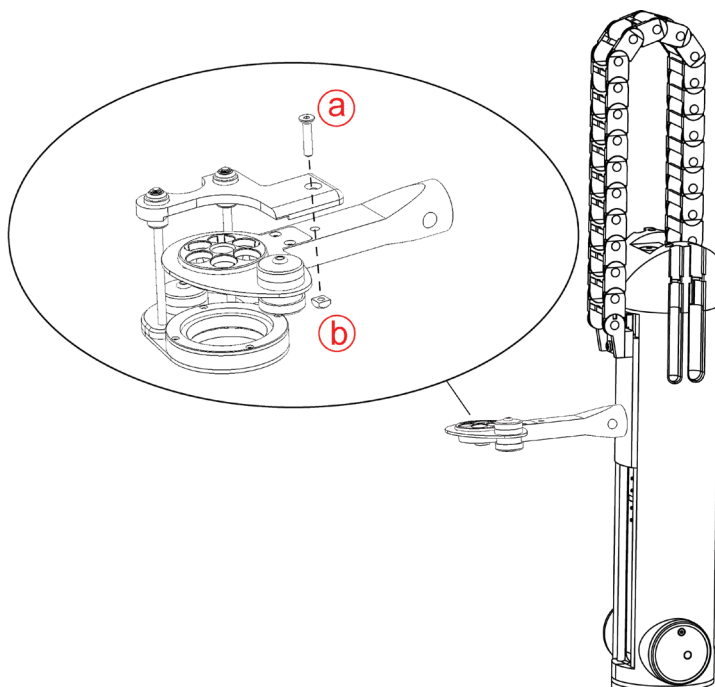
The Beaker cover module for SAC950, X16T016 is used to remove metal covers from any beakers (*) of the SAC950 turntable. The Beaker cover module cannot work without the Dynamic rinse module.

Another application of the Beaker cover module is the automatic detection (by the beaker detector) of beakers with solid samples, powder samples or liquid samples with heights of liquid lower than the limit of detection (10 mm).


Switch on the sample changer and move the arm at its upmost position:
press key  (see chapter "Presentation - Keypad", key **(6)**).

Switch off the sample changer.

Fix the module to the arm using the bolt (a), nut (b) and Allen key supplied.

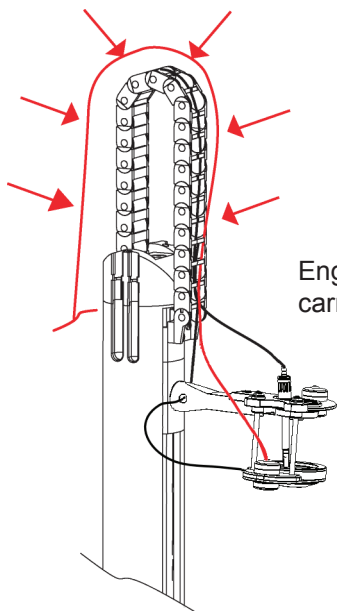
(*) Metal covers are listed in chapter "Accessories", table "Beaker and beaker covers".


Engage the electromagnet cable in the cable carrier from the 3rd chain link.

Delicately tip the sample changer onto the side. While doing this operation, do not pick up the instrument by the electrode holder or by the arm.

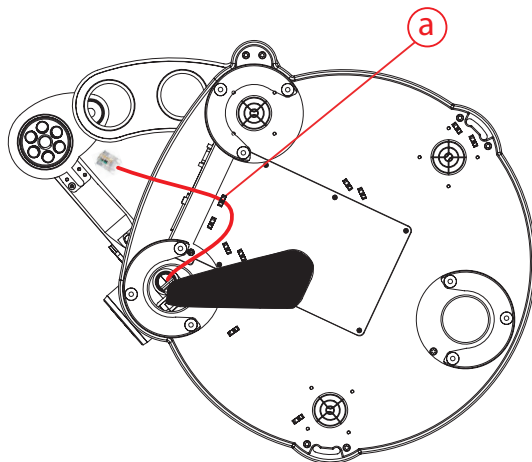Slot the electromagnet cable in holder **(a)**.

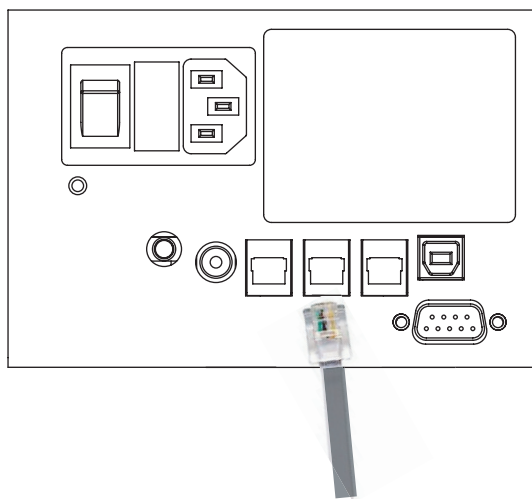Put the sample changer back in place. While doing this operation, take care to not pinch the cables and pump tubings.

Connect the electromagnetic cable to the corresponding socket of the SACxxx connection panel.

4. Use

Switching on

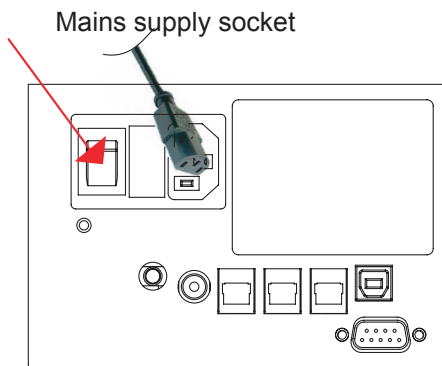
1. Connect the mains socket of the SAC 850/SAC950 sample changer and the workstation (TIMxxx or IONxxx) to the mains supply using the 3-lead power cord provided.
In the USA or Canada, use a UL listed power cord only.

Important:

The sample changer and the workstation must be connected to an earthed mains socket for safety reasons. Efficient grounding is essential for reliable measurements and security.

The power supply connector on the rear panel must remain accessible and close to the user (2 m maximum) so that you can quickly disconnect the cables in case of emergency.

2. Switch on the sample changer (O/I power switch on the rear panel set to "I", see below) then the workstation.

The sample changer runs its autotest routine:

- Lifts/lowers the arm to search for the arm upmost position,
- Starts the turntable rotation motor (SAC850 and SAC950) and start the arm rotation motor (SAC950 only) to search for reference positions of both elements.
- Dips the electrodes into the Park beaker. If no Park beaker detected, leaves the electrodes above the Park beaker (Service position).

Running sequence with a workstation

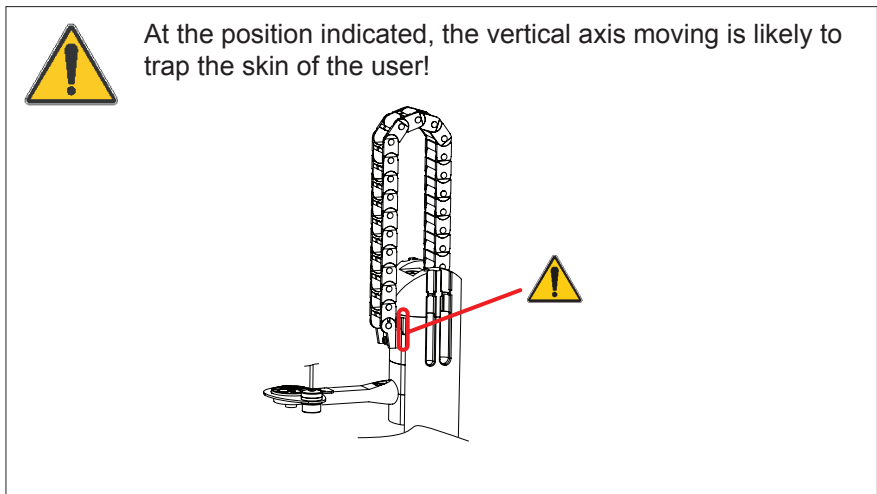
After having started a sequence from the workstation or TitraMaster 85, the sample changer performs the following operations if 1 to 9 dynamic rinses and if 2 static rinses R1 and R2 have been programmed (*):

1. Moves to Park beaker (**), and starts rinsing dynamically the electrodes. At the end, the Park beaker is left empty.
See paragraph "Dynamic rinsing".
2. Moves to R1 beaker (***), dips the electrodes into that beaker and rinses under stirring during a preset delay. *See paragraph "Stirring".*
3. Moves to R2 beaker (***), dips the electrodes into that beaker and rinses under stirring during a preset delay. *See paragraph "Stirring".*
4. Moves to beaker no.1, prepares sample/standard running the sample preparation number attached to the method. During this time, the workstation displays the message "Sample changer processing in beaker 1" message.
See paragraph "Sample preparation".
5. At the end of analysis, starts rinsing dynamically the electrodes. At the end, the beaker no.1 is left nearly empty with a small amount of

solvent that was used to flush the electrodes and delivery tips.

- (*) The number of static rinses (0, 1 (R1) or 2 (R1 and R2)), the duration of static rinses and the number of dynamic rinses (0 to 9) are programmed in the Setup > Configuration parameters of the workstation (refer to the Reference Manual of the workstation used).
- (**) A dynamic rinse before first beaker analysis can occur in the Park or in the R1 beaker if 0 or 1 static rinse has been programmed. Selection between Park and R1 is made in the Setup > Configuration parameters of the workstation (refer to the Reference Manual of the workstation used).
- (***) R1: Static rinse beaker no.1. R2: Static rinse beaker no.2. The position of these beakers is given in chapter "Setting up - Placing beakers".

6. Repeats steps 2 and 3.
7. Dips the electrodes in beaker no. 2, prepares sample/standard running the sample preparation number attached to the method.
8. At end of the sequence (end of analysis in the last beaker of the sequence):
 - starts rinsing dynamically the electrodes,
 - repeats steps 2 and 3,
 - moves the electrodes above the Park beaker,
 - left the electrodes as is or dip the electrodes into the Park beaker depending on the selection made in the Setup > Configuration parameters of the workstation (refer to the Reference Manual of the workstation used).

Dynamic rinsing

1 to 9 dynamic rinses (or shower rinse/emptying operations) can be performed between 2 sample runs. To improve rinsing efficiency and to save on rinse solution, the rinse process is a pulsed rather than a continuous operation.

A dynamic rinse is carried out in the just analysed sample beaker or in the Park beaker (*). At the start of a sequence, the dynamic rinse can be performed either in the Park beaker or in R2 (static rinse no.2) beaker (*). Dynamic rinses never take place in calibration beakers. At the end of a calibration beaker analysis, the dynamic rinse is performed in the Park or R2 (static rinse no.2) beaker (*).

A dynamic rinse performed in a Park or Rinse beaker consists of the following operations:

1. The electrodes are positioned above the Park or Rinse beaker.
2. The electrodes are dipped into the beaker. The beaker is emptied to a waste in the same time. At the end, the electrodes are located in the emptied beaker at their downmost position.
3. The electrodes are washed with rinse solution (usually demineralised water) then start to move up under rinsing. At the end, the electrodes are located above a beaker filled with rinse solution and some remaining impurities.

Steps **2** and **3** can be repeated up to 8 times as up to 9 dynamic rinses can be programmed. Steps **2** and **3** are performed under stirring.

- (*) Depending on the selection made in the Setup > Configuration parameters of the workstation (refer to the Reference Manual of the workstation used).

A dynamic rinse performed in a Sample beaker consists of the following operations:

1. The electrodes are dipped into the Sample beaker. The beaker is emptied to a waste in the same time. At the end, the electrodes are located in the emptied beaker at their downmost position.
2. The electrodes move up and are washed in the same time with rinse solution (usually demineralised water). At the end, the electrodes are located above a beaker filled with rinse solution and some remaining impurities.
3. The sample beaker is emptied to a waste.
4. A last rinsing of the electrodes and delivery tips is carried out. At the end, the electrodes are located above a nearly emptied beaker containing a little solvent that was used to flush the electrodes and delivery tips.

Steps **1** to **3** can be repeated up to 8 times as up to 9 dynamic rinses can be programmed. Steps **1** to **3** are performed under stirring.

Stirring

A magnetic and a propeller stirrer can be used to stir the solutions. When the propeller stirrer is connected (see chapter "Setting up"), it will be automatically used in place of the built-in magnetic stirrer. The magnetic or propeller stirrer speed is set in the Cell menu of the workstation (value adjustable between 100 and 1100 rpm by steps of 50 rpm).

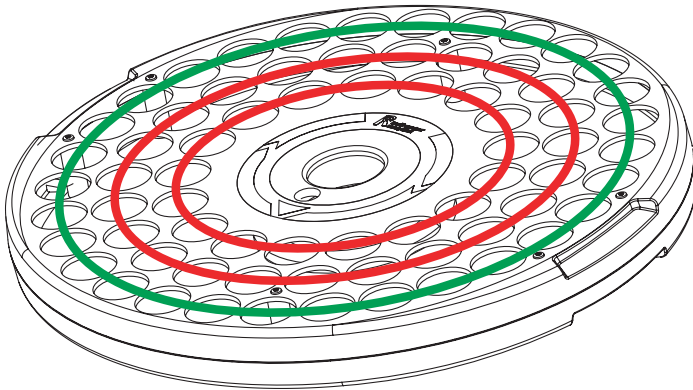
Dynamic and static rinses are performed under stirring.

Magnetic stirring with a SAC950

When using a 2 or 3-radii turntable with a SAC950, you must use a propeller stirrer to stir the solutions that are placed on the smallest row (2-radii turntable) or on the 2 smallest rows (3-radii turntable).

Green positions : magnetic and propeller stirring possible

Red positions : propeller stirring only

Sample preparation

A SAC850 or SAC950 can run up to 10 different sample preparations in a sequence. Sample preparations are created and edited from TitraMaster 85 PC software.

Each element of a sample changer sequence is associated to a sample preparation. A sample preparation is a succession of up to 10 elementary actions required by the sample changer to prepare each sample beakers to be run thereafter.

On running a sequence, the workstation sends to the sample changer all sample preparations to be runned for the sequence. The compatibility between the sample preparation and the sample changer used is tested and an error message is displayed in case of failure (example: use of a peristaltic pump that is not installed yet).

See *TitraMaster 85 on line help, topics "Handling sample preparations"*.

Important:

When starting a sample preparation, if you have a solid or powder sample in the beaker or if you have, for some beakers, less liquid than a minimum height limit (see chapter "Accessories - Beakers"), then you must disable the "beaker presence detection" option before running the sequence. If you keep that option activated, then the sample changer will not prepare the sample in that beaker and will jump to the next beaker considering that the current position is empty.

To disable the "beaker presence detection" option in TitraMaster 85, see TitraMaster 85 on line help, topics "Editing an application - Configuration".

With a SAC950 sample changer, you can detect beakers containing solid samples, powder samples or liquid samples with heights of liquid lower than the limit of detection (10 mm). In that case, you must install the Beaker cover module on the sample changer and cover the beakers with the appropriate metal covers (see chapter "Accessories", table "Beakers and beaker covers").

Actions that can be done during a sample preparation:

Control peristaltic pump.

Start/stop electrode rinsing.

Start/stop beaker emptying.

Control stirring.

Set delay.

Control stirrer.

TTL command (wait for a TTL IN for a preset time).

5. Troubleshooting

A number of tests are automatically performed when the sample changer is switched on and during calibration and sample batch measurements. The errors reported can be divided in two groups:

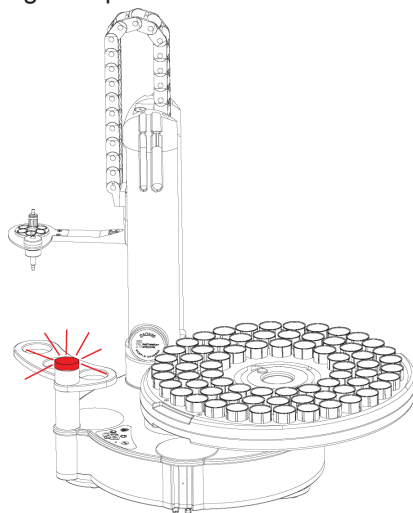
- Autotest errors occurring on switching on the sample changer and errors occurring while operating the sample changer alone (manual mode without workstation).
- Errors displayed by the workstation.

Autotest and manual mode errors

If an error is detected, the sample changer stops the current action and the status lamp lights in red.

Switch off the sample changer, identify the causes of the error (generally, arm or turntable blocked or arm or turntable movement obstructed), correct the error then switch on the instrument.

It should be noted that when no movement is in progress, the simple fact of manually moving the arm or the turntable is interpreted by the apparatus like a movement forced resulting in an error with a lit red status lamp.

Important: Never, at any time during setup or use of the instrument, manually move the arm or turntable even if blocked. Manual movement may cause damage to the motor axis, which in turn can cause excessive play resulting in positioning errors.

Electrode head not centered correctly above the beaker

Adjust the reference position of the tray and arm rotations.

See chapter 6 "Maintenance", paragraph "Adjusting the mechanics reference positions".

Errors displayed by the workstation

Messages	Error/Action
Communication failure	<p>Data transmission between the sample changer and the workstation cannot be performed properly.</p> <p>Check the cable connections and make sure that the sample changer is switched on and connected to the workstation via the RS232 serial cable.</p> <p>Press the workstation key 3 or Stop and restart the sequence.</p> <p>It is not possible to continue the sequence from the point it stopped.</p>

Messages	Error/Action
<p>Missing beaker</p>	<p>Beaker not detected. Solve the problem and restart the sequence from the beaker it stopped (key 1 Resume analysis).</p> <p>Warning! <i>Do not change the turntable before restarting the sequence from the beaker it stopped (key1) or before restarting the sequence from the next beaker (key 2). The sample changer identifies a turntable only when a new sequence is initialized (equivalent to a keystroke on 3 End of sequence followed by 1 Run sequence).</i></p> <p>Causes:</p> <ol style="list-style-type: none"> 1. There is no beaker at the dedicated position and/or there is not enough liquid in the beaker and the beaker cannot be detected. <i>Refer to chapter "Accessories - Beaker" to know the minimum height of liquid required for detection.</i> Add solution in the beaker or deactivate the beaker automatic detection by clearing the option "Beaker detection" in TitraMaster 85 (refer to TitraMaster 85 on-line help, topics "Editing an application - Configuration"). 2. Using a level, check that the sample changer is installed on a flat and horizontal bench surface. 3. If points 1 and 2 have failed, adjust the beaker detector position (see chapter 6 "Maintenance").

Messages	Error/Action
<p>SAC arm obstructed or Turntable blocked</p>	<p>The arm or turntable has been blocked and cannot function properly. Remove the obstruction and press Resume analysis (key 1) to continue the sequence from the point it stopped.</p> <p>Warning! <i>Do not change the tray before restarting the sequence from the beaker it stopped (key1) or before restarting the sequence from the next beaker (key 2). The sample changer identifies a turntable only when a new sequence is initialized (equivalent to a keystroke on 3 End of sequence followed by 1 Run sequence).</i></p>
<p>Tray missing</p>	<p>On starting a sequence, the sample changer cannot read the RFID tag of the turntable. Check that a turntable is correctly mounted on the sample changer. Check the model of turntable used and see chapter "Accessories".</p> <p>In particular, note that a SAC850 turntable can be used on a SAC950 but a SAC950 turntable cannot be used on a SAC850. Solve the problem then restart the sequence.</p>
<p>SAC switch Off/On or ERR#32</p>	<p>Either the data transmission between sample changer and the workstation cannot be performed properly, in which case you should check the cable connections, or the measurement stopped due to a movement error. End the sequence by pressing key Stop on the sample changer keypad or key Stop of the workstation, and check that nothing is obstructing sample changer movements.</p>

Messages	Error/Action
SAC option missing	<p>The SAC option addressed is missing and not installed on the sample changer.</p> <p>Examples:</p> <p>1) Dynamic rinses are programmed in the sequence and no dynamic rinse module is installed and the sample changer. Review the Setup > Configuration parameters and/or install missing pump (in particular, check the electrical connection between the pump and the sample changer).</p> <p>2) A reagent addition is programmed in the sample preparation and no peristaltic pump is installed on the sample changer. Review the Edit sequence parameters, parameter Sample preparation no. and/or install missing peristaltic pump (in particular, check the electrical connection between the pump and the sample changer). <i>For TitraMaster 85 users only, check and, if necessary, edit the sample preparation routine.</i></p>
Wrong type	<p>The sample changer connected does not correspond to the one programmed in the Setup > Configuration menu.</p> <p>Example:</p> <p>A SAC850 is connected to the workstation, SAC socket and you have declared a SAC950 in the Setup > Configuration menu. Review the Setup > Configuration parameters and/or connect the correct model of sample changer to the SAC socket of the workstation.</p>

6. Maintenance

Cleaning

The sample changer requires minimum maintenance. The exterior surface should be cleaned regularly with tepid water or a mild detergent and wiped dry with a soft damp cloth.

Transporting the instrument

Always use the packaging supplied by the manufacturer.

IMPORTANT!

Remove the tray before transporting the instrument.

Never pick-up or carry the instrument by the electrode head holder or by the arm.

Servicing

DO NOT ATTEMPT TO SERVICE THIS PRODUCT YOURSELF, except as noted in this User's Guide.

To ensure optimum maintenance, it is recommended to have the instrument serviced at least once a year by a qualified Radiometer Analytical service technician. Please contact your service representative at the address given below:

RADIOMETER ANALYTICAL SAS
72, rue d'Alsace
69627 Villeurbanne CEDEX - France
Tel.: +33 (0) 4 78 03 38 38
Fax: +33 (0) 4 78 03 38 27
E-mail: radiometer@nalytical.com

or your local service representative:

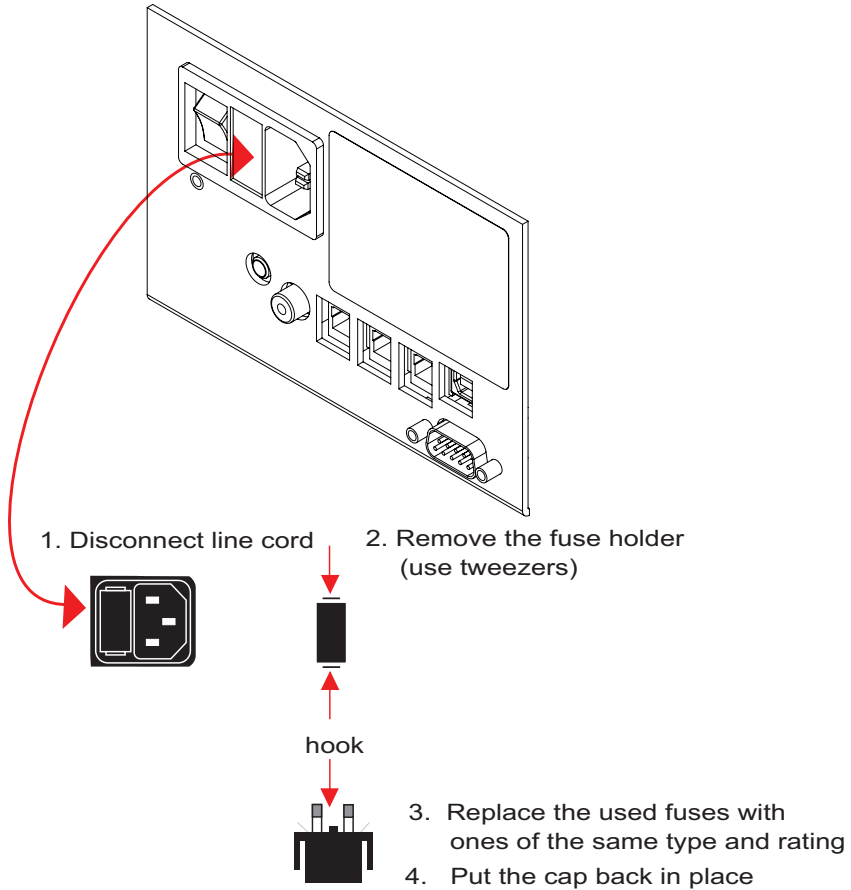
Replacing the fuses

For continued protection replace the fuse with one of a high interrupting capacity, same type and rating:

2 x fuses, slow blow, T2.5A L250V (5 x 20 mm), part no. X16T017.


To replace the fuses:

Locate the fuse housing on the rear panel of the sample changer.

Adjusting the beaker detector position

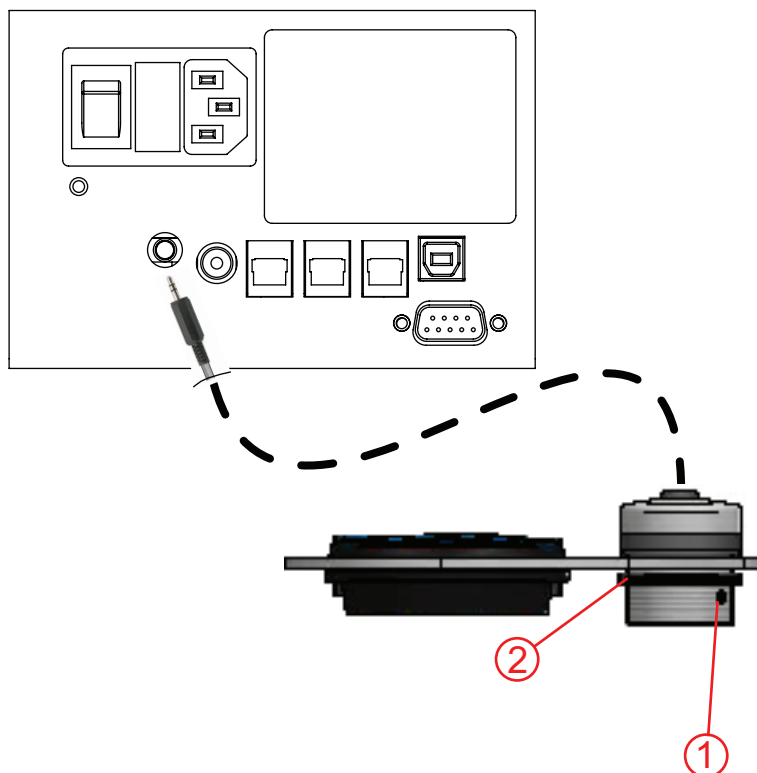1. Switch off the sample changer
2. Unplug the beaker detector.


3. Switch on the sample changer. Wait for the end of instrument initialization. The electrode/beaker detector holder is now above the Park beaker at the upmost position (Service position).
4. Pour a 10 mm height of water in the Park beaker.
5. Connect the beaker detector to the sample changer.

6. Press successively and calmly on keys ,  and  and .

Warning!

Take care to press these keys in the order indicated!

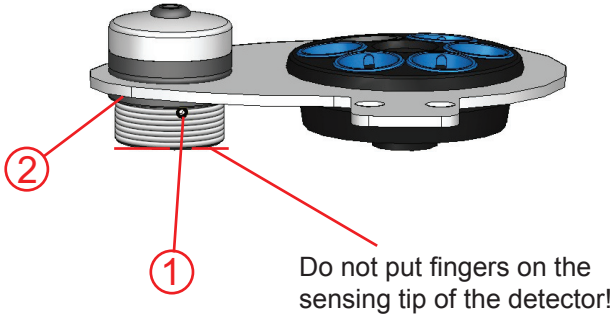7. Check the colour of the status lamp:
 - Blue (difficult to reach): optimal adjustment, detection
 - Purple: correct adjustment, detection
 - Red: wrong adjustment.
8. If relevant, adjust the position of the beaker detector:
 - using the Allen key supplied with the sample changer, screw or unscrew the side screw (1),
 - and/or unscrew the ring (2), rotate the detector and screw ring (2).
 - check the colour of the status lamp that must turn to purple or blue.

Note: In case of problems, it is recommended to turn the detector in order to direct the side screw (1) towards the back of the instrument then to screw/unscrew the side screw (1).

9. Once the adjustment is done, leave the procedure by pressing on

Adjusting the mechanics reference positions

1. Switch off the sample changer.
2. Check the connection of the Ultrasonic transducer plug (3.5 mm jack stereo) to the 'U.SENSOR' socket on the rear panel of the instrument.
3. Place a beaker on **Park** and, in the case of a SAC950, place another beaker on **position 1**. See chapter "Setting up", paragraph "Placing beakers" for more information.
4. Pour a 10 mm height of water in these 2 beakers.
IMPORTANT: The beakers and water must be clean.

Hints:

- To facilitate the adjustment operation, we recommend you to mount the stirrer and one/two electrodes on the electrode head.
 - Pay attention to center correctly the beakers, use blotting paper to eliminate any play of the beaker in its location.
5. Switch on the sample changer. Wait for the end of instrument initialization. The electrode/beaker detector holder is now above the Park beaker at the upmost position (Service position).

6. Press successively and calmly on keys , ,  and  and .

Warning!

Take care to press these keys in the order indicated!

7. For a SAC850:
The electrode head moves from Service down to a position above the Park beaker. The turntable starts rotating by small steps clockwise.

For a SAC950:

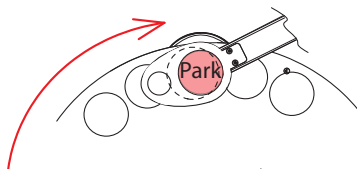
The electrode head moves from Service down to a position above the Park beaker. The arm starts rotating by small steps anti-clockwise.

In both cases, during this phase, evaluate when the stirrer and electrodes are accurately centered above the Park beaker.


When OK (see the figures on the next page for a SAC850), confirm the

position by pressing key .


Too early. Wait a little.

Adjustment correct. Press key .

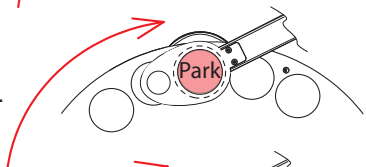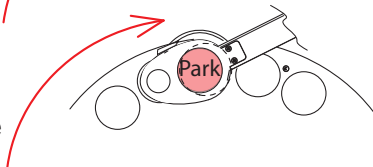Too late. Press key , switch off the sample changer then restart the procedure from step 5.

8. For a SAC850:

The adjustment is complete and the electrodes are dipped into the Park beaker.

For a SAC950:

The electrode head moves to a position above sample beaker no.1. The turntable starts rotating clockwise by small steps.

During this phase, evaluate when the stirrer and electrodes are accurately centered above the Sample beaker no.1.

When OK, confirm the position by pressing key .

The adjustment is complete and the electrodes are dipped into the Park beaker.

7. Accessories

Dynamic rinse module for SACxxx

Dynamic rinse/waste module for SACxxx, part no. X91T107 comprising:

- | | | |
|---|---------|---|
| 1 | X31T166 | Shower head for SACxxx |
| 1 | X16T019 | Pump membrane assembly for SACxxx |
| 1 | X31T154 | Silicon crystal tube, d=4/7 mm, L= 15 m |
| 1 | | OI-Tri, Dynamic rinse module for SACxxx |

Reagent addition module for SACxxx

Reagent addition module for SACxxx, part no. X91T096 comprising:

- | | | |
|---|---------|--|
| 1 | X16T020 | Peristaltic pump assembly for SACxxx |
| 1 | X31T156 | Maprene ® tube, d=4.8/8 mm, L= 0.3 m |
| 1 | X31T154 | Silicon crystal tube, d=4/7 mm, L= 15 m |
| 1 | X31T157 | Addition Tip |
| 1 | | OI-Tri, Reagent addition module for SACxxx |

Cables for SACxxx

Cables for SACxxx:

- | | | |
|---|---------|---|
| 1 | A95X501 | Serial Cable, SubD9/SubD9 |
| 2 | X16T017 | Fuse, slow blow, 2.5A (5x20mm) |
| 2 | A95A126 | Cable, BNC/1m/BNC with Adapter BNC f/f |
| 1 | A95A127 | Cable, Banana-m/1m/Banana-m with Adapter Banana f/f |
| 1 | A95A128 | Cable, CINCH-f/1m/CINCH-m |

Turntables (SAC850 and SAC950)

Kit 25-position turntable, SAC850, 50 ml glass low form, 22-45 ml & 8-25 ml PP beakers, X91T080 comprising:

- 1 25-position turntable, beakers d= 44 mm
- 1 904-489 Titration Vessel PP, 22-45 ml (50 pcs)

Kit 22-positions, turntable, SAC850, glass 100ml tall form beakers, X91T082 comprising:

- 1 22-position turntable, beakers d= 48 mm

Kit 20-position turntable, SAC850, glass 150ml low form beakers, X91T105 comprising:

- 1 20-position turntable, beakers d= 62 mm

Kit 20-position turntable, SAC850, 40-100 ml PP beakers, X91T084 comprising:

- 1 20-position turntable, beakers d= 60 mm
- 1 904-490 Titration Vessel PP, 40-100 ml (50 pcs)

Kit 20-position turntable, SAC850, GOSSELIN 180 ml PP beakers, X91T086 comprising:

- 1 20-position turntable, beakers d= 54 mm

Kit 18-position turntable, SAC850, glass 250ml tall form beakers, X91T088 comprising:

- 1 18-position turntable, beakers d= 60 mm

Kit 15-position turntable, SAC850, glass 250ml low form beakers, X91T090 comprising:

- 1 15-position turntable, beakers d= 70 mm

Kit 15-position turntable, SAC850, glass 400ml tall form beakers, X91T092 comprising:

- 1 15-position turntable, beakers d= 70 mm

Turntables (SAC950 only)

Kit 70-position turntable, SAC9xx, 50 ml glass low form, 22-45 ml & 8-25 ml PP beakers, X91T081 comprising:

- 1 70-position turntable, beakers d= 44 mm
- 2 904-489 Titration Vessel PP, 22-45 ml (50 pcs)

Kit 42-position turntable, SAC9xx, 100 ml glass tall form beakers, X91T083 comprising:

- 1 42-position turntable, beakers d= 48 mm

Kit 36-position turntable, SAC9xx, glass 150ml low form beakers, X91T106 comprising:

- 1 36-position turntable, beakers d= 62 mm

Kit 36-position turntable, SAC9xx, 150 ml glass tall form, 40-100 ml PP beakers, X91T085 comprising:

- 1 36-position turntable, beakers d= 60 mm
- 1 904-490 Titration Vessel PP, 40-100 ml (50 pcs)

Kit 36-position turntable, SAC9xx, GOSSELIN 180 ml PP beakers, X91T087 comprising:

- 1 36-position turntable, beakers d= 54 mm

Kit 30-position turntable, SAC9xx, 250 ml glass tall form beakers, X91T089 comprising:

- 1 30-position turntable, beakers d= 60 mm

Kit 25-position turntable, SAC9xx, 250 ml glass low form beakers, X91T091 comprising:

- 1 25-position turntable, beakers d= 70 mm

Kit 25-position turntable, SAC9xx, 400 ml glass tall form beakers, X91T093 comprising:

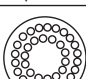
- 1 25-position turntable, beakers d= 70 mm

Turntables and beakers (SAC850 and SAC950)

Beaker part no.	Volume min./max.	Turntable kit to be used	Dynamic rinsing - Shower head skirt	Propeller size
400 ml tall form * d = 70 mm	300 - 360 ml	 X91T092 15 positions		18 mm 
250 ml low form * d = 70 mm	130 - 225 ml	 X91T090 15 positions		18 mm 
250 ml tall form * d = 60 mm	180 - 225 ml	 X91T088 18 positions		18 mm 
180 ml Gosselin, d = 54 mm, PP, X31V005 (50pcs)	100 - 160 ml	 X91T086 20 positions		18 mm 
40 - 100 ml PP 904-490 (50 pcs) d = 60 mm	55 - 90 ml	 X91T084 20 positions		18 mm 
150 ml low form * d = 62 mm	80 - 135 ml	 X91T105 20 positions		18 mm 
100 ml tall form * d = 48 mm	50 - 90 ml	 X91T082 22 positions		18 mm 
50 ml low form * d = 44 mm	30 - 45 ml	 X91T080 25 positions		18 mm 
904-489 (50 pcs) d = 44 mm	22 - 45 ml			10 mm 
904-488 (50 pcs) d = 44 mm	10 - 22 ml			

* Standard glass beakers specified in accordance with DIN12331 ISO3819.

Turntables and beakers (SAC950 only)

Beaker part no.	Volume min./max.	Turntable kit to be used	Dynamic rinsing - Shower head skirt	Propeller size
400 ml tall form * d = 70 mm	300 - 360 ml	 X91T093 25 positions		18 mm 
250 ml low form * d = 70 mm	130 - 225 ml	 X91T091 25 positions		18 mm 
250 ml tall form * d = 60 mm	180 - 225 ml	 X91T089 30 positions		18 mm 
180 ml Gosselin, d = 54 mm, PP, X31V005 (50pcs)	100 - 160 ml	 X91T087 36 positions		18 mm 
40 - 100 ml PP 904-490 (50 pcs) d = 60 mm	55 - 90 ml	 X91T085 36 positions		18 mm 
150 ml low form * d = 62 mm	80 - 135 ml	 X91T106 36 positions		18 mm 
100 ml tall form * d = 48 mm	50 - 90 ml	 X91T083 42 positions		18 mm 
50 ml low form * d = 44 mm	30 - 45 ml	 X91T081 70 positions		18 mm 
904-489 (50 pcs) d = 44 mm	22 - 45 ml			10 mm 
904-488 (50 pcs) d = 44 mm	10 - 22 ml			10 mm 

* Standard glass beakers specified in accordance with DIN12331 ISO3819.

Beakers and beaker covers

Beaker type Ø x H (mm)	Volume min./max.	Detection min. height **	Beaker Part no.	Beaker cover***
250 ml tall form * 60 x 120	180 - 225 ml	10 mm (30 ml)		X31T129 (36 pcs.)
180 ml, Gosselin, PP, 54 x 102	100 - 160 ml	10 mm (25 ml)	X31V005 (50 pcs.)	X31T162 (36 pcs.)
150 ml low form * 62 x 81	80 - 135 ml	10 mm (30 ml)		X31T129 (36 pcs.)
40-100 ml, PP 60 x 80	55 - 90 ml	10 mm (30 ml)	904-490 (50 pcs.)	X31T129 (36 pcs.)
100 ml tall form * 48 x 80	50 - 90 ml	10 mm (20 ml)		X31T161 (42 pcs.)
50 ml low form * 42 x 60	30 - 45 ml	10 mm (15 ml)		X31T160 (70 pcs.)
22-45 ml, PP 44 x 70	22 - 45 ml	10 mm (15 ml)	904-489 (50 pcs.)	X31T160 (70 pcs.)
8-25 ml, PP 44 x 70	10 - 22 ml	10 mm (15 ml)	904-488 (50 pcs.)	X31T160 (70 pcs.)

* Standard glass beakers specified in accordance with DIN12331 ISO3819.

** When starting an analysis, if you have less liquid sample than this specified value, then the beaker automatic detection may not work.

When working with a workstation, you must clear the option "Beaker detection" in TitraMaster 85 (refer to the TitraMaster 85, on-line help, topics "Editing an application - Configuration").

You have also the possibility, with a SAC950 and the Beaker cover module, to cover beakers with the suitable metal covers so that the automatic control of the beakers functions.

*** Beaker metal covers that can be uncovered using the beaker cover module, part no. X16T016

Adapters for Bayonet Electrode Head

Bayonet adapters & Delivery tubings for SACxxx:

- | | | |
|---|---------|--|
| 1 | X31T082 | Bayonet Adapter 4xD4362 |
| 1 | X31T015 | Bayonet Adapter, L=120mm FX Elect. |
| 2 | X31T018 | Bayonet Adapter, 12/7,5 mm |
| 2 | 842-302 | Delivery Tubing Long, 3,25m, for burette stand \geq 5 ml |

Spare parts and consumables not included in kits

- | | |
|---------|--|
| X31T130 | Set of propeller, 10 & 18 mm, for 24Vdc Stirrer |
| 842-151 | Delivery Tubing Long, 160 cm, for burette stand $<$ 5 ml |
| A94L336 | CL336 Cable FX/COAX/3m/MAB6 |
| 842-152 | Delivery Tubing Long, 160 cm, B505/510/525 |

8. Technical specifications

Functions

Easy removable turntables, automatic identification of turntable model by RFID:

- SAC850: 1 radius
- SAC950: 1, 2 or 3 radii

Dip rinse and/or reconditioning: 0 to 2 positions.

Service position (electrodes above the park beaker).

Automatic detection of empty position by ultrasonic transducer.

Stirring: magnetic or propeller stirring. Magnetic stirring if propeller stirrer not connected. 0 to 1100 rpm by steps of 50 rpm.

Built-in keypad: lift electrode head up/down, move to next/previous beaker, start/stop dynamic shower rinse, start/stop emptying to waste, start/stop adding a reagent, stop action.

Embedded programmable sample preparation sequences.

Unused electrodes and delivery tips storage: 4 positions Ø16 mm, on the top of the sample changer arm.

Optional modules

Dynamic shower rinse & Beaker emptying: Shower rinse and emptying operations controlled by 2 PTFE membrane/PE body pumps, available in any beaker size. Flow rate: up to 600 ml/min. Pumps can be controlled manually and individually through keypad.

Addition of a reagent with a peristaltic pump to prepare sample. The pump can be controlled manually through keypad. The pump can be programmed individually as part of sample preparation sequence.

Beaker cover option (SAC950 only): Remove beaker covers using an electro-magnet. Covers are disposed beside the sample changer in a dedicated location once removed.

Inputs/outputs

USB-B (Slave) connection to computer for sample preparation programming sequence.

RS232C serial connection to workstation unit.

TTL I/O. Start external peripherals (RJ11 socket). Wait for external signal to continue the sample preparation sequence.

Electro-magnet input for beaker cover option (RJ11 socket).

Propeller stirrer output (RJ11 socket).

Temperature input (CINCH socket) for use as input in a sample preparation sequence.

Ultrasonic transducer input for beaker detection module (Jack socket 3.5 mm).

Casing, keyboard

PETP – PVC and/or PE

Protected against projections – For use in a laboratory environment

Dimensions (H x W x D)

70 x 58 x 68 cm.

Weight

27 kg (SAC950), 26 kg (SAC850).

Power requirements

100-240 Vac - 50-60 Hz. 76 VA

Fuses: 2 mains fuses : slow blow T2.5A L250V (230 V)

Environmental operating conditions

Interior use only

5 to 30°C temperature.

20 to 80% relative humidity.

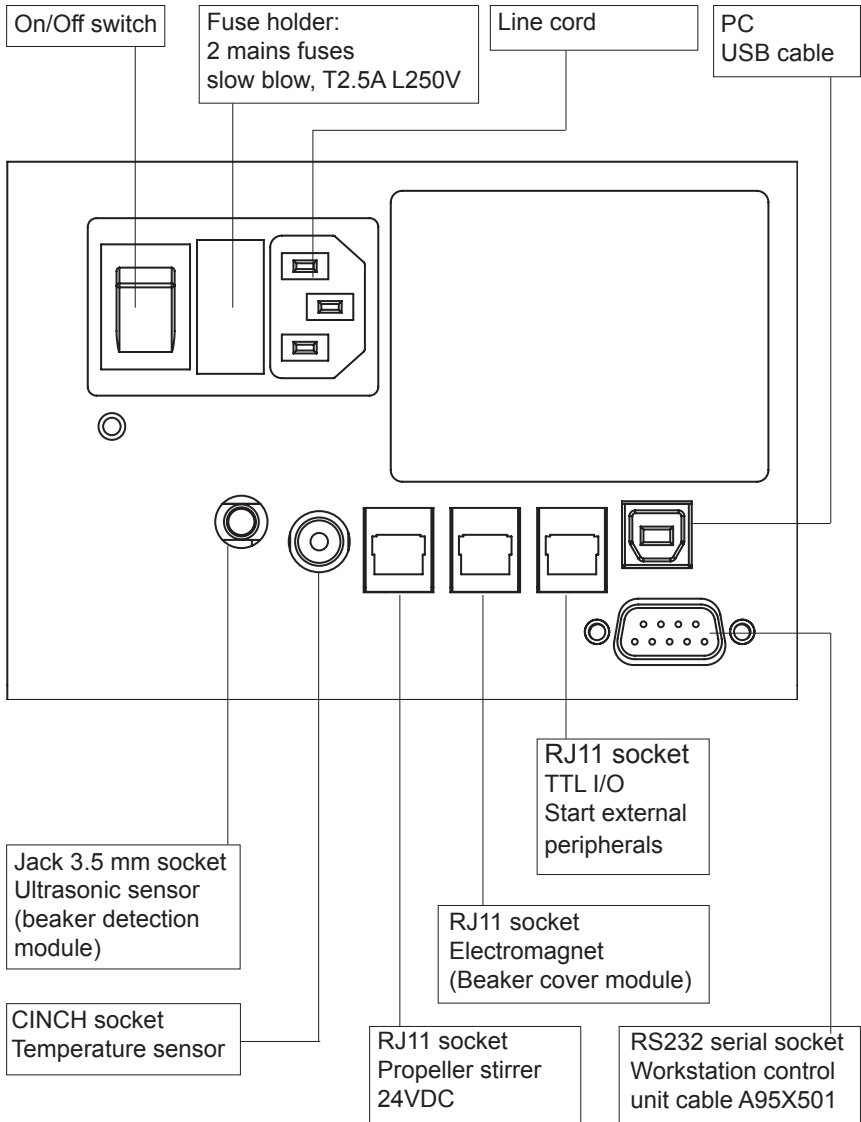
Altitude \leq 2000 m.

Variations of the mains voltage do not exceed \pm 10% of the nominal value.

Level of pollution: 2.

Transitory overvoltage: level II.

Rear panel

International Standards

Radio and Telecommunication Directive (99/5/EEC)

Standards :

EN300330-1

en301489

EMC Directive (2004/108/EEC)

The sample changer complies with following standards:
Class A equipment for laboratory use, according to the standard
EN 61326-1.

Emission:

EN 55022

EN 61000-3-2

EN 61000-3-3

Immunity:

EN 61000-4-2

EN 61000-4-3

EN 61000-4-4

EN 61000-4-5

EN 61000-4-6

EN 61000-4-11

Low Voltage Directive (2006/95/EEC)

The sample changer complies with the following standard:
Reference standard: EN 61010-1.

The sample changer complies with the following standards:

UL 61010-1

CAN / CSA C22 2 N° 61010-1

The sample changer complies with part 15 of the FCC Rules. Operation is subject to the following two conditions:

- (1) This device may not cause harmful interference, and
- (2) this device must accept any interference received, including interference that may cause undesired operation.

See Chapter 1, "Compliance with FCC rules, part 15 - Information to the user".

The sample changer complies with IC Rules (FCC Rules applied to Canada)

European RoHS directive 2002/95/EC

The following products:

Part no.	Description
R41T043	SAC850, Sample changer
R41T044	SAC950, Sample changer
X91T095	Dynamic rinse module
X91T096	Reagent addition module
X16T016	Beaker cover module
X61T007	Propeller Stirrer

comply with the European RoHS directive 2002/95/EC.

Electrical and electronic equipment marked with this symbol may not be disposed of in European public disposal systems after 13 August of 2005. In conformity with European local and national regulations (EU Directive 2002/96/EC), European electrical equipment users must now return old or end-of life equipment to the Producer for disposal at no charge to the user.

Note: For equipment produced by "Radiometer Analytical", please talk to your local contact.

Important document. Retain with product records.

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