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
For current price information, technical support and ordering assistance, contact the Hach office or distributor serving your area.

In the United States, contact:

HACH COMPANY
P.O. Box 389
Loveland, Colorado 80539-0389
U.S.A.
Telephone: 800-227-4224
Fax: 970-669-2932
E-mail: orders@hach.com
www.hach.com

In Europe, the Middle East, and
Mediterranean Africa, contact:

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
DOC02.98.00729 Sept.05 Printed by Hach Lange Germany All rights reserved

Hach DataTrans

Data Collection software for spectrophotometers

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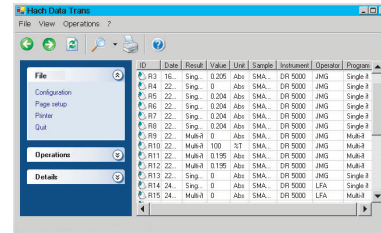Introduction

Hach Data Trans is a Windows-based PC program dedicated for use with the DR 5000 / DR 2800 / DR 4800 spectrophotometer.

With the Hach Data Trans, the method is edited and started from the instrument. Measurements are collected, displayed and saved by the software.

Hach Data Trans is a powerful data archiving tool, which enables you to access your stored results and curves in a flash via user-programmable requests. You can consult the database and filter and sort the archived results by date, program type etc. providing the information you need at a glance.

Minimum configuration recommended

Operating system: W2000 Pro (SP2) or XP Pro (SP2).

Microsoft® .NET Framework Version 1.1 and Microsoft SQL Server Desktop Engine (*).

- Pentium 1 GHz processor or equivalent, 256 MB RAM,
- SVGA 1024 x 768 screen resolution,
- 1 available USB 1.1 or USB 2.0 port,
- 1 CD-ROM drive,
- 1 PS2 or USB mouse type,
- 1 Windows compatible printer (optional),
- Hard disk, 150 MB of free space for software installation, 300 MB required for using properly the software with your database. The result database size is limited to 2 GB.

(* If .NET Framework Version 1.1 and/or Microsoft SQL Server Desktop Engine are not installed on your PC, you can ask the Hach Data Trans setup software to install them.

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Software installation

Before starting installation

Refer to system requirements and make sure that your PC complies with the minimum configuration recommended (see previous chapter).

Installation procedure

- Close all applications.
- Insert the CD-ROM into the CD-ROM drive. Installation starts automatically (AUTORUN). If it does not, use your Windows Explorer to locate Setup.exe and run it.
- You are prompted to install Microsoft® .NET Framework Version 1.1. Click **Yes** then accept the license agreement for installing .NET Framework Version 1.1 or click **No** if .NET Framework Version 1.1 is already installed on your PC.
- The Hach Data Trans Setup window is displayed. The Setup software checks if the Microsoft SQL Server Desktop Engine (MSDE) is installed and updated on your PC. Click "Hach Data Trans Prerequisites" to install/update MSDE if prompted to do so, otherwise click "Hach Data Trans". If you click **Exit**, you quit the installation program and Hach Data Trans is not installed.
- At the end of MSDE installation, click "Hach Data Trans".
- Hach Data Trans installation starts with the Welcome dialogue box. Click the **Next** button.
- Accept the installation directory proposed by default clicking **Next** or change it clicking the **Browse** button.
- Review and confirm your installation settings. Click **Next**.
- The installation program creates the destination directory, transfers the files into the directory and creates an icon for starting the software in the "Start > All Programs > Hach Data Trans" folder.

10. Click the **Close** button at the end of installation.
11. Click **Exit** in the Hach Data Trans Setup window.

The software is installed in the language used by Windows (English, French, German, Spanish or Italian). The software is installed in English for all other languages.

If you change your Windows operating system on your PC, you must uninstall then reinstall Hach Data Trans.

It is highly recommended to use Microsoft Windows Update to check for any critical updates to the .Net Framework product.

Getting help

Click  to gain access to the Hach Data Trans online help. You can also press F1 or select the command ? > **Contents** to display the contents menu of the help file. From here, select the item of interest or search for a certain topic.

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Starting up

Connect the USB port socket on the rear panel of the instrument (for example, the DR 5000) to an USB port on the PC.

If necessary, connect a printer to a USB port on the PC.

Connect the instrument to the mains using the line cord supplied.

Switch on the instrument (**O/I** switch on the rear panel set to **I**).

Install the instrument driver:

- Start windows on the PC. A new hardware is detected automatically.
- In the Found New Hardware Wizard window, select "Install from a list or specific location" and click **Next**.
- Insert the Hach Data Trans CD-ROM supplied into a CD drive.
- Under "Search for the best driver in these locations", select "Include this location in the search" and browse the Hach Data Trans CD to select the driver installation directory: \USB Drivers.
- Click **OK**.
*You may have a Windows warning message telling that your hardware has not passed a Windows Logo Testing. Do not pay attention to this message and click **Continue anyway**.*
- Click **Finish** at end of installation.

Start Hach Data Trans, select **Start > Programs > Hach Data Trans > Hach Data Trans**.

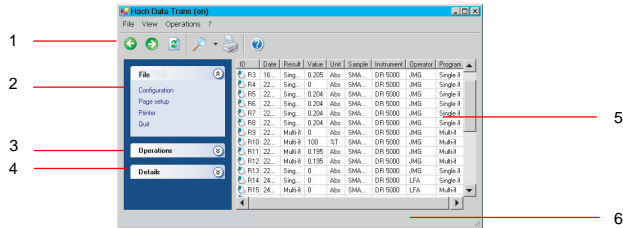
Hach Data Trans scans the PC's port(s) and detects the instrument connected and on-line.

The Main window is displayed.

You are now ready to run a program and collect measurements!

Main window

- Function toolbar:** comprises icons (see next page). To display or hide the tool bar, use the **View > Tool Bar** command.
- File menu:** gives access to a series of commands:
Configuration: to configure printouts and define the data separation character when exporting results to a LIMS.
Page setup: to open the Page setup window where you select a printout format.
Printer: to open the Printout window where you select a printer.
Quit: to quit Hach Data Trans.
- Operations menu:** gives access to a series of commands:
Search: to consult the database and filter and sort the archived results.
Export to LIMS: to export results to a text format file ("txt" file).
Export to Excel: to export results to an Excel format file ("xls" file).
Preview: to preview a printout before starting printing.
Print: to start printing the results currently displayed.
Show raw data: display the coordinates of the measurement points.
Show curve: display the curve of the selected scan method.
- Details menu:** Displays information about the results currently displayed (description of the request criteria selected, etc.).
- Results area:** This area displays all the results of the database or all the results of the database which satisfy the search criteria of a selected request.
- Status bar:** displays information about the software working status. To display or hide the status bar, use the **Display > Status Bar** command.

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Icons

Previous: icon of the Main window. Displays the last viewed results screen.

Next: icon of the Main window. Displays the next viewed results screen.

Refresh list: icon of the Main window. Runs the last search request and refreshes the result screen.

Search: icon of the Main window. Selects and runs a request or displays the Search results window. A request is used to filter and sort the results of the database in order to retrieve those of interest to you.

Print: icon of the Main window. Prints the results currently displayed.

Help: Gives access to online help.

Modify: icon of the Search results window. Gives access to the edit request parameters window. When in the edit mode you are able to modify the searching criteria of a selected request.

Rename: icon of the Search results window. To change the name of a request.

Create: icon of the Search results window. To create a request.

Delete: icon of the Search results window. To delete a request.

Uninstalling

1. Select "Start > Settings > Control Panel > Add/Remove Programs" command in Windows.
2. Select "Hach Data Trans" and click **Add/Remove**.
3. The Hach Data Trans Setup window is displayed. Click "Hach Data Trans". If you click **Exit**, you quit the uninstallation program and Hach Data Trans is not uninstalled.
4. Select "Remove Hach Data Trans" in the next window and click **Finish**. All the files that have been created and copied by the Hach Data Trans setup program are removed. The Hach Data Trans icon placed in the "Start" folder is also removed.
5. At the end of uninstallation, click **Close**.
6. The Hach Data Trans Setup window is displayed. You can now uninstall Microsoft SQL Server Desktop Engine (MSDE) by clicking "Hach Data Trans Prerequisites", otherwise click **Exit**.

Click **Exit** if you want to keep MSDE installed; the Hach Data Trans result database is not deleted in this case.

To delete the Hach Data Trans database, uninstall Hach Data Trans (steps 1 to 5 above) then click "Hach Data Trans Prerequisites" to uninstall MSDE.

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