FG7001014

User Instructions

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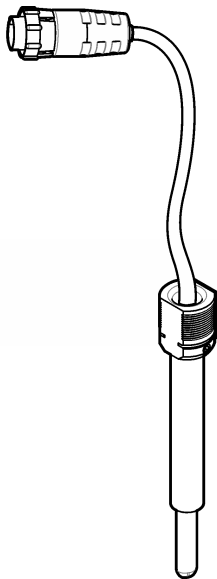
www.hach.com

PT-BR

Perigo

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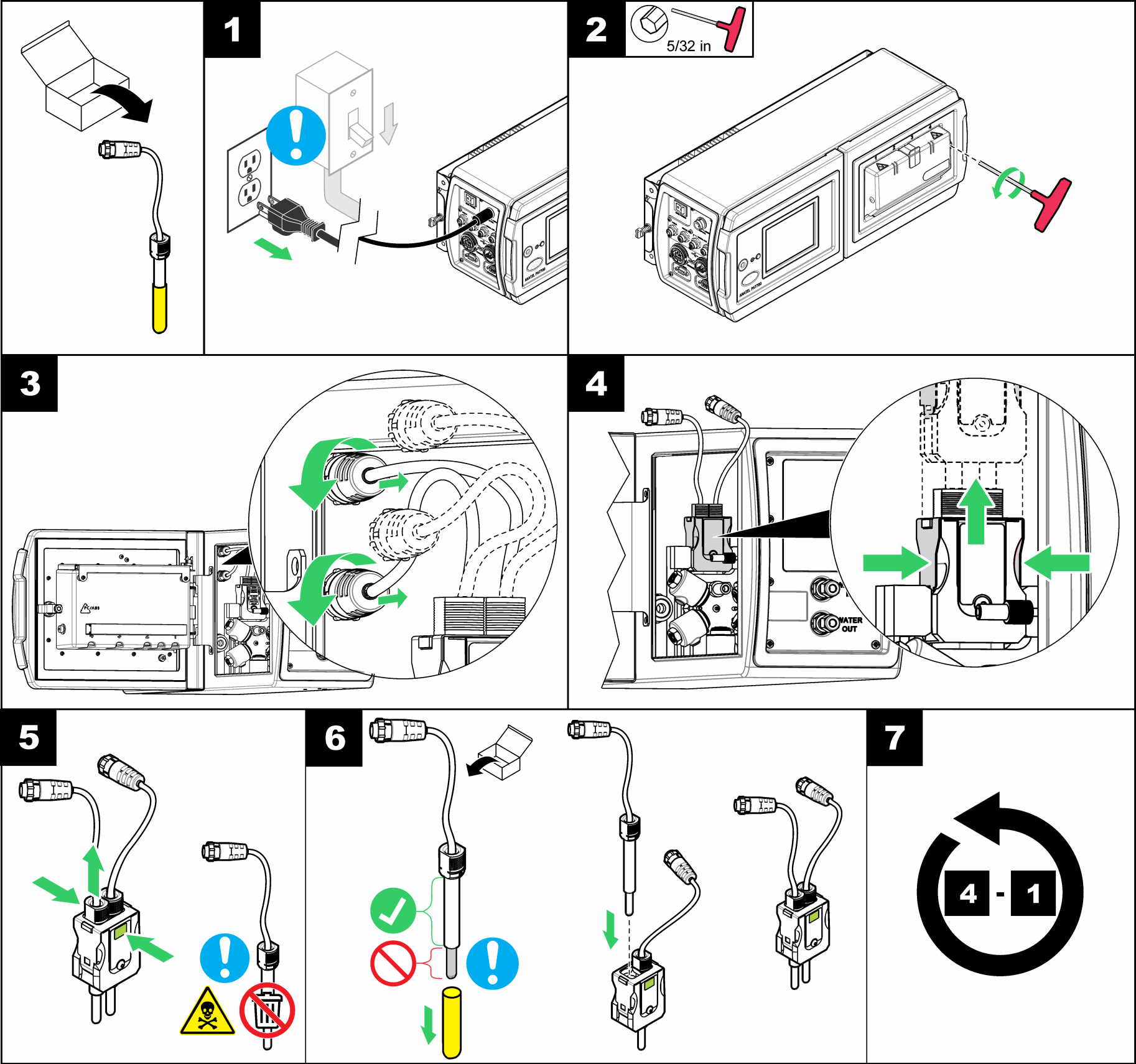
EN

Danger
⚠ **Electrocution hazard.** Remove power from the instrument before this procedure is started.

Caution
⚠ **Ultraviolet (UV) light exposure hazard.** Exposure to UV light can cause eye and skin damage. Protect eyes and skin from direct exposure to UV light.

⚠ **Burn hazard.** Turn off power and allow the instrument to cool before this procedure.

⚠ **Chemical hazard.** UV lamps contain mercury. Consult the correct local regulatory agency for safe handling and current lamp disposal requirements.

⚠ **Chemical hazard.** Never use a vacuum cleaner to clean up broken lamps. All mercury-contaminated debris must be cleaned up as hazardous waste. Obey local regulations and guidelines for the removal and disposal of mercury waste.

Notice
Do not touch the glass surface of the lamp because oil from skin will cause damage. Clean the glass surface of the lamp with isopropyl alcohol and a lint-free cloth as necessary.

- Before power is removed:**
1. Push << to show the sliding menu.
 2. Push **DIAGNOSTICS**, then select the General tab.
 3. Push **REPLACE LAMP**, then select the lamp to replace.
- Select **Run Wizard** for guided lamp replacement.
 - Select **Replace Now** to bypass the wizard.

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