9377700, 9394900

EN: NOTICE! Replace the trays in the instrument at intervals of 6 to 12 months depending on individual tray usage to prevent measurement errors. If frequent error or warning messages show on the display, replace the identified tray.

Items to collect: Cross-slotted (Phillips) screwdriver, cotton tipped applicator, isopropyl alcohol and mild detergent

The SL1000 instrument has four Chemkey trays and the SL250 instrument has one Chemkey tray. The Chemkey trays are all the same. Refer to the illustrations in the *Maintenance* section of the SL250/SL1000 user manual and the steps that follow to replace the Chemkey tray(s).

1. Remove the Chemkey tray(s) from the instrument. 2. Discard the used tray(s) in accordance with local, regional and national regulations. 3. Clean the light shield with water, mild detergent and a cotton tipped applicator. 4. Clean the sample detector and access cover with isopropyl alcohol and a cotton tipped applicator. 5. Clean the Chemkey slot(s) with isopropyl alcohol. Make sure to move the cotton tipped applicator in one direction only. 6. Install the new Chemkey tray(s) in the correct orientation and assemble the instrument.

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P.O. Box 389, Loveland, CO 80539-0389 U.S.A.
(800) 227-4224 (U.S.A. only)